

Acidic Herbicides in Drinking Water Using Xterra Column

Waters Corporation



This is an Application Brief and does not contain a detailed Experimental section.


Abstract

This application brief demonstrates analysis of acidic herbicides in drinking water.

Introduction

The compounds used in this study are –

1. Picloram
2. Chloramben
3. 4-nitrophenol (non-linear above 500)
4. Bentazon (non-linear above 300)
5. 2, 4-D
6. MCPA
7. 2, 4, 5-TP
8. Dichloroprop
9. MCPP
10. Dichlorobenzoic
11. Acifluorfen (non-linear above 300)
12. 2, 6, 5 -TP
13. 2, 4-DB
14. Dinoseb (non-linear above 200)
15. Pentachlorophenol

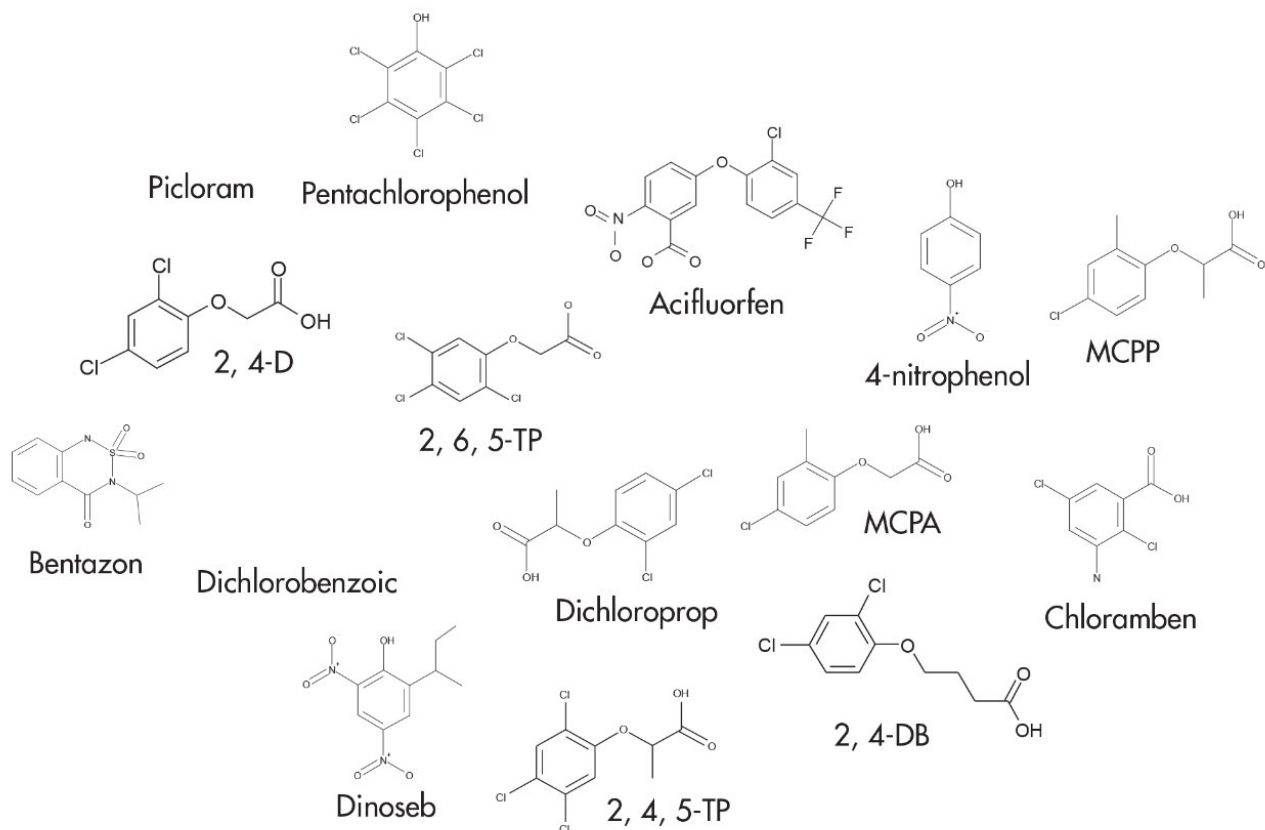

Compounds.

Experimental

Conditions

Column:	Xterra MS C ₁₈ 2.1 x 100 mm, 3.5 μm
Part number:	186000404
Mobile phase A:	15 mM NH ₄ COOH, pH 3.4
Mobile phase B:	ACN
Flow rate:	0.2 mL/min to MS

Injection volume: 20 µL

Detection: MS ESI-


Instrument: Alliance 2695, Micromass ZQ

Gradient


Time (min)	Profile	
	%A	%B
0.0	75	25
9.0	40	60
14.0	40	60
30.0	10	90

Oasis® MAX SPE Method
for Acidic Herbicides
Conditions for Oasis® MAX Cartridge, 6 cc, 150 mg
Part Number 186000369

Sample is first hydrolyzed at pH 12 for 60 min. Then, pH is adjusted to approx. neutral with HCl before SPE.

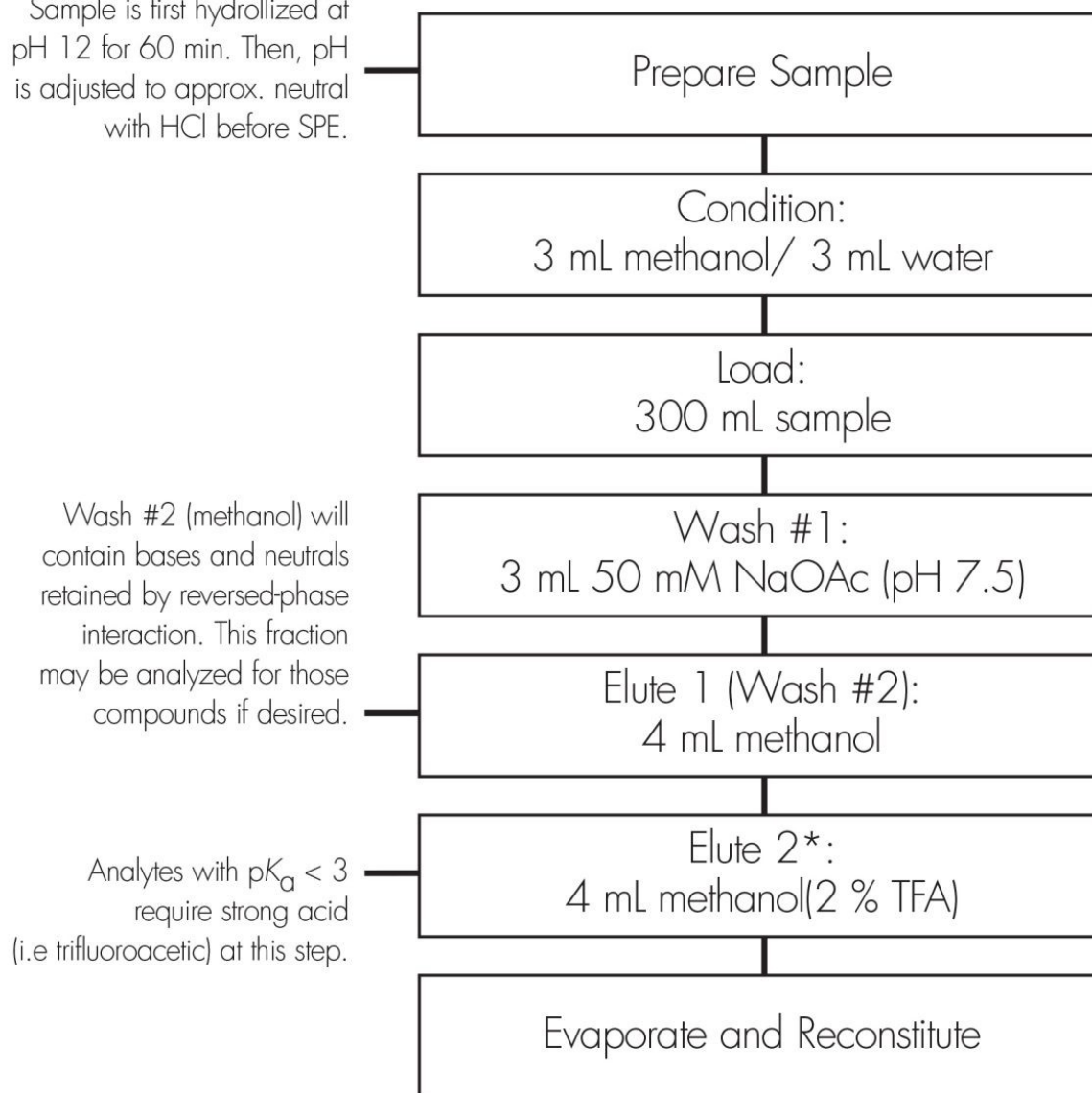

Results and Discussion

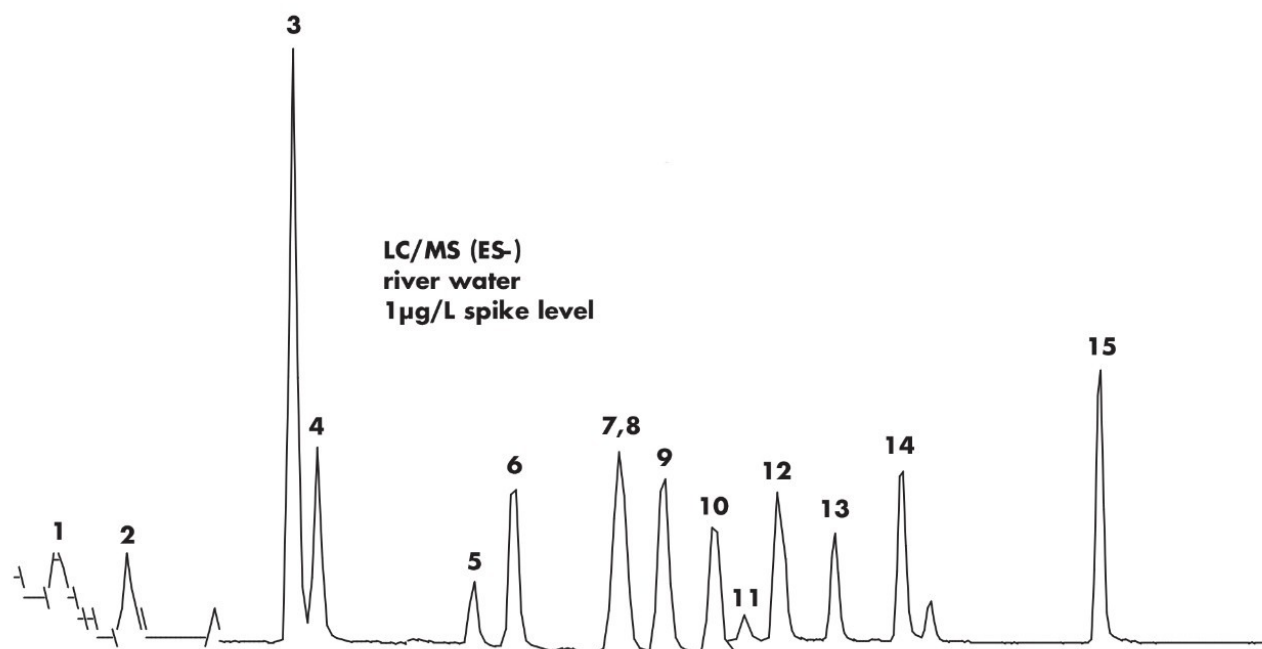

Featured Products

· [Alliance HPLC System <https://www.waters.com/534293>](https://www.waters.com/534293)

WA20738.004, June 2002



© 2021 Waters Corporation. All Rights Reserved.