

Atenolol

Waters Corporation




This is an Application Brief and does not contain a detailed Experimental section.

Abstract

This application brief demonstrates analysis of atenolol.

Introduction

The compound analyzed in this study is atenolol.

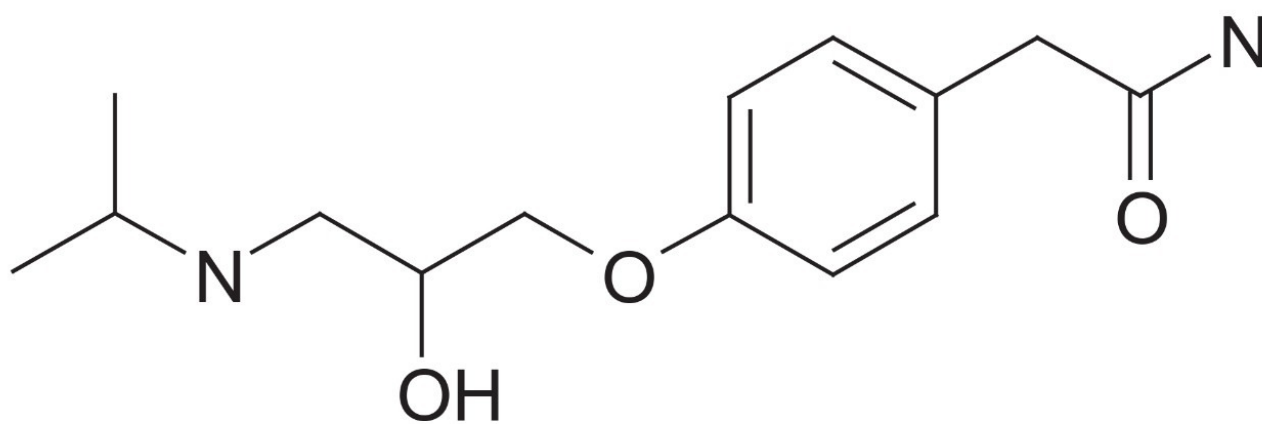

Atenolol


Experimental

Conditions

Column:	Xterra RP ₁₈ 4.6 x 50 mm, 5 μm
Part number:	186000492
Mobile phase:	pH 3.0: H ₂ O/ACN/100 mM NH ₄ COOH, pH 3.0 75:15:10 pH 7.0: H ₂ O/ACN/100 mM NH ₄ HCO ₃ , pH 7.0 80:10:10 pH 10.0 H ₂ O/ACN/100 mM NH ₄ HCO ₃ , pH 10.0 78:12:10

Flow rate:	1.0 mL/min
Injection volume:	5 µL of 250 µg/mL
Temperature:	30 °C
Detection:	UV @ 270 nm
Instrument:	Alliance 2695, 2996 PDA
Mobile Phase pH	USP Tailing
3.0	1.21
7.0	1.38
10.0	1.06

Results and Discussion

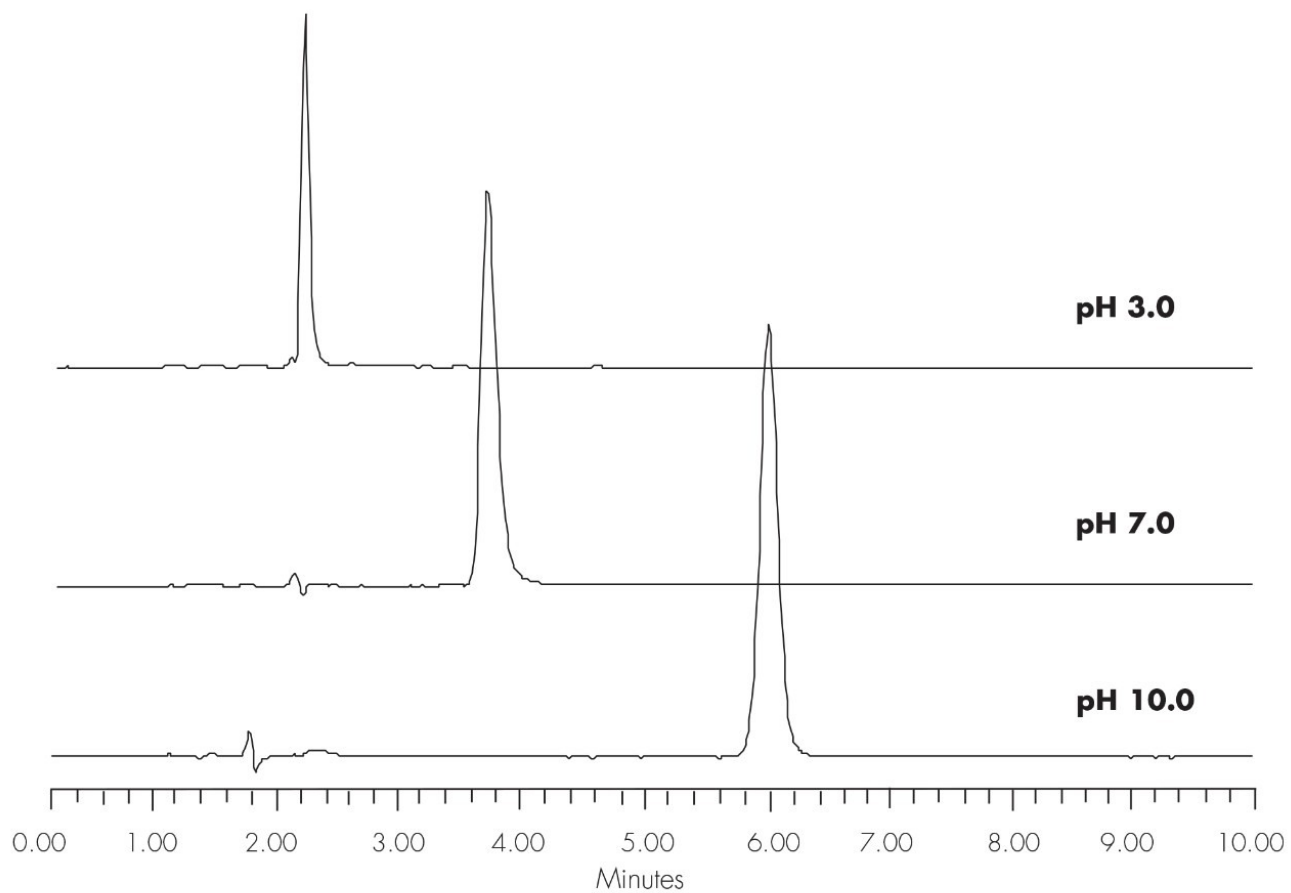

Featured Products

· [Alliance HPLC System <https://www.waters.com/534293>](https://www.waters.com/534293)

WA20738.012, June 2002



© 2021 Waters Corporation. All Rights Reserved.