

Nota applicativa

Warfarin in Rat Plasma

Waters Corporation




This is an Application Brief and does not contain a detailed Experimental section.

Abstract

This application brief highlights the analysis of warfarin in rat plasma using XTerra MS C₁₈ columns.

Introduction

Warfarin in rat plasma has been analyzed in this application brief.

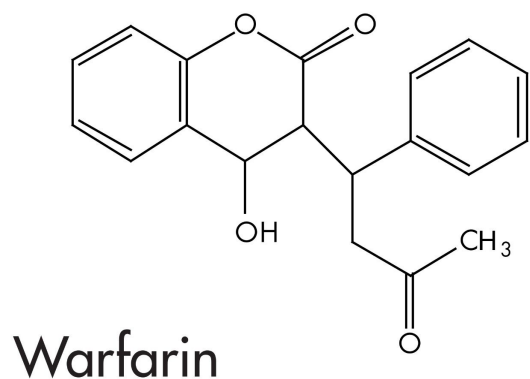

Experimental

HPLC Conditions

Column:	XTerra MS C ₁₈ 2.1 x 30 mm, 3.5 μm (p/n: 186000398)
Mobile phase A:	0.5% HCOOH
Mobile phase B:	ACN
Flow rate:	0.2 mL/min
Isocratic mobile phase composition:	40% A; 60% B
Injection volume:	20 μL
Detection:	MS ESI-

Instrument:	Alliance 2790, Micromass Quattro Ultima
-------------	---

MS Conditions

Ion source:	ESI-
-------------	------

Source temp.:	150 °C
---------------	--------

Gas cell:	1.5×10^{-3} mbar, 30 eV
-----------	----------------------------------


Desolvation temp.:	350 °C
--------------------	--------

Cone gas flow:	150 L/hr
----------------	----------

Drying ga flow:	600 L/hr
-----------------	----------


Cone voltage:	30 V
---------------	------

Oasis® MAX Extraction Method
Oasis® MAX Extraction Plate, 10 mg/ 96-well
Part Number 186000375

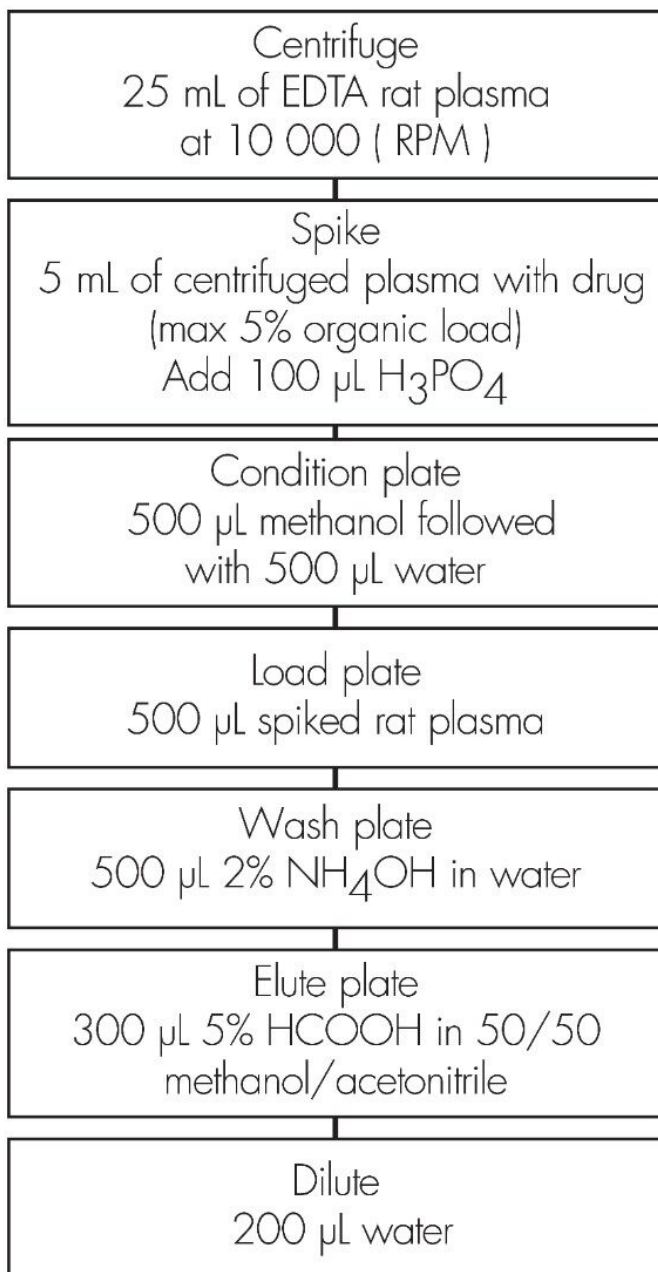

Results and Discussion

CID mass spectra

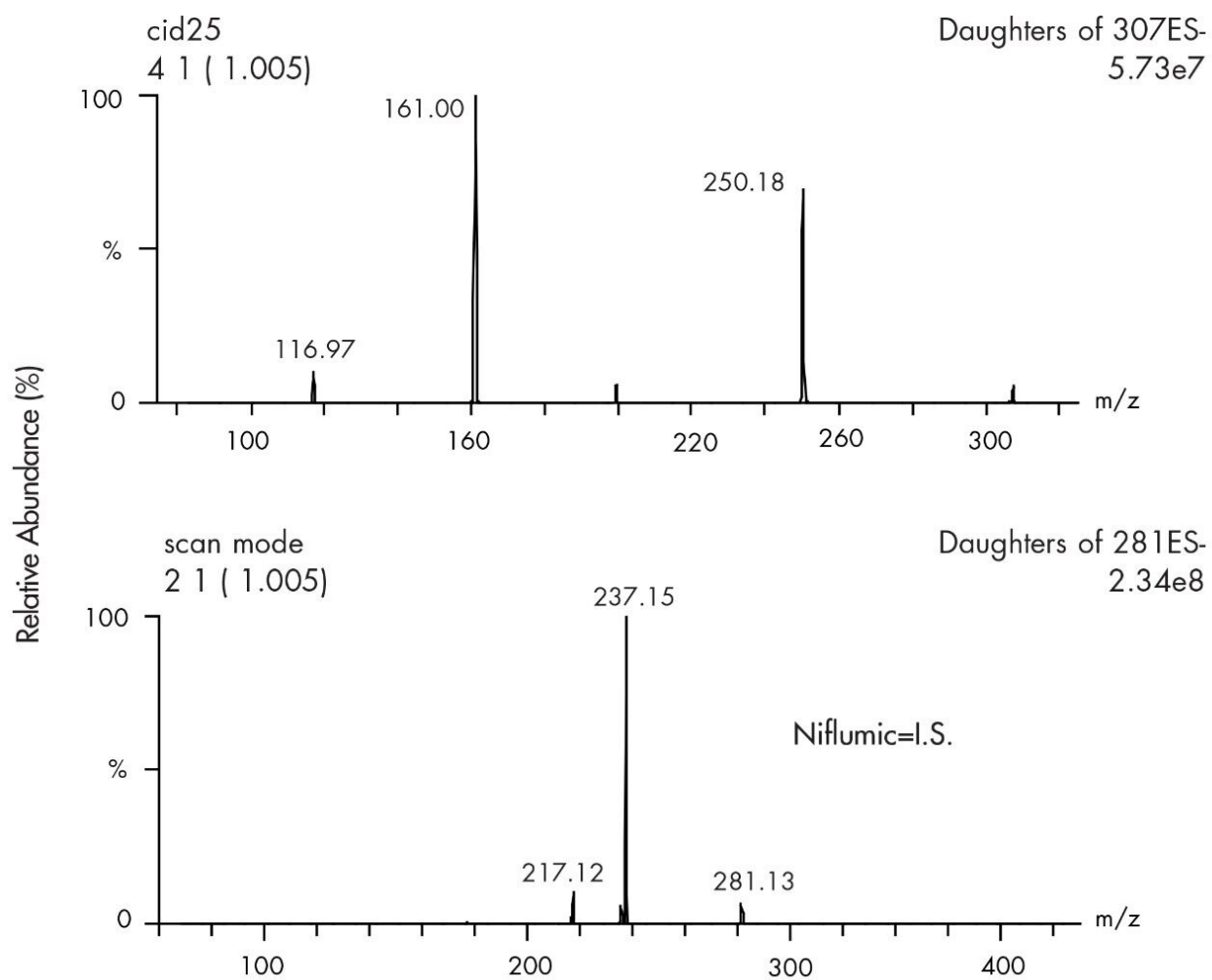

Featured Products

Alliance HPLC <<https://www.waters.com/514248>>

WA20738.118, June 2002

