

On-Line Protocol for Basic Drugs in Cell Culture Lysate Using LC-MS/MS

Waters Corporation



This is an Application Brief and does not contain a detailed Experimental section.


Abstract

This application brief highlights the analysis of basic drugs in cell culture lysate using Oasis HLB on-line


protocol.

Introduction

Basic drugs has been analyzed in cell culture lysate using LC-MS/MS in this application brief.

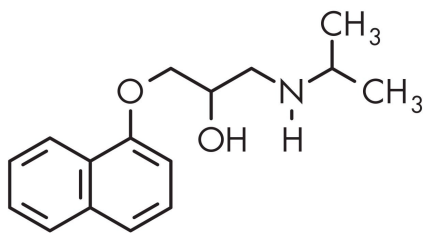

PROPRANOLOL

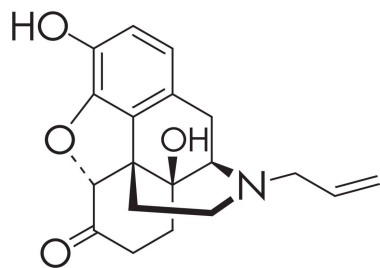

NALOXONE

Experimental

HPLC Method

Column:

XTerra MS C₁₈, 2.1 x 30 mm, 3.5 μm (p/n:
186000398)

Mobile phase A:

Acetonitrile + 0.5% Formic Acid

Mobile phase B:	Water + 0.5% Formic Acid
Flow rate:	1.0 mL
LC ₁ :	Alliance 2690–0.4 mL/min
LC ₂ :	Waters 515–4.0 mL/min
Loading mobile phase:	100 % water
Eluting mobile phase:	1 minute gradient 5% ACN to 95% ACN
Eluting mobile Phase additive:	0.5 % Formic acid
Extraction column temp.:	40 °C
Switching valve:	Rheodyne LabPro 10 ports, 2 position
MS:	Waters Micromass Quattro Ultima
Source:	Electrospray positive
Sorce temp.:	150 °C
Desolvation gas:	600 L/hr
Gas cell:	1.5e ⁻³ mbar
Cone voltage:	20 volts
Collision energy:	20

Gradient

Time (min)	Profile	
	%A	%B
0.0	5	95
0.5	5	95
1.5	95	5
4.40	95	5
4.50	5	95
6.0	5	95

Compound 2 name: Propranolol

Coefficient of Determination: 0.998650

Calibration curve: $-1.52393e^{-5} * x^2 + 0.00390583 * x + 0.000102086$

Response type: Internal Std (Ref 1), Area * (IS Conc. / IS Area)

Curve type: 2nd Order, Origin: Exclude, Weighting: 1/x², Axis trans: None



Compound 3 name: Naloxone

Coefficient of Determination: 0.995263

Calibration curve: $-2.47393e^{-5} * x^2 + 0.00892751 * x + -5.00665e^{-5}$

Response type: Internal Std (Ref 1), Area * (IS Conc. / IS Area)

Curve type: 2nd Order, Origin: Exclude, Weighting: 1/x², Axis trans: None

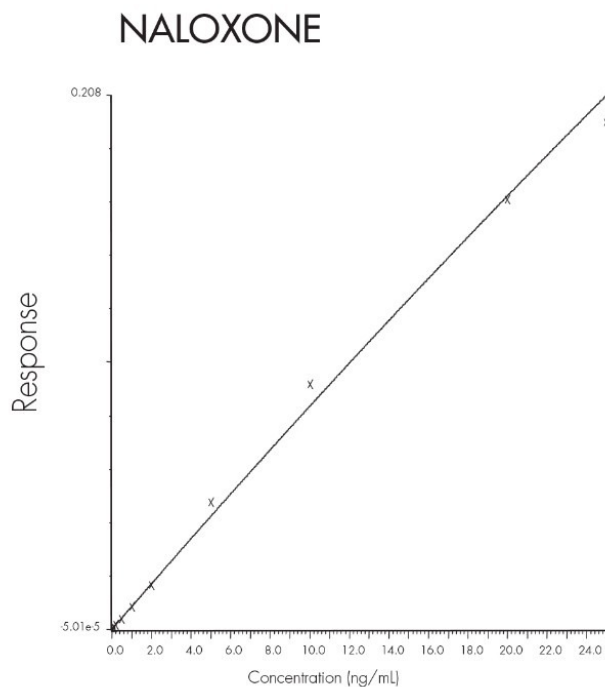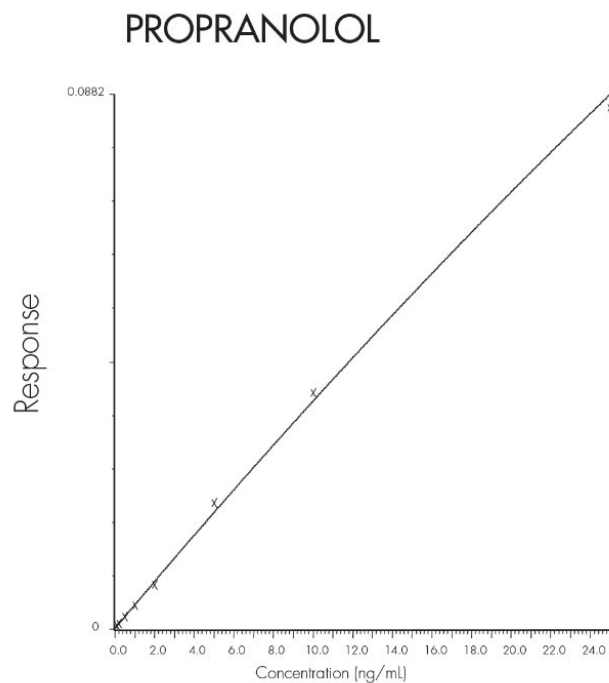


On-Line Protocol

Oasis HLB 2.1 x 20 mm, 25 µm Extraction Column Part Number [186000706](#) <


<https://www.waters.com/nextgen/xg/en/shop/sample-preparation--filtration/186000706-oasis-hlb-online-column-80a-25--m-21-mm-x-20-mm-1-pk.html>>

ON-LINE PROTOCOL

Oasis® HLB 2.1 x 20 mm, 25 µm Extraction Column
Part Number 186000706

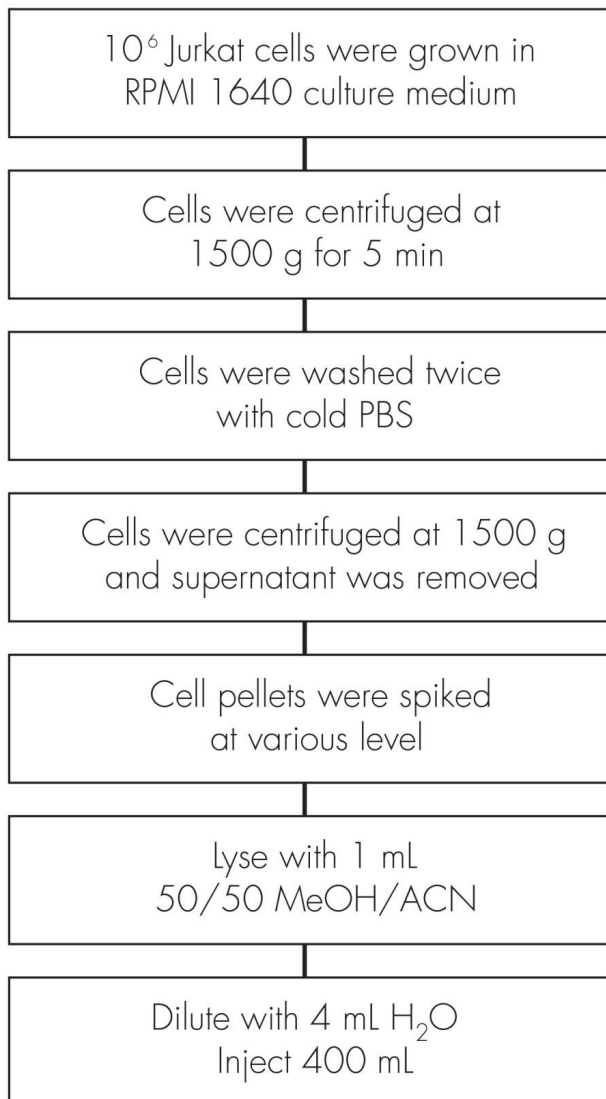


Conc. ng/mL n=6	Average	Standard deviation	RSD%
Propranolol calibration curve			
0.1	0.09	0.05	4.8
0.5	0.49	0.03	5.9
1.0	1.03	0.04	3.9
5.0	4.93	0.26	5.3
10.0	9.76	0.34	3.5
20.0	20.21	0.89	4.4
25.0	25.21	0.69	2.7
Naloxone calibration curve			
0.1	0.10	0.04	4.2
0.5	0.48	0.02	5.0
1.0	0.99	0.06	6.8
5.0	5.25	0.25	4.8
10.0	10.43	0.41	3.9
20.0	19.98	0.54	2.7
25.0	24.44	1.12	4.6

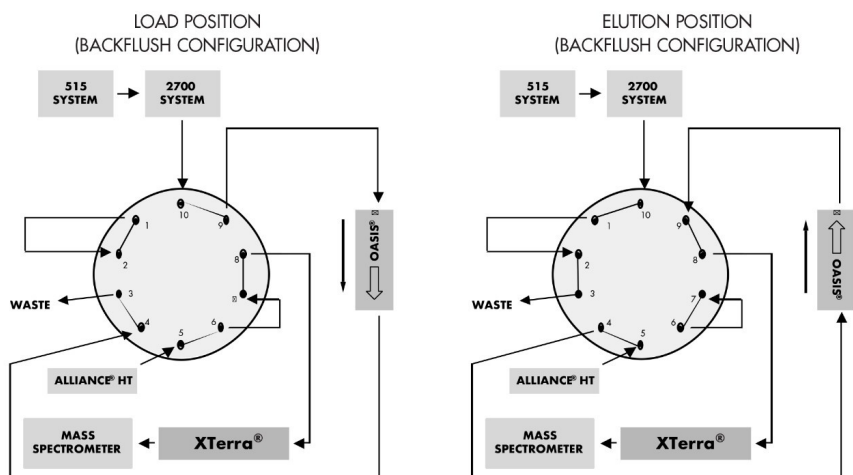


Results and Discussion

CID Mass Spectra

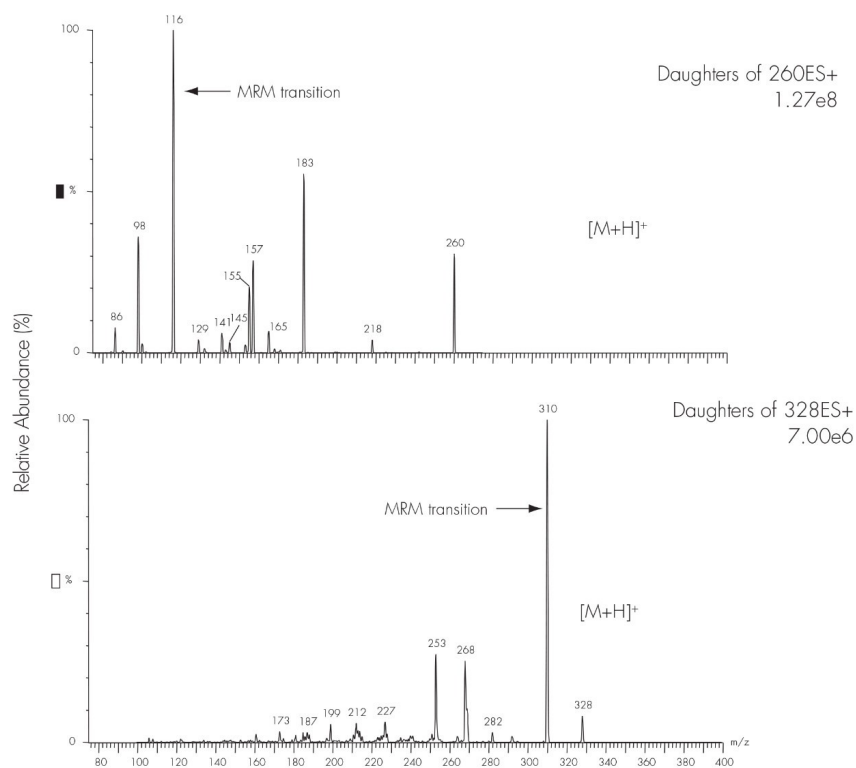


50/50 MEOH/Acetonitrile Cell Lysing

PROPRANOLOL AT 0.1 ng/mL AND 25 ng/mL

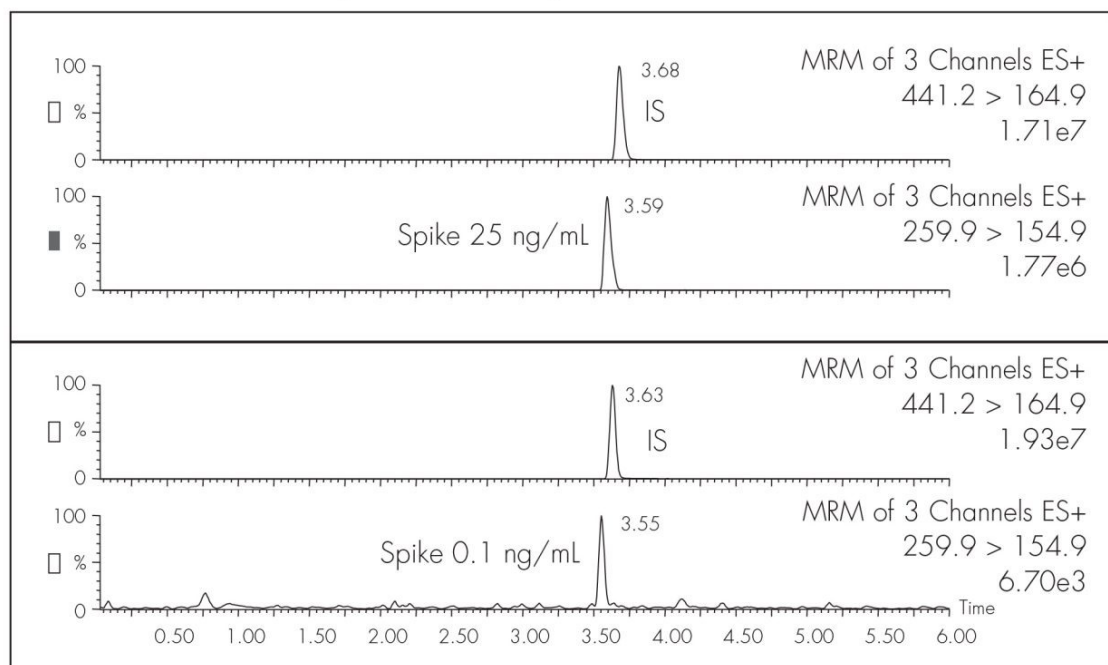

NALOXONE AT 0.1 ng/mL AND 25 ng/mL

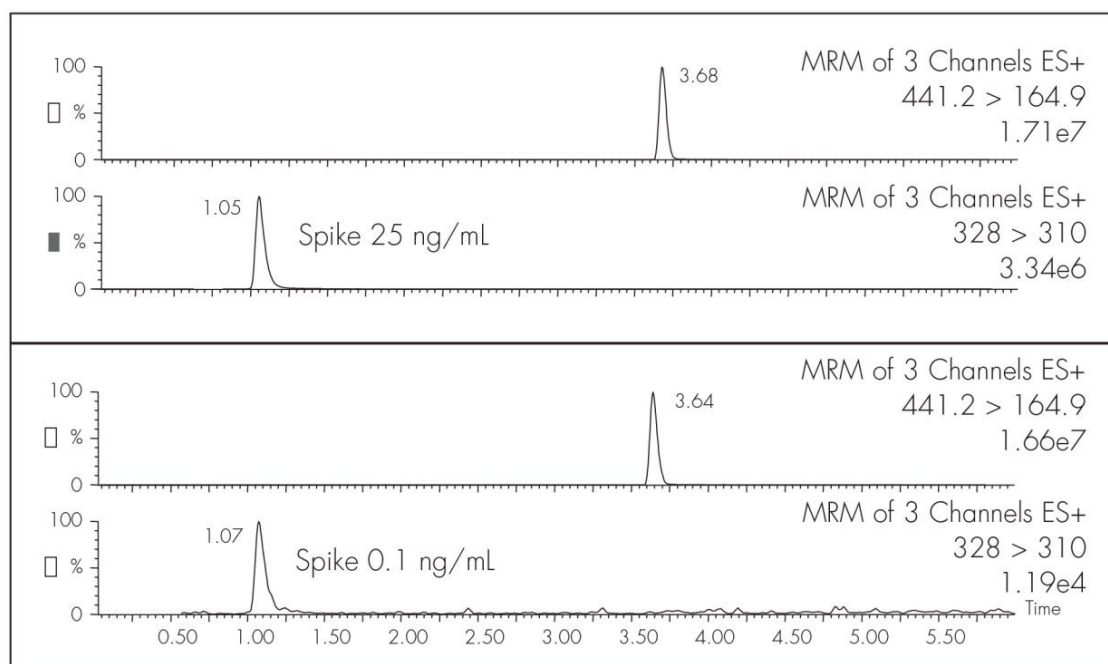

Featured Products

· [Alliance HPLC <https://www.waters.com/514248>](https://www.waters.com/514248)

WA31764.112, June 2003



© 2021 Waters Corporation. All Rights Reserved.