

## Propranolol in Rat Plasma by Mixed-Mode SPE and LC-MS/MS (Low Elution Volume)

---

Waters Corporation



This is an Application Brief and does not contain a detailed Experimental section.

---


### Abstract

This application brief demonstrates the analysis of Propranolol in Rat Plasma by LC-MS/MS.

---

## Introduction

The compound analyzed in this study is Propranolol.

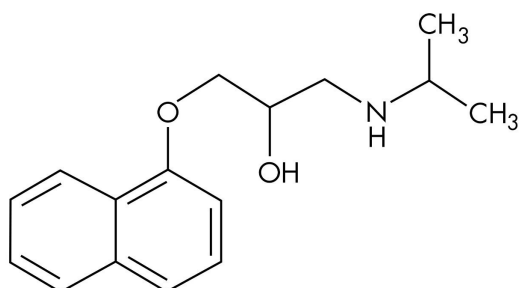

PROPRANOLOL

---

## Experimental

### HPLC Method

Column:	XTerra MS C <sub>18</sub> , 2.1 x 30 mm, 3.5 μm
Part number:	186000398
Mobile phase A:	Water + 0.5 % NH <sub>4</sub> OH
Mobile phase B:	ACN + 0.5 % NH <sub>4</sub> OH
Flow rate:	0.2 mL/min
Temperature:	Ambient
LC instrument:	Alliance 2795

### Gradient

Time (min)	Profile
------------	---------

	%A
--	----

0	5
---	---

1	95
---	----

## MS Conditions

MS instrument:	Micromass Quattro Triple Quadrupole
----------------	-------------------------------------

Ion source:	ESI+
-------------	------

Source temp.:	150 °C
---------------	--------

Gas cell:	2.0 e <sup>-3</sup> bar Argon
-----------	-------------------------------

Desolvation temp.:	350 °C
--------------------	--------

Drying gas flow:	500 L/hr
------------------	----------

Cone gas flow:	50 L/hr
----------------	---------


Cone voltage:	25 V
---------------	------

Collision energy:	20
-------------------	----

Capillary voltage:	3.5 Kv
--------------------	--------

MRM transition:	Metoclopramide (IS) <i>m/z</i> 299.8 → 226.7
	Propranolol <i>m/z</i> 259.9 → 154.9
	Amitriptyline <i>m/z</i> 278.1 → 232.9
	Nortriptyline <i>m/z</i> 263.9 → 190.8


# NEW! OASIS® $\mu$ ELUTION PLATE

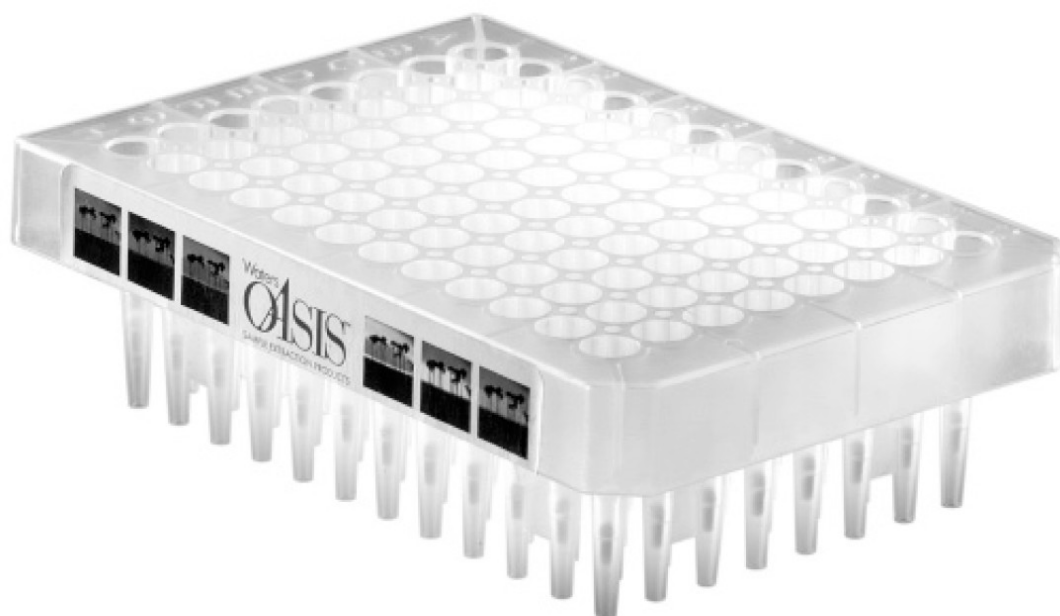

## OASIS® MCX GENERIC EXTRACTION PROTOCOL

Conditions for Oasis® MCX  $\mu$ Elution Plate



Part Number 186001830

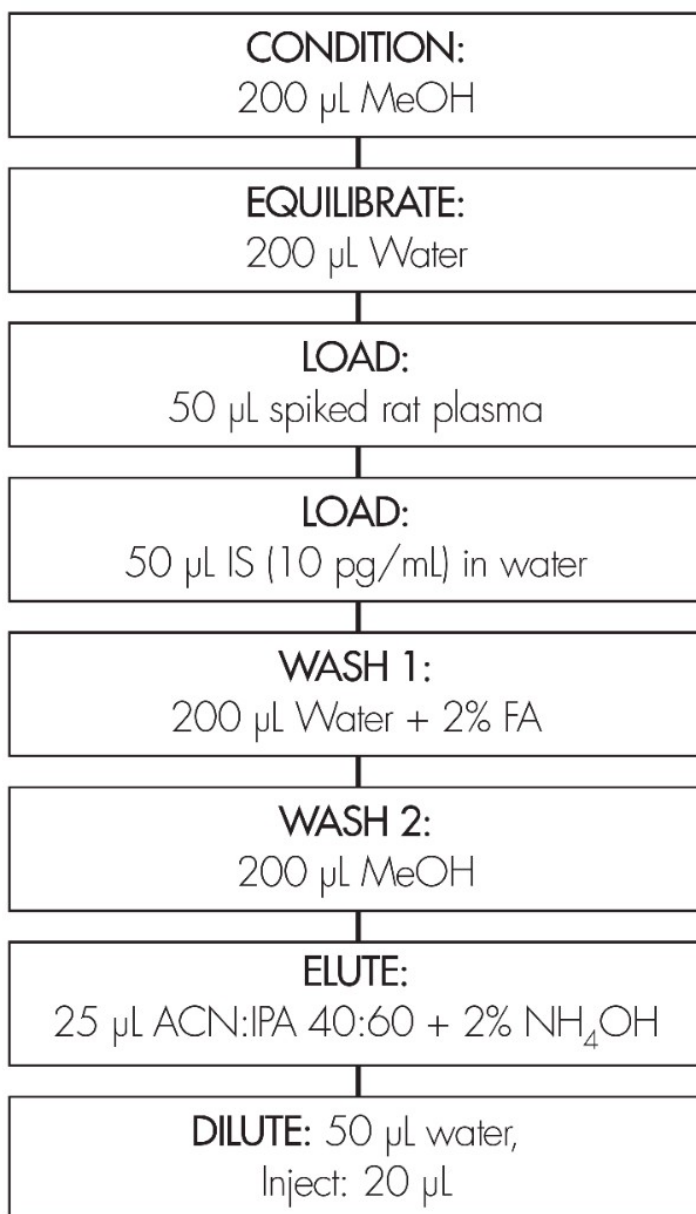

---

## Results and Discussion

# CALIBRATION CURVE

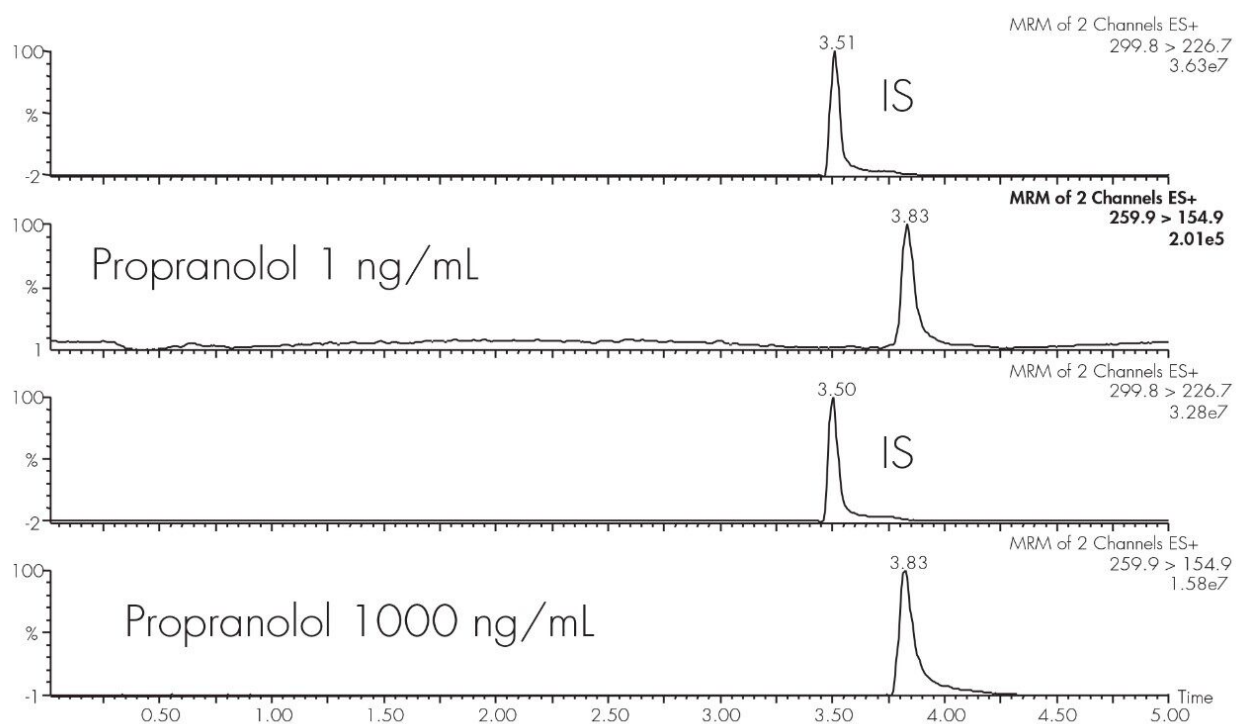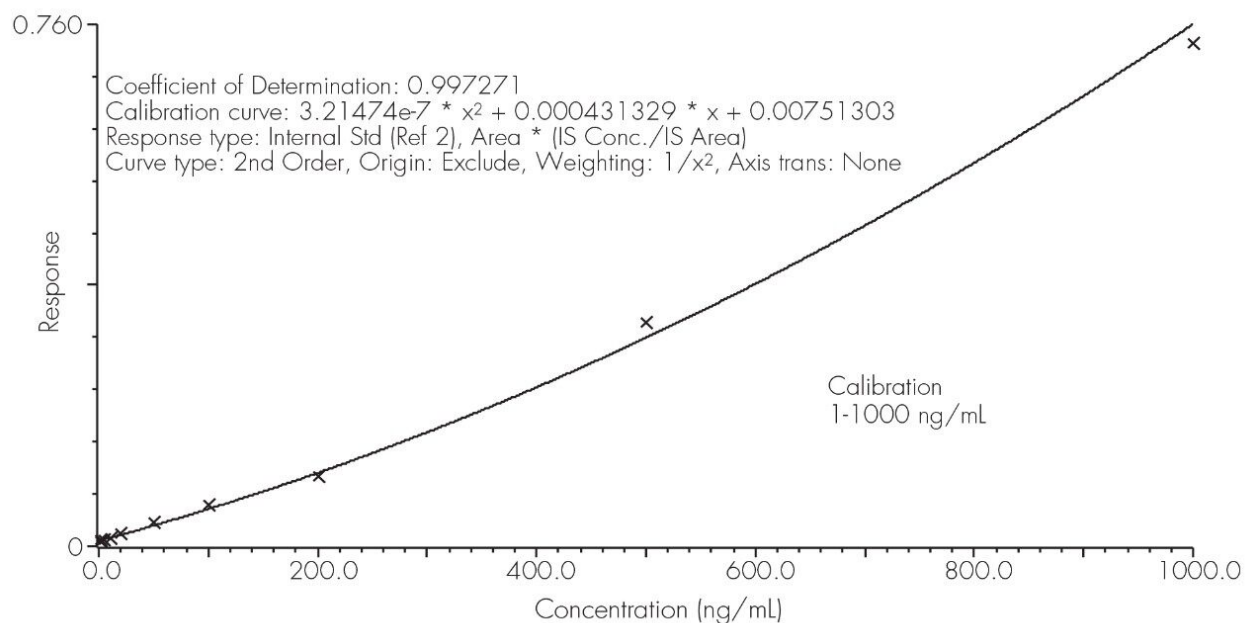

---

## Featured Products

WA31764.132, June 2003

© 2021 Waters Corporation. All Rights Reserved.