

## Application Note

# Barbiturates in Human Urine using Oasis Sample Preparation Products

---

Waters Corporation

*For forensic toxicology use only.*

This is an Application Brief and does not contain a detailed Experimental section.

---

## Abstract

This application brief demonstrates analysis of barbiturates in human urine.

---


## Introduction

The compounds used in this study are –


1. Phenobarbital
  2. Butabarbital
  3. Butalbital
  4. Amobarbital (I.S.)
-

5. Mephobarbital


6. Secobarbital

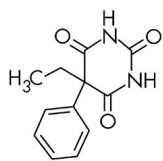


PHENOBARBITAL

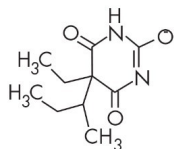


BUTABARBITAL

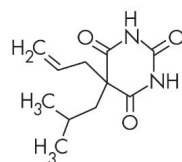


BUTALBITAL

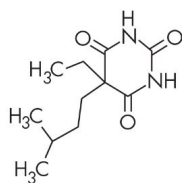

AMOBARBITAL (I.S.)

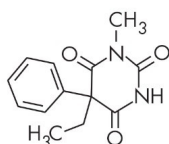

MEPHOBARBITAL

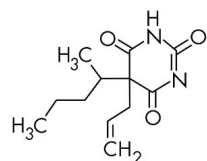

SECOBARBITAL

---

## Experimental

### HPLC Method

Column:

Symmetry Shield RP18, 2.1 x 150 mm, 5 µm

Guard column:

Symmetry Shield RP18, 3.9 x 20 mm, 5 µm

Part numbers:

Column - 186000111, Guard - 186000107


Mobile phase:

50 mM potassium phosphate, pH 7.0/acetonitrile  
71:29

Flow rate:	1 mL/min
Injection volume:	80 µL urine extract
Temperature:	30 °C
Detection:	UV @ 214 nm (0.350 AUFS)

## OASIS® HLB EXTRACTION METHOD


Oasis® HLB Extraction Plate, 30 mg/96-well  
Part Number WAT058951

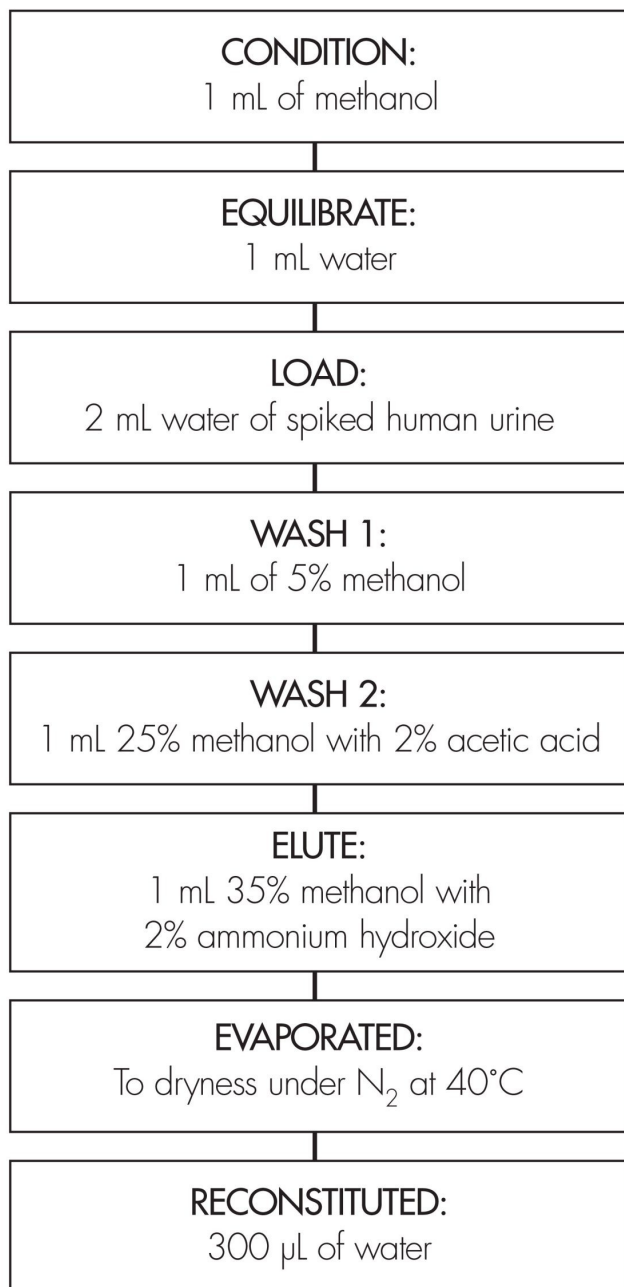

## Results and Discussion

Compound	%Recovery (n=8) 0.2 µg/mL	(%RSD) (n=8) 1.0 µg/mL
Phenobarbital	114.3 (1.7)	106.5 (0.5)
Butabarbital	95.7 (1.3)	105.5 (0.7)
Butalbital	109.5 (0.9)	104.2 (0.9)
Amobarbital		86.3 (1.7)
Mephobarbital	92.5 (3.6)	92.4 (1.7)
Secobarbital	101.5 (5.2)	94.8 (2.2)

### Chromatogram of A) Blank Urine, B) Spiked Urine

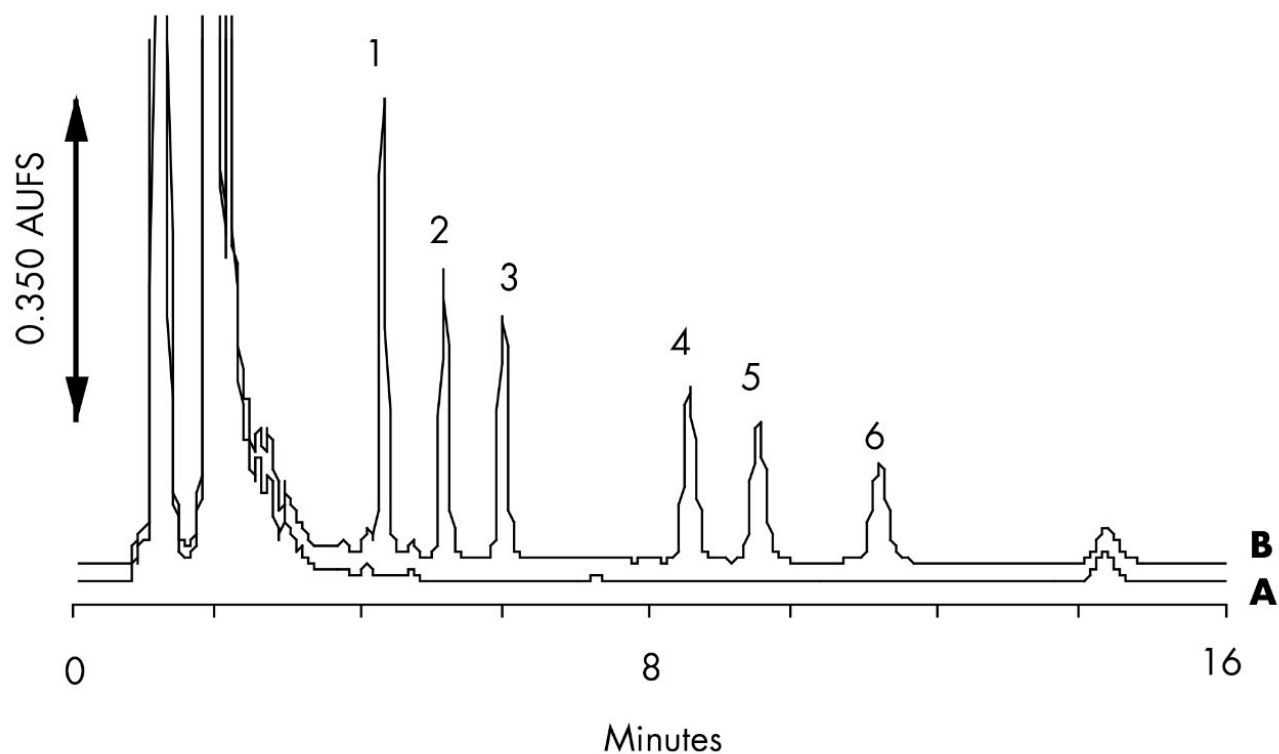

---

## Featured Products

WA31764.48, June 2003

© 2022 Waters Corporation. All Rights Reserved.