Waters[™]

ASTM D5501- 04 for Denatured Fuel Ethanol: Automating Calculations and Reports with Empower 2 Software

Larry Meeker, Alice J. Di Gioia

Waters Corporation



Abstract

This application note briefly describes the use of Empower 2 Software for denatured fuel ethanol analysis and reporting.

Benefits

Easy-to-use to analyze and report fuel ethanol according to ASTM D5501-04

Introduction

Ethanol is becoming an increasingly important alternative fuel on a worldwide scale due to social, environmental, political and economic factors. Production from fermentation alone in the U.S. has increased from 1.6 billion gallons in 2000 to 3.4 billion gallons in 2004.¹ To ensure use as a fuel product, the ethanol is denatured by the addition of alkanes before shipment. ASTM2 method D5501-04, "Determination of Ethanol Content of Denatured Fuel Ethanol by Gas Chromatography," details the analysis.

Producer labs and analysts are busy monitoring the lengthy ethanol fermentation process. Using Waters Empower 2 Software to streamline the final denatured ethanol analysis can save time. When an analyst can automate GC column conditioning and equilibration, data processing and calculations according to ASTM D5501-04 recommendations and print a report by filling in a small amount of information and pressing a Run button, that analyst is available to do other work in the lab or plant. Record keeping is easier with Empower as well. The sample identifiers that a lab uses, such as, sample name, batch, lot number, are trackable, traceable and searchable.

This application note briefly describes the use of Empower 2 Software for denatured fuel ethanol analysis and reporting. The analytical system was comprised of an Agilent 6850 GC with an FID detector and a client server running Empower 2.

Experimental

Standards were from a commercial source for the ASTM method.

Conditions, Software and System

System:	Agilent 6850 GC with FID
Software:	Empower 2
Column:	DB-1 100 m x 0.25 µ film
Conditions:	ASTM D5501-04

Results and Discussion

Setting Parameters

With Empower 2 Software anyone in the lab can easily automate the calculation and reporting of all the required values outlined in Section 12 of ASTM D5501- 04. The GC must be set up and run with parameters to comply with the ASTM method. These parameters are entered on individual tabs in the Empower instrument method (Figures 1–6).

Pa D5501 ii	n Larry GCW-5	501 2004 as	System/Admi	nistrator - Ins	trument Met	hod Editor	
<u>Eile E</u> dit <u>V</u> i	iew <u>H</u> elp						
		5 66850					
Instrumen	t Oven Inlet	Column Det	ector Channel	Injector Events	Valves Auxil	iary	-
IJ ∪ 	Oven Enable Oven Setpoints Initial Tim Initial Temperatur	e 15.00 e 60 Rate (*C/min)	C tyo Enable Min E *C Final Temp (*C)	e quilibration Time Maximum Temp. Hold Time (min)	Temp F 0.00 350 Total Time (min)	Ramp Enable Min °C	?
	1 2 3 4 5	30.00	230	15.00	35.67		
	Cryogenic Options	Cooling Enable It Detection Ena etection Enable	ble	Ambient Temp Timeout Value	20	°C Min	
• Ready							

Figure 1. Oven Parameters.

🐴 D5501 in Larry GCW-5501 2004 as System/Administrator - Instrument Method Editor	
Eile Edit Yiew Help	
	•
Instrument Oven Inlet Column Detector Channel Injector Events Valves Auxiliary	
- Settings	?
Inlet Type Split/Splitless Mode Split	
Temperature	
I I Initial Temp	
Gas Saver	
Enable On Time 0.00 min Flow 15.0 ml/min	
Split/Splitless Purge	
Split Ratio 200.0 to 1 Flow 0.0 ml/min	
Split Flow 200.0 ml/min 0.00 min	
Pulse Time 0.00 min	
Pulse Pressure 00 psi	
	_
٠	
Ready	11.

Figure 2. Inlet Parameters.

D5501 in Larry GC\D-5501 2004	1 as System/Ad	ministrator -	Instrument Me	thod Editor	
<u>Eile E</u> dit <u>V</u> iew <u>H</u> elp					
Instrument Oven Inlet Column	Detector Channe	el Injector Eve	ents Valves Aux	iliary	<u>*</u>
Column Information					?
Length (m) 150.0	De	scription DB-1	100m×.251uflm		
Diameter (u) 250	_ Serial	Number			
Film Thickness (u) 1 00	-		-		
1.00					
Inlet Connection Inlet	Detector Co	nnection Detec	tor 🗾		
Vacuum Correct Pro	essure Correct	Pressure	Correction Value	0.00	
Inlet Settings					
Mode Constant Flow	-				
Initial Time					
1.00 min	Rate	Final	Hold Time	Total Time	-
Initial Flow	Trate	T N KM	(min)	(min)	_
1.0 ml/min 2	-				
Initial Pressure 3					
1.0					
<u> </u>					
Ready					1.

Figure 3. Column Parameters.

a D5501 in Larry GCID-5501 2004 as System/Administrator - Instrument Method Editor	
Eile Edit View Help	
Instrument Oven Inlet Column Detector Channel Injector Events Valves Auxiliary	<u> </u>
Settings	?
Source Detector	
Sampling Rate 100.0	
Durinting Inc.	
Description FID	
Analog Output Settings	
Enable Fast Peaks Range 0	
Zero Enable Offset 0.0 Attenuation Attenuation Value 0	
l • I Ready	_

Figure 4. Detector Parameters.

🐴 D5501 in Larry GCID-5501 2004 as System/Administrator - Instrument Method Editor 💦 📕	
<u>File Edit View H</u> elp	
Instrument Oven Inlet Column Detector Channel Injector Events Valves Auxiliary	
General Detector Type FD General Detector Type FD General Temperature Fenable Setpoint 250 *C Makeup Gas Flow Mode Constant Plus Column Gas Type N2 Flo	?
Image: Flame Enable Fuel Flow 40.0 ml/min Lit Offset 2.0 pA Image: Flame Enable Oxidizer Flow 400.0 ml/min	
TCD Filament Enable Negative Polarity Off Reference Flow 1.0 ml/min	_
Ready	

Figure 5. Channel (Sampling Rate) Parameters.

🚹 D5501 in Larry GCW-5501 2004 as System/Administrator - Instrument Method Editor 👘 🔲 🔲
Elle Edit View Help
Instrument Oven Inlet Column Detector Channel Injector Events Valves Auxiliary
Solvent A Pre-Washes 3 Skim Enable
Solvent B Pre-Washes 3 Skim Depth 100 (x 0.1 mm)
Sample Washes 3 TSlow Plunger Enable
Injector Pumps 6 Pre Dwell Time 0 (x 0.01 min)
Viscosity Delay 0 sec Post Dwell Time 0 (x 0.01 min)
Solvent A Post-Washes 6
Solvent B Post-Washes 6
4
Ready

Figure 6. Injector Parameters.

Running Samples

To run samples, the next area of the software to use is the RUN SAMPLES page (Figure 7). When a sample template is used, the lab analyst only enters a minimum of information reducing the data entry burden. Data fromtwo other required tests can be entered, stored, searched and plotted. Ds is the sample density determined using ASTM method D 1298 or D 4052 and % water is calculated using ASTM D1364 or E 1064. These data are needed to calculate the final % Mass of alcohols. If the sample density and % water are unavailable when the GC analysis is run, the values can be entered later and the final result calculated.

C 6850 in Larry GCD-5501 2004 as System/Administrator - Run Samples										
Elle Edit View Inject Actions Customize Help										
🔥 🗞 🖶 🥥 🔌 🌭 🧼 🗰 🔜 🗶 🖹 📰 🕺 Run and Process 💌 Reinject on Fault						•				
					5	Sample	Set Method: test			
ø	Vial	SampleName	Ds	Percent_Water	Inj Vol (uL)	#of Injs	Function	Run Time (Minutes)	s	
1							Condition Column	30.00		
2							Equilibrate	30.00		
3	1	Ethanol STD	0.7940	0.0	1.0	1	Inject Standards	30.00		
4	2	sample 1	0.7883	0.5	1.0	3	Inject Samples	30.00		
5	3	check STD	0.7940	0.0	1.0	1	Inject Controls	30.00		
6							Clear Calibration 💌			
7							Calibrate			
8							Quantitate			
9							Report			
		Cingle) Sa	mplac	Comple Cate	/ Du	nning				
1	17.0	Sindle V24	inples /	V cample sets	ARU	miniq		-		
	Instrument Method:									
	D5501									
	Edit Monitor Setup									
For Help, press F1 System Idle 🕅 🕅 🖋 🐙 🖉						* # C/				

Figure 7. Running Samples.

Automatic Monitoring and Processing

Critical to success in daily operation is ensuring that unattended systems generate quality data. With Empower 2 you can do more than just set up initial conditions and wait for temperatures to meet set points. The software has features that allow easy automation and monitoring of system setup, column conditioning, system equilibration and calibration, running samples, performing required calculations, and printing a custom report. This interactive process ensures the analysis begins, runs properly and generates data. Control values can also be set using the System Suitability option. This feature analyzes the data from eachinjection relative to defined parameters, such as retention time or response. If the value is outside tolerance, the user can set the system to take corrective action, for example, re-inject of the sample or stopping the run completely.

Automatic Calculations and Reporting

After the analysis has been run, Empower 2 can be set up to automatically process the data, performcalculations and report results. There are three key benefits to these features. First is the ability to smooth the GC FID data with software as opposed to filtering the data in the hardware.

In the processing method the user can define both the smoothing type and by how much the raw data should be smoothed (Figure 8). This is important with high speed data typically collected today. The data shown in Figure 11 were collected at 100 Hz. Second, performing calculations to meet the needs of Section

12 of ASTMD5501 can be done automatically with Empower 2 instead of using a calculator or spreadsheet.

😸 D5501sample B in Larry GCW-5501 2004 as System/Administr 🔳 🗖 🔀
Eile Edit View Plot Process Navigate Options Window Help
46 26 28 28 36 28 28 28 28 28 28 28 28 28 28 28 28 28
Integration Smoothing/Offset Components Default Amounts Named Groups Tim
Concentration
Time Offset (min) 0.000
Smoothing Type Smoothing Level
Mean 💌 31 💌
None Savitsky-Golay Mean
For Help, press F1 Meth Set: Untitled

Figure 8. Smoothing Parameters.

A unique feature of Waters Empower 2 Software is Custom Fields, a tool that lets users define equations for their needs in standard math syntax not a programming code (Figure 9). The interface is as easy to use as writing an equation on a piece of paper. These custom fields meet the needs of Section 12.

Pro	Project Properties - 'Larry GC\D-5501 2004'							
Ge	General Custom Fields Access Integrity Processing							
	3 9	Name	Formula	CField Id	All or 1			
	1	Adj_area	Area*CConst1	103	Г			
	2	Dilution		97	Г			
	3	Ds		108	Г			
	4	Mi	RMi*(100-Percent_VVater)/100	102	Г			
	5	Percent_Water		100	Г			
	6	RMi	Adj_area*100/total_adj_area	101	Г			
	7	SampleName		99	Г			
	8	SampleWeight		98	Г			
	9	Vi	Mi*Ds/CConst3	107	Г			
1	10	total_adj_area	SUM(Adj_area)	104	Г			
	<				>			
	New Edit Delete Lock/Unlock Save to Project							
	OK Cancel Help							

Figure 9. Equations for ASTM 5501 Calculations.

The equation for the custom field, %mass or Mi, was created using simple math. Empower 2 is easier to use than writing an equation on a piece of paper (Figure 10).

Edit Custom Field	- Mi
Field Type C Sample	Name: Mi
C <u>R</u> esult	Width : 12 Translation Definition:
• Peak • Sample Set	Precision : 3
C Component	Min.: -399999993.399
	Mag. : 100000000.000
Data Type	Default Value : 🔲 User Entry Required 🔲 Custom Field Locked
C Integer (0)	
Real (0.0)	Calculation Criteria
C Lext	Search Order: Use As:
C Date	Result Set First All or Nothing
C B <u>o</u> olean	Sample Type: All
C Enu <u>m</u>	Peak Type: All 💽 Missing Peak
Data Source	Calculated Field Formula
C Keyboard	RMi*(100-Percent_Water)/100
C External	
• <u>C</u> alculated	E dit <u>F</u> ormula
	OK Cancel <u>H</u> elp

Figure 10. Creating equations using simple math.

The third key benefit of Empower 2 software is the reporting capability. After setting up templates and fields one time, the system can automatically print a report. The report is fully customizable. Figure 11 is an example of such a report.



Figure 11. Empower 2 report.

Besides sample information, Empower 2 software can be used to generate reports for tracking and trending data. Figures 12 and 13 are summary plots of 14 ethanol batches showing the % Mass (Mi) for methanol, ethanol and the heptanes. The Limit bars in Figure 13 were set as part of the system suitability.



Figure 12. Methanol.



Figure 13. Ethanol and Heptane.

Conclusion

Empower 2 Software is easy-to-use to analyze and report fuel ethanol according to ASTM D5501-04. The software allows use of a template that includes instrument parameters such as conditioning and equilibration, data processing instructions and report instructions. The equations needed to calculate required values, such as % mass, can be built-in one time with the Custom Field feature and re-used. All the analyst must do is load the sample and a template, enter minimum information and hit run. After the run is finished, the analyst can be confident that quality data are printed. Using interactive system suitability, runs can be re-injected or stopped based on user pre-defined limits. With the System Suitability option, more advanced trending and tracking data features are available.

References

- 1. BioFuels Journal, posted 3/7/2005, Annual and Monthly U.S. Ethanol Production.
- 2. www.astm.org D5501-04 Standard Test Method for Determination of Ethanol Content of Denatured Fuel Ethanol by Gas Chromatography.

Featured Products

Empower Chromatography Data System https://www.waters.com/513188

720001840, July 2006

© 2021 Waters Corporation. All Rights Reserved.