

Synthetic Reaction Monitoring Using UPLC-MS

Marian Twohig, Darcy Shave, Paul Lefebvre, Robert S. Plumb

Waters Corporation

Abstract

This application note demonstrates the advantages of speed and ease of use that self-service UPLC with photodiode array (PDA)/evaporative light scattering (ELS)-MS detection brings to reaction monitoring studies.

Introduction

Once a chemical hit is found through a library screening process and is verified, optimization of the compounds' desired properties takes place. This step involves an iterative process of synthesis and reactivity measurement of the new compounds to further develop drug candidates into the lead phase.

Because these reactions may take a long time, chemists need to know as soon as possible if their syntheses are proceeding as desired. This means utilizing measurement capabilities that require minimal sample preparation and provide a fast response giving low detection limits. Another advantageous property might be the ability to measure multiple parameters simultaneously.¹

High throughput approaches can provide important time savings in the optimization of process parameters. Open access LC-MS is replacing TLC as a reaction monitoring tool.² Sample preparation of reaction mixtures can be as minimal as filtering and dilution before injecting into the LC-MS system. This allows fast turnaround of results to allow the chemist to advance to the next step.

The purpose of this application note is to demonstrate the advantages of speed and ease of use that self-service UPLC with photodiode array (PDA)/evaporative light scattering (ELS)-MS detection brings to reaction monitoring studies.



Figure 1. The ACQUITY SQD for synthetic reaction monitoring.

Experimental

Chromatographic separations were carried out using an ACQUITY UPLC System coupled to an ACQUITY SQ Mass Detector. PDA and ELS signals were collected simultaneously. Samples were analyzed using gradients less than 1 minute. For chromatographic flexibility, a column selection module was added.

LC conditions

LC system:	Waters ACQUITY UPLC System
Column:	ACQUITY UPLC BEH C ₈ Column 2.1 x 30 mm, 1.7 μ m
Column temp.:	45 °C

Flow rate:	800 μ L/min
Mobile phase A:	0.1% Formic acid in water
Mobile phase B:	0.1% Formic acid in acetonitrile
Gradient:	5 to 95% B/0.7 min

MS conditions

MS system:	Waters SQ Detector
Ionization mode:	ESI positive/ESI negative
Capillary voltage:	3.0 KV
Cone voltage:	20 V
Source temp.:	150 °C
Desolvation temp.:	450 °C
Desolvation gas:	L/Hr
Cone gas:	50 L/Hr
Acquisition range:	100 to 1300 <i>m/z</i>
Scan speed:	10,000 amu/sec

Note: A low volume micro-tee was used to split the flow to the ELS and SQ.

ELS conditions

Gain:	500
N ₂ gas pressure:	50 psi
Drift tube temp.:	50 psi
Sampling rate:	20 points/sec

PDA conditions

Range:	210 to 400 nm
Sampling rate:	20 points/sec

Results and Discussion

During the compound optimization stage of a discovery cycle, medicinal chemists are not only interested in determining the key structural features responsible for activity and selectivity, but also what structural changes need to be made to improve these characteristics. Because the reactions necessary to bring about these changes may take many steps, chemists need to be sure they are progressing as expected during the course of the reaction synthesis.

To illustrate the functionality of such a system, the synthesis of atenolol (Figure 2) was used as a reaction model. The increase in the formation of atenolol was monitored, as was the decrease in the intermediate 4-hydroxyacetamide³ (Figure 3). A reaction by-product 4-hydroxyphenylacetic acid was also observed.

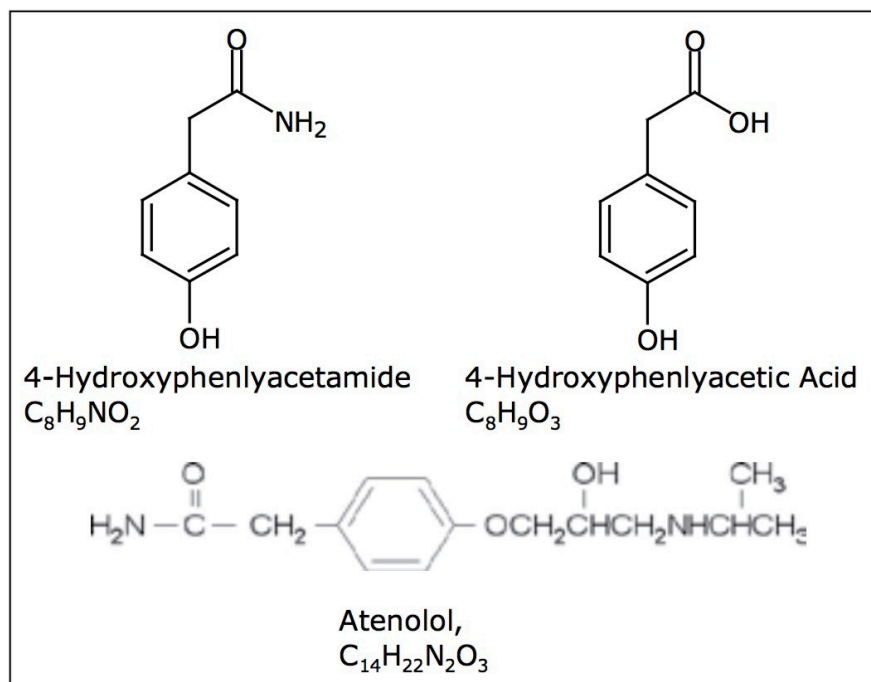


Figure 2. Structures of atenolol and 4-hydroxyacetamide.

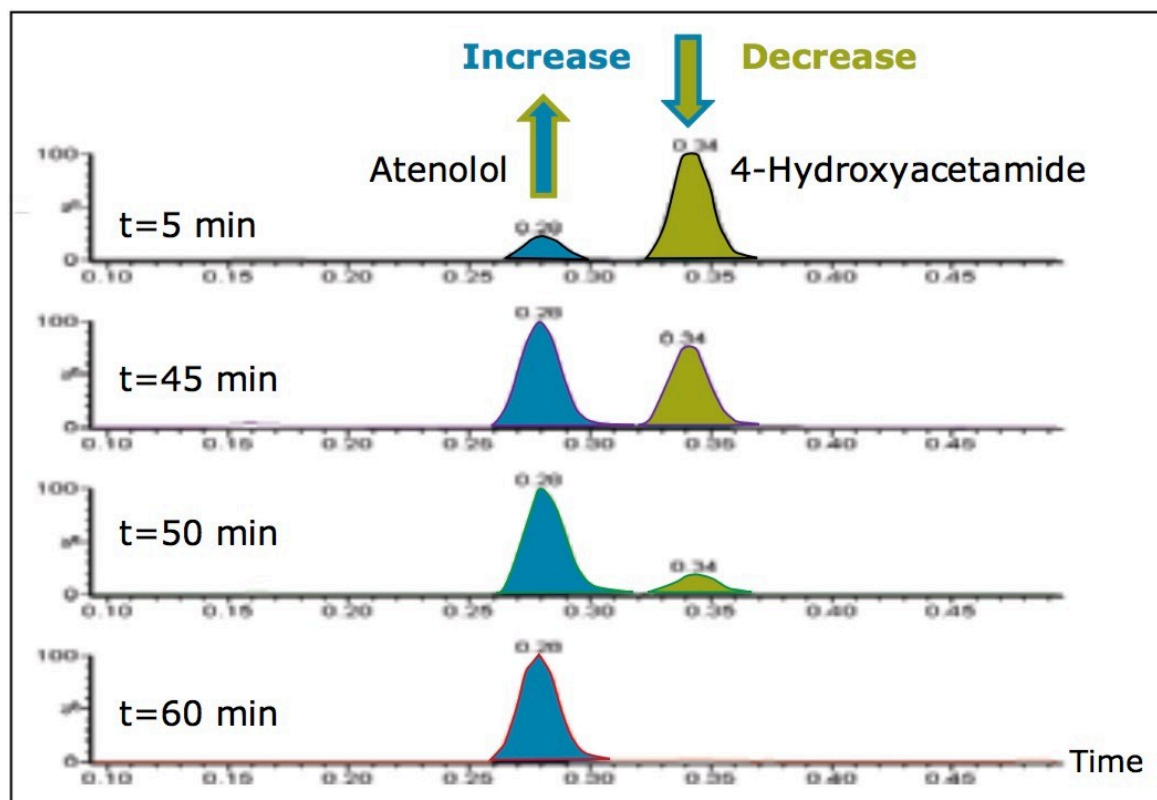


Figure 3. UPLC-MS chromatograms. The reaction mixture was sampled at various time points.

The ACQUITY SQD is capable of scan speeds of up to 10,000 amu/sec. Consequently, it is possible to employ a

large number of scan functions in a single run while still maintaining adequate peak characterization. The fast scan speed is essential for this functionality, as peak widths of 1 second or less are common with the use of UPLC. Scanning multiple functions allows confirmation of compound synthesis to be obtained on reaction components whether they ionize in positive ion mode or negative ion mode, ESI or APCI. The total cycle time of the method was 1 minute 20 seconds, facilitating increased sample throughput.

A single run can also provide UV spectral information and an estimation of compound purity at low wavelengths. ELS detection is based upon the degree to which solute particles scatter light. It has been known to give rise to similar responses for related compounds.⁴ The signal can give a tentative estimation on the relative quantities of the components present (Figure 4). It is also an alternative detector to UV, which depends on the presence of a chromophore. As can be seen from Figure 4, atenolol ionizes in ESI positive ion mode (retention time 0.28 min). The reaction intermediate 4-hydroxyphenylacetamide ionizes in both positive and negative ion mode (Rt 0.34 min) and 4-hydroxyphenylacetic acid (Rt 0.39 min) only ionizes in negative ion mode.

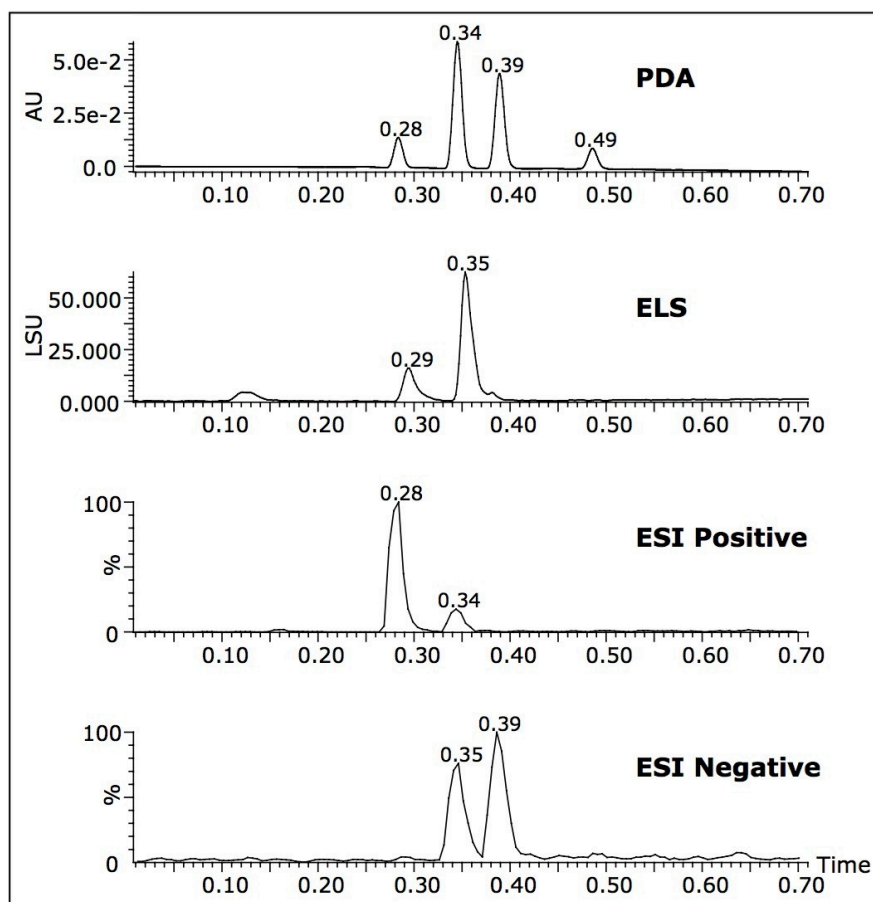


Figure 4. UPLC/PDA/ELS-MS chromatograms.

The OpenLynx Open Access Application Manager, part of MassLynx Software, allows chemists to walk up to a

terminal and log in samples while entering the minimum information required to run the samples.

The OpenLynx OA Login screen shown in Figure 5 allows the administrator to set up the system such that the user only needs to input the information requested, and then upon completion, select the Login Samples button. This will either tell the user the designated autosampler position, or confirm the position that the user has chosen, and ask for confirmation of position before it will run the sample. In addition to a simplified sample submission process, the OpenLynx Application Manager can then process data automatically and produce a summary report that can be emailed or printed as desired.

OA Login

Login Information

Your Name: E.L. Job ID: 9-26_18

Method: C1_waters/EndcapH_F18_P01_Meg Description: LC/MS method with broad range (5-95%) gradient over 1.7 minute at a flow rate of 0.3mL/min. Acids Mobile Phase B: 1.75M TFA. Performance: none

☐ E-Mail Results to Address:

Sample Holder: 04AUC148Vial2ndHolder (Single Shot - 40 wells/walk)

Holder Details: 8 Rows x 8 Columns, 32.0mm depth, 12.0mm diameter

Sample Details

Number of Samples: 1

SampleID: RM

Mass: 286

Wavelength: 254

Press button to login samples. There are no jobs waiting

Figure 5. The OpenLynx single page login.

The information contained in the summary report is viewed via the OpenLynx browser shown in Figure 6. It clearly defines what components are found or not found. Chromatograms and spectra are generated based on

the processing parameters set up by the administrator in the OpenLynx method.

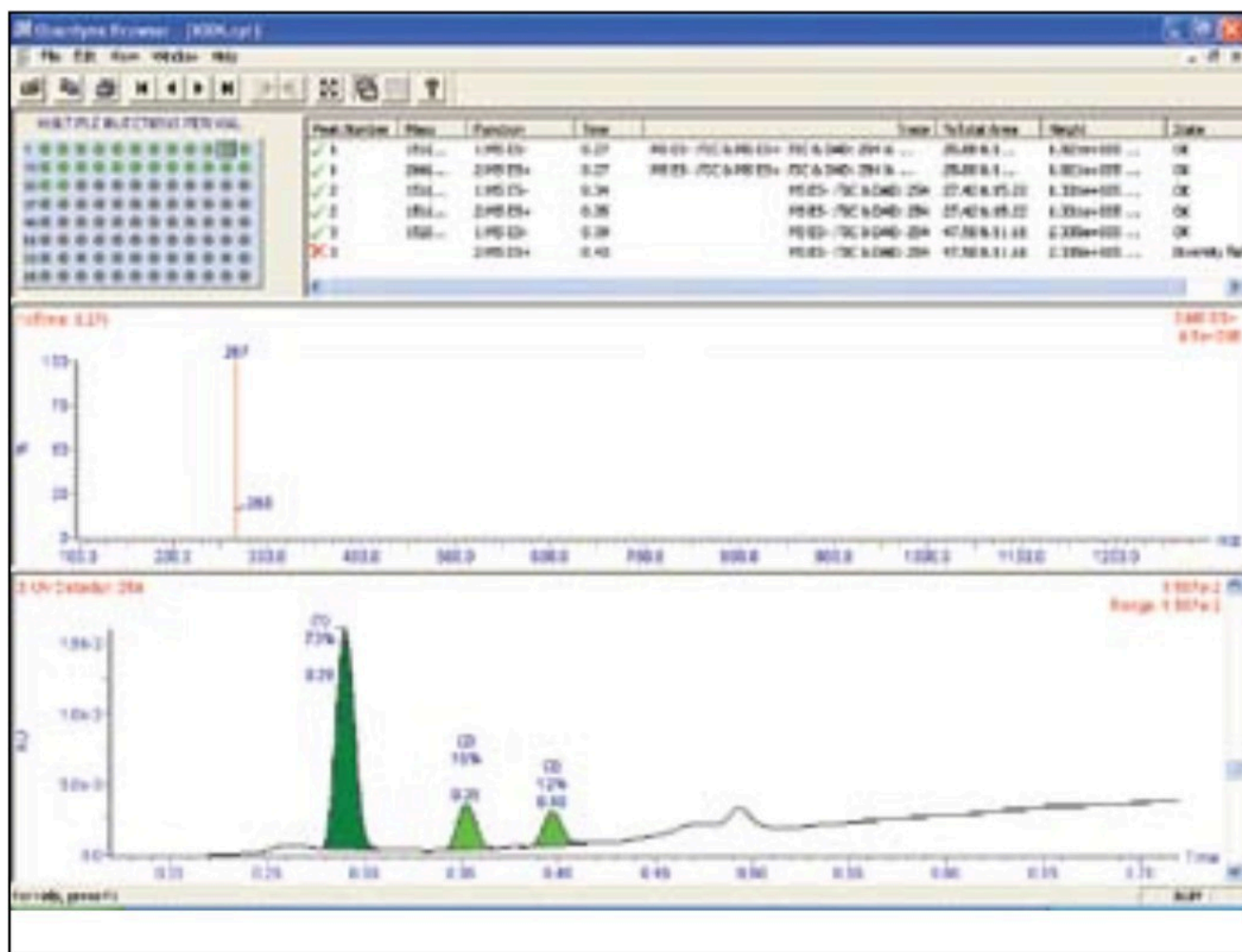


Figure 6. The OpenLynx Application Manager browser.

Conclusion

During the compound optimization stage of a discovery cycle, medicinal chemists are not only interested in determining the key structural features responsible for activity and selectivity, but also what structural changes need to be made to improve these characteristics. Because the reactions necessary to bring about these changes may take a long time, chemists need to be sure they are progressing as expected.

By using a walk-up UPLC-MS system, chemists were able to quickly and easily monitor their reactions, noting the relative amounts of starting materials and products. They were also able to note the formation of any side products and make necessary alterations to their reaction protocol to minimize these.

The described system and software combination can autonomously evaluate large numbers of samples, with a cycle time of 1 minute 20 seconds. Data can then be automatically processed and a summary report can be generated. The scan speed capabilities of the ACQUITY SQD make it possible to better characterize narrow chromatographic peaks. This has become a necessity since the advent of sub-2 μm particle technology, where chromatographic peaks can be 1 second wide or less. The fast scan speed allows the chemist to extract as much data as possible per injection by switching between APCI and ESI as well as positive and negative ion modes.

Open Access gives the chemist a walk-up system that is flexible for analytical data acquisition. It runs as a complete system, from sample introduction to end results.

The use of the fast-scanning MS along with the throughput of UPLC technology allows the chemist to obtain high quality and comprehensive data about their compounds in the shortest possible time. This combined with intelligent open access software allows informed decisions to be made faster, thus supporting the needs of the modern drug discovery process.

References

1. Analysis and Purification Methods in Combinatorial Chemistry, Wiley-Interscience. (5): 87-123.
2. LC/MS Applications in Drug Development, Wiley-Interscience. 96-106.
3. A Synthesis of Atenolol using a Nitrile Hydration Catalyst. Organic Process Research and Development. 1998; 2: 274-276.
4. Kibbey, C.E. Mol. Diversity. 1995; 1: 247-258.

Featured Products

- [ACQUITY UPLC System <https://www.waters.com/514207>](https://www.waters.com/514207)
- [SQ Detector 2 <https://www.waters.com/134631584>](https://www.waters.com/134631584)
- [OpenLynx Open Access <https://www.waters.com/10008851>](https://www.waters.com/10008851)

720002258, June 2007



©2019 Waters Corporation. All Rights Reserved.