


Analysis of Cephalosporin Antibiotics Using XBridge C₁₈

日本ウォーターズ株式会社

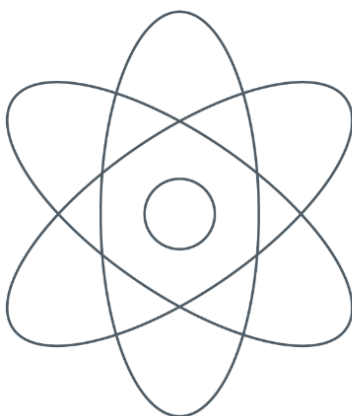


This is an Application Brief and does not contain a detailed Experimental section.


Abstract

This application brief highlights the analysis of cephalosporin antibiotics using XBridge C₁₈ Columns.


Introduction

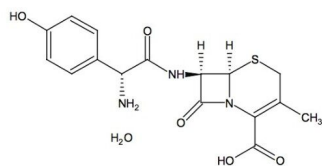


1. Cefadroxil

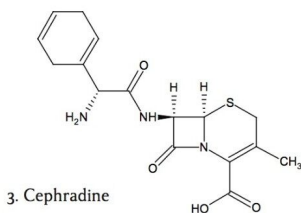

3. Cephadrine

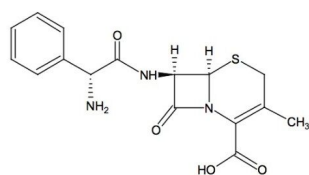

2. Cephalexin

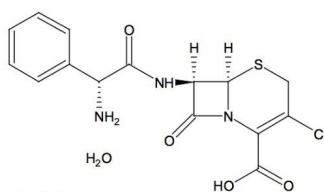

4. Cefaclor

Compounds used in this study.

Experimental


Test Conditions

Column:	XBridge C ₁₈ , 5 µm, 4.6 x 150 mm p/n: 186003116
Isocratic Mobile Phase:	0.5% acetic acid/acetonitrile (85:15)
Flow Rate:	1.0 mL/min
Injection Volume:	10 µL
Sample Concentration:	20 µg/mL in water
Temperature:	25 °C
Detection:	UV @ 254 nm

Instrument:

Alliance 2695 with 2996 PDA

Results and Discussion

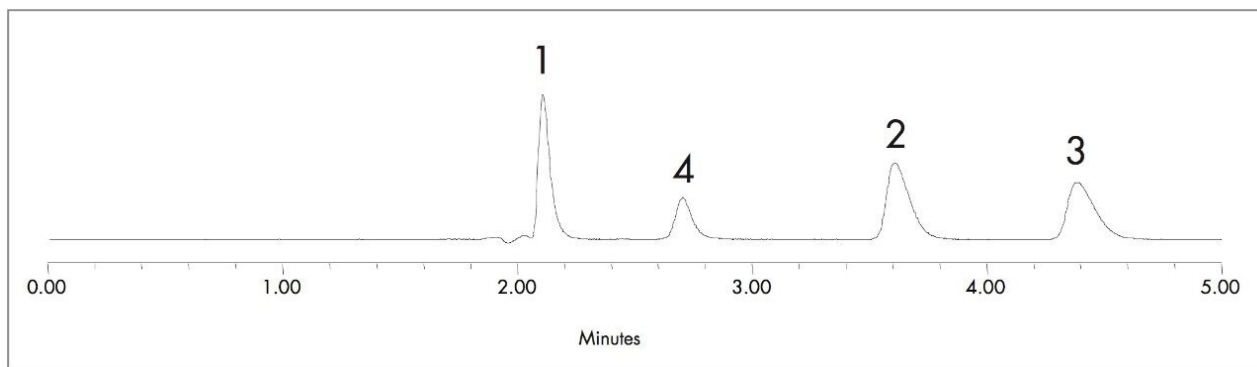

1. Cefadroxil 2. Cephadrine 3. Cephalexin 4. Cefaclor.

Featured Products

Alliance HPLC <<https://www.waters.com/514248>>

WA60165, April 2007