

An Automated Application Template for Content Uniformity Studies in SDMS Vision Publisher

Wolfgang Lemmerz, Chris L. Stumpf, Thomas Schmidt, Phil Kilby

Waters Corporation



This is an Application Brief and does not contain a detailed Experimental section.

Abstract

SDMS Vision Publisher serves as a reporting and authoring module incorporated as standard for NuGenesis SDMS and provides templates for Content Uniformity testing; whereas, NuGenesis SDMS acts as a data repository for analytical raw data as well as printed reports. A brief example will illustrate a Content Uniformity template for a tablet HPLC based assay. Other forms of Content Uniformity testing are also possible.

Introduction

Releasing a pharmaceutical drug product to market may require a Content Uniformity assessment to ensure that the downstream consumer receives the recommended dosage of active pharmaceutical ingredient (API). Content Uniformity serves as one of several tests (other tests include appearance, average mass, dissolution, etc.) that are performed during manufacturing and that adhere to GMP guidelines. As defined by the USP, content uniformity tests may be performed on coated tablets, transdermal systems, soft gel capsules, packaged inhalations, single packaged solids, and suppositories¹⁻² A typical Content Uniformity workflow is described in Figure 1 and consists of six primary steps. Each of these steps requires accurate documentation due to the adherence to GMP guidelines. Fortunately, the Content Uniformity workflow is straightforward and can make ready use of templates.

Content Uniformity studies employ a straightforward workflow that may be readily captured in a standardized template by using electronic report and authoring tools (see Figure 1 for workflow example). A standardized template both guides the analyst through the workflow and ensures reliable data capture without skipping necessary steps.³ Further benefits to recording Content Uniformity studies electronically include: searchable reports, electronic interface to analytical data systems (e.g., CDS, LIMS, balances), and electronic records that are 21 CFR Part 11 Compliant Ready.³

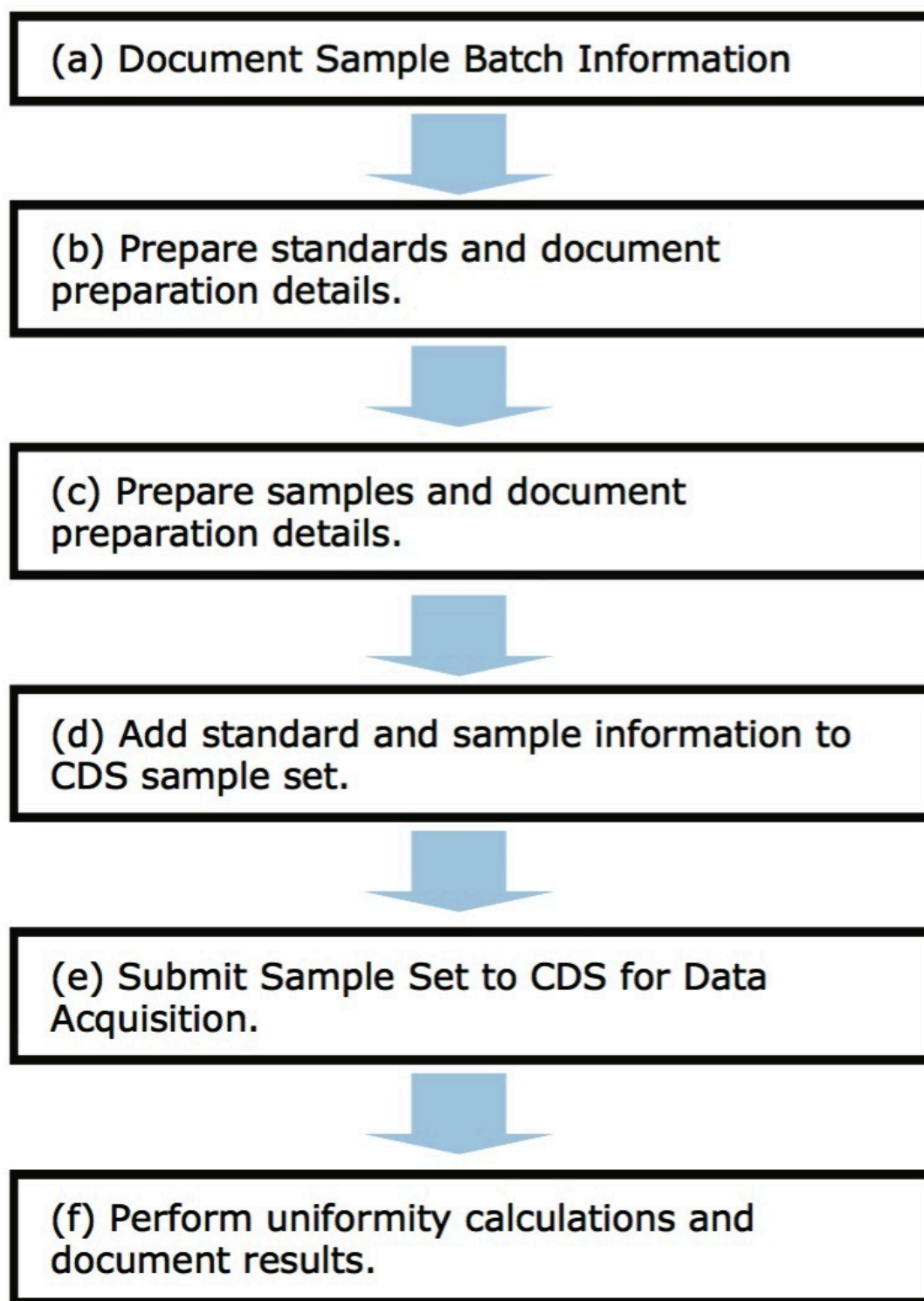


Figure 1. Typical

Content Uniformity workflow when using HPLC.

SDMS Vision Publisher serves as a reporting and authoring module incorporated as standard for NuGenesis

SDMS and provides templates for Content Uniformity testing; whereas, NuGenesis SDMS acts as a data

repository for analytical raw data as well as printed reports. A brief example will illustrate a Content Uniformity

template for a tablet HPLC based assay. Other forms of Content Uniformity testing are also possible.

Results and Discussion

Template Description

A typical tablet based content uniformity HPLC assay document requires four primary sections that include the reference standard preparation, the test sample preparation, the chromatographic system setup, along with the results and conclusions. The template highlighted in Figure 2 captures the aforementioned information as well as provides a convenient mechanism to automatically populate the Empower 2 sample set with reference standard and sample batch information, as well as automatic submission of the completed sample set to an Empower 2 System for acquisition. Subsequently, the results are then pulled back into the document from Empower 2 and the results are summarized.

Waters Vision Publisher - [080130_HPLC_0000(2)]

File Edit Document Tools Workspace Services Administration Options Window Help

Message Board Documents New Document Templates Workspaces New Workspace External Systems SDMS Project Browser

Append Delete Replace Properties History Sign Options Reload Save New Transfer Print

Filter: [Icons]

Title: Content Uniformity Template Example Document ID: 080130_HPLC_0000

Nr	Title	Ver
1	Batch Information	
2	Reference Standard	
3	Sample Preparation	
4	CDS Sample Set	
5	CDS Results	
6	Content Uniformity Plot	
7	Conclusions	
8	Status Report	

Batch Information

Method	Acetaminophen, Aspirin, and Caffeine Tablet- USP Monogram
Project (Drug)	
Study	
Analytical Run Batch	
Comment	

Description	Work sheet #	Lot #	Exp.Date
See Standard solutions preparation sheet:			

API's to monitor:

	1
--	---

Standard	Conc.	#	Component	Weight

Tablet masses:

Item Number	Balance Reading
Table 1:	
Table 2:	
Table 3:	
Table 4:	

080130_HPLC_0000(2) Jim Waters on wli-eu-demo Wednesday, J

Figure 2. An example Content Uniformity template. This template captures the reference standard preparation, sample preparation, Empower sample set, and results in a single electronic document.

Automated Record Keeping

An electronic document provides a convenient means to find data and records by allowing online searching. It also empowers the user to do more than what was possible with a paper document because it can interface with other electronic data sources such as balances, CDS, LIMS, etc.³ For example, it can automatically record balance measurements during the standard and sample preparation stage; hence, it reduces the risk of transcription errors (Figure 3). Additionally, assay method reports can be stored in NuGenesis SDMS and can be easily retrieved and added to the content uniformity document, reducing the analyst's documentation burden (Figure 3). Finally, the HPLC sample set can be automatically populated, uploaded to Empower 2, and the results returned again to the Content Uniformity document (this is also possible with other Chromatography Data Systems).

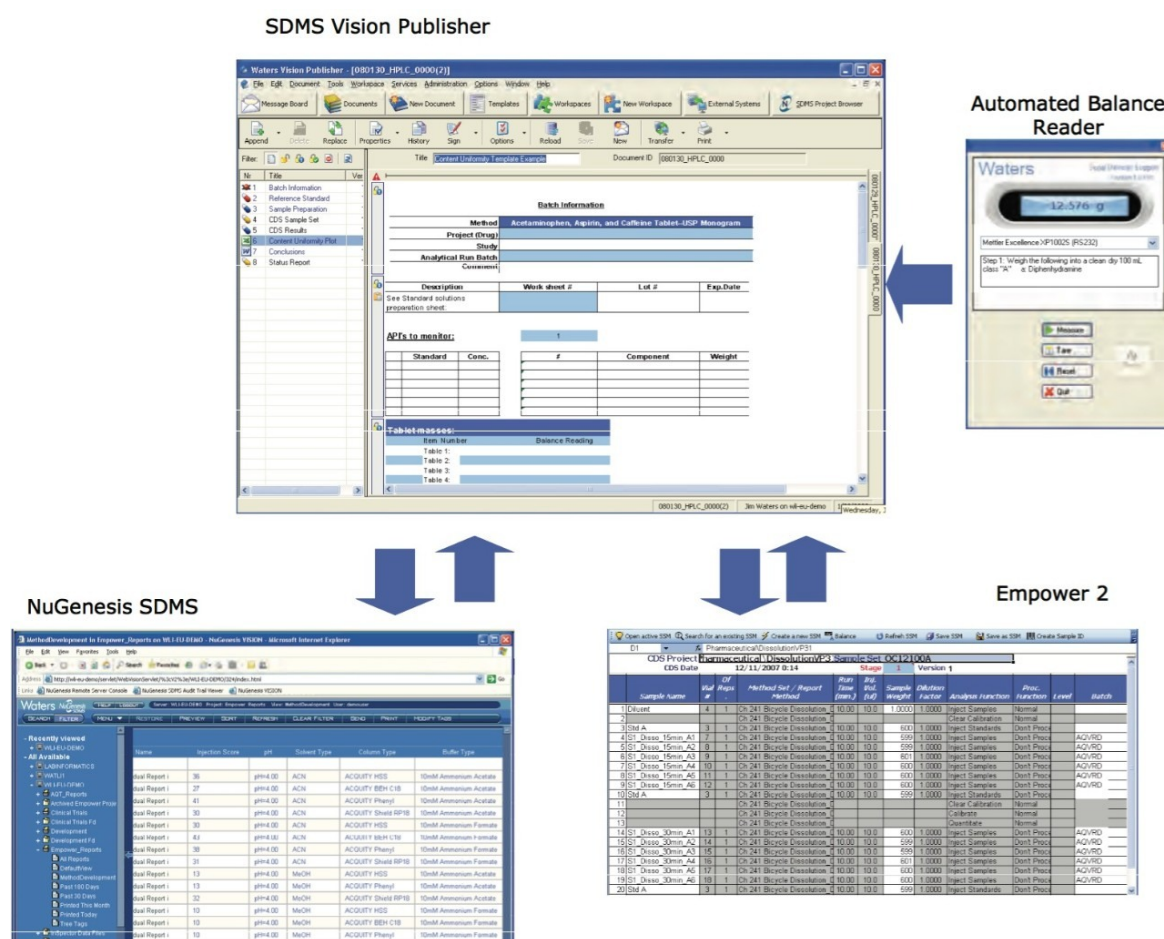


Figure 3. The Content Uniformity document automatically interfaces with external systems such as an automated balance reader, Empower 2, and NuGenesis SDMS. The arrows show the flow of data into, or out of the document.

Conclusion

Content Uniformity is a critical test required before releasing a drug product to market. The process comprises a multi-step workflow where accurate documentation of results is an absolute necessity. The Waters Informatics suite provides an integrated solution that can dramatically streamline workflow, increase laboratory productivity, and facilitate product development and release.

References

1. 2008 USPC Official Monographs: Acetaminophen, Aspirin, and Caffeine.
2. 2008 USPC, 905: Uniformity of Dosage Units.
3. Ping Du, *et al J Assoc Lab Autom.* 12(3): 157–165, 2007.

Featured Products

- [NuGenesis Lab Management System <https://www.waters.com/10067099>](https://www.waters.com/10067099)
- [Empower Chromatography Data System <https://www.waters.com/513188>](https://www.waters.com/513188)

720002554, April 2008



© 2021 Waters Corporation. All Rights Reserved.