

## Note d'application

# Multi-Residue Analysis of Pesticides in Oranges Using AOAC QuEChERS Method by GC-MS

---

Waters Corporation

This is an Application Brief and does not contain a detailed Experimental section.

---

## Abstract

This application brief highlights about Multi-Residue analysis of pesticides in oranges using AOAC QuEChERS method by GC-MS

---

## Experimental

### Extraction Procedure

1. Add 15 mL 1% acetic acid in acetonitrile into the 50 mL DisQuE extraction tube 1.
2. Add 15 g of homogenized sample into the 50 mL tube.
3. Add any internal standards and standard mixture.
4. Shake vigorously for 1 minute and centrifuge > 1500 rcf for 5 minute.

5. Transfer 1 mL of the acetonitrile extract into the clean-up tube 2.
6. Shake for 30 seconds and centrifuge >1500 rcf for 1 minute.
7. Transfer 0.5 mL extract into a tube.
8. Add any post-extraction internal standards.
9. Add 0.25 mL toluene.
10. Evaporate at 50 °C with N<sub>2</sub> to < 0.1 mL.
11. Bring volume up to 0.2 mL with toluene.
12. Transfer to vial with insert for analysis.


## GC Conditions

Instrument:	Agilent 6890N GC
Column:	RTX-5MS, 30 x 0.25 mm, (0.25 µm film)
Carrier Gas:	Helium
Flow Rate:	1.0 mL/min
Temp. Program:	Initial 100 °C, hold 1 min, then 10 °C/min to 320 °C, hold for 7 minutes
Injection Volume:	2 µL splitless

## MS Conditions

Instrument:	Waters Quattro micro GC-MS
Ionization:	Electron Impact (70 eV)
Acquisition:	Single Ion Recording (SIR)

## Results and Discussion

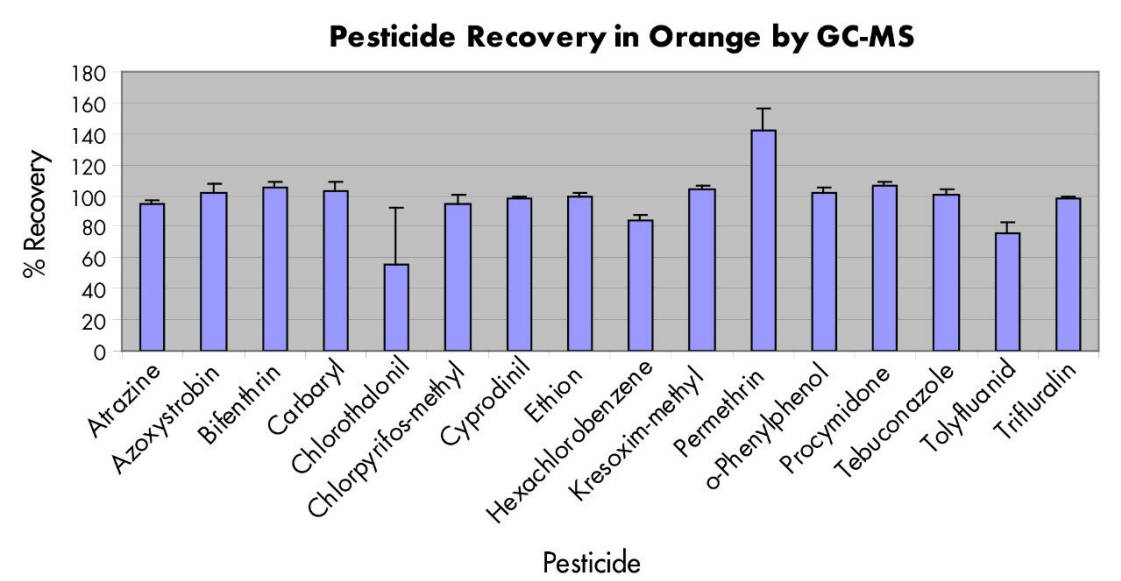

*Pesticides in Oranges by GC-MS*

## Featured Products

©2019 Waters Corporation. All Rights Reserved.