

Nota de aplicación

Improving Resolution Using eXtended *P*erformance (*XP*) Columns

Mia Summers, Kenneth D. Berthelette, Kenneth J. Fountain

Waters Corporation



This is an Application Brief and does not contain a detailed Experimental section.

Abstract

This application brief demonstrates improved resolution of *XP* 2.5 μm Columns over traditional HPLC particle size columns for challenging separations.

Benefits

Improving the resolution of a related-compounds separation using eXtended Performance (*XP*) columns.

Introduction

It is widely accepted that transferring methods to smaller particle sizes can result in faster analysis time. By transferring a method directly to a smaller particle size, there may also be improvements in resolution. As the particle size gets smaller, however, the back pressure across the column will increase. While the use of sub-2- μm columns may necessitate the use of a UPLC System, HPLC users can still realize significant benefits in resolution by transferring their HPLC methods to a eXtended Performance (*XP*) 2.5 μm Column. This may be particularly beneficial for the separation of complex mixtures, where the added resolution from a smaller particle size column may help to identify impurities or target compounds without resorting to increasing column length and run times.

An example of improved resolution using an *XP* 2.5 μm Column is demonstrated using a related compounds mixture of abacavir. Abacavir is a nucleoside reverse-transcriptase inhibitor that is used in anti-HIV therapy. The mixture of related compounds contains five compounds, including the main component, abacavir, shown here in glutarate form (Figure 1). The separation of abacavir from its *trans*-isomer is particularly challenging. Here, the overall improved separation of abacavir from its related compounds is demonstrated, comparing the use of a 3.5- μm column to a high efficiency *XP* 2.5 μm Column.

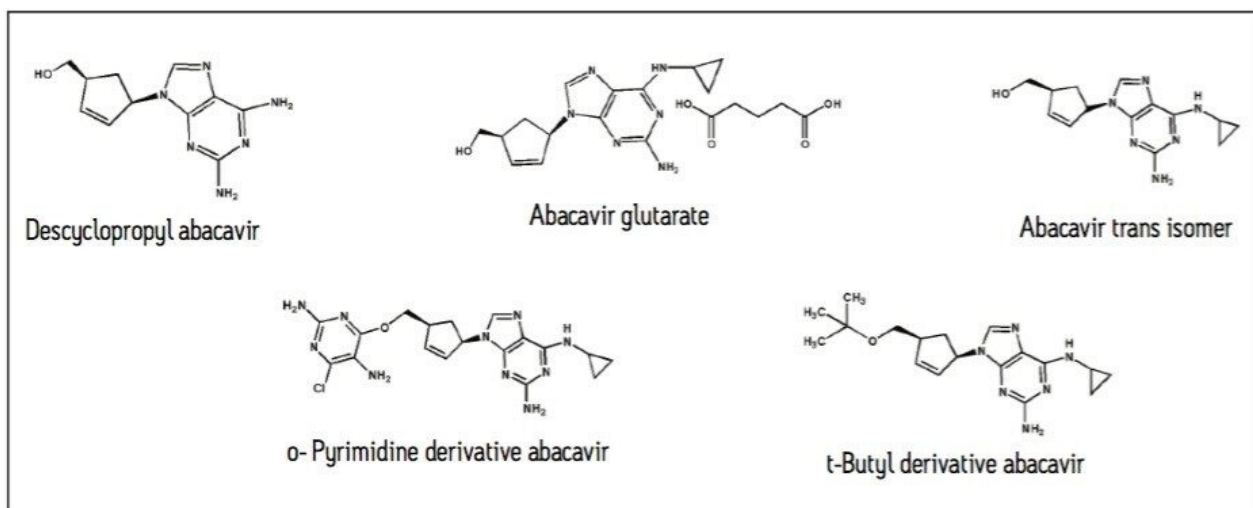


Figure 1. Abacavir components in the USP related compounds mixture.

Results and Discussion

To properly separate related compounds while minimizing extensive method development in HPLC, a highly efficient column with higher resolving power should be used. *XP* Columns contain 2.5- μm particles packed at high pressures in UltraPerformance hardware. The back pressure allowances of the *XP* 2.5 μm particle column still allow for use on an HPLC system.

To demonstrate the improvement in performance using *XP* Columns, the related compounds mixture for abacavir was tested on a 100-mm 3.5- μm , XSelect CSH C_{18} Column. The same method was run using the same column chemistry and dimensions with *XP* 2.5 μm Columns. The comparative separations are shown in Figure 2.

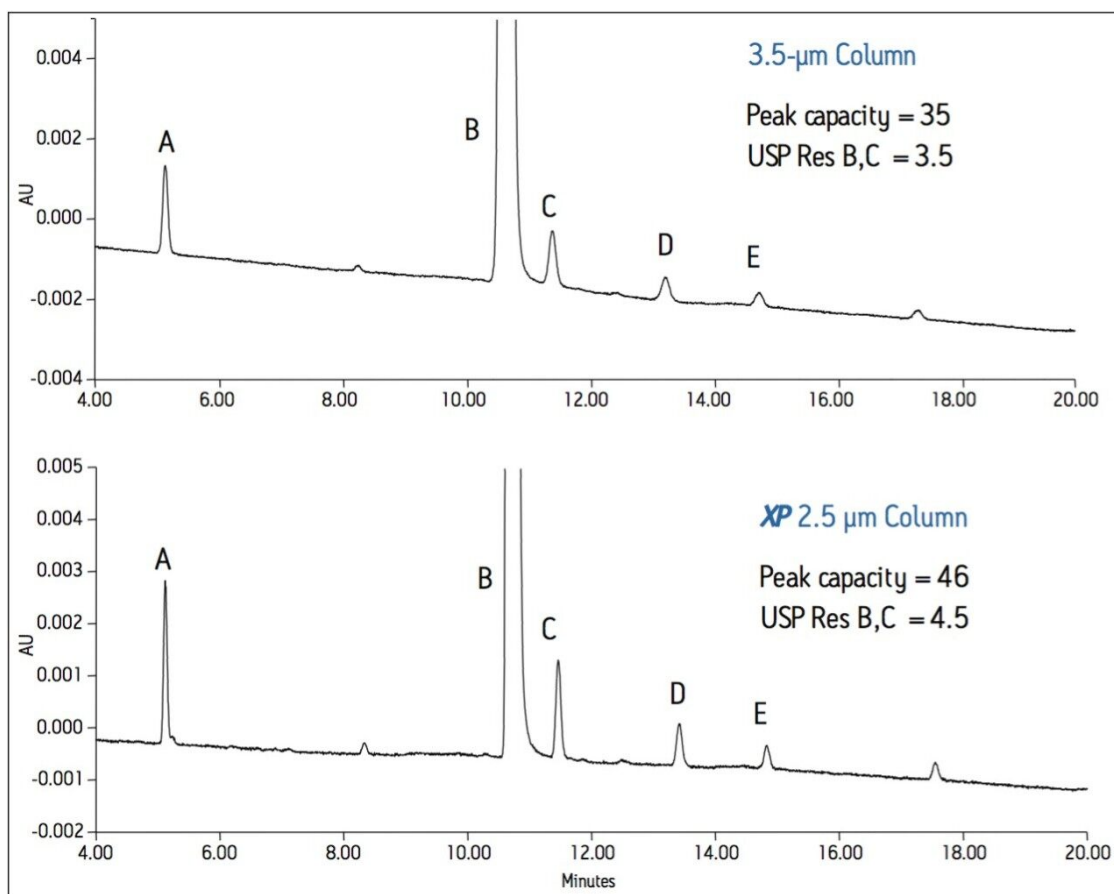


Figure 2. Separation of abacavir related compounds, demonstrating improved assay performance using an eXtended Performance (XP) Column. A) Descyclopropyl abacavir, B) Abacavir glutarate, C) 1R,4R Trans abacavir, D) o-Pyrimidine abacavir, E) t-Butyl abacavir.

In this case, by simply changing the 3.5-μm column to an XP 2.5 μm Column, significant performance improvements are seen as the overall peak capacity for the separation increases 31%, and the peak heights increase up to 42%. A 28% increase in resolution demonstrates the improved separation between closely eluting compounds abacavir and *trans*-abacavir. This example illustrates the capability to increase sensitivity and resolution by using an XP Column, which can result in more accurate identification and quantification of target compounds such as impurities.

Conclusion

By transferring HPLC methods to an *XP* 2.5 μm Column, improvements in resolution and sensitivity can be achieved. This was demonstrated with a related compounds method for abacavir, in which an overall improvement of 31% in peak capacity, up to 42% in peak height, and 29% in resolution were observed by changing the 3.5- μm column to an *XP* 2.5 μm Column. The use of *XP* Columns allows HPLC users to maximize the separation performance on their HPLC systems, lessening the need for further method development, and promoting cost-effective asset utilization.

720004706, May 2013

©2019 Waters Corporation. All Rights Reserved.