


## UPC<sup>2</sup>-MS for Characterization of Complex Oligomeric Materials

---

Baiba Cabovska, Michael O'Leary

Waters Corporation

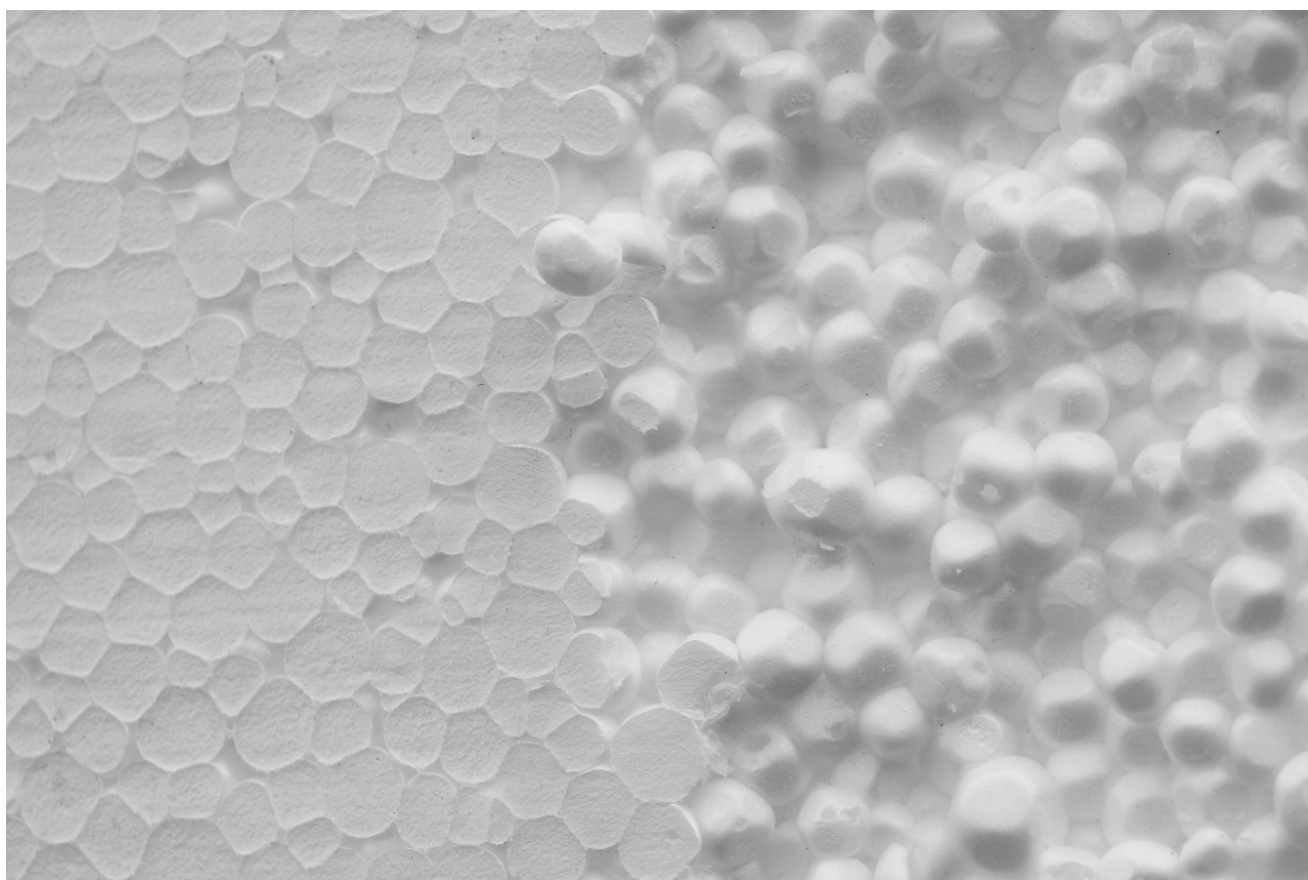

### Abstract

In this application note, we investigate various addition polymers, polystyrenes (PS) and

---

polymethylmethacrylates (PMMA)s, to evaluate the separations range of UltraPerformance Convergence Chromatography. This knowledge was then applied to the analysis of condensation co-polymers, bisphenol A-formaldehyde condensation polymer (PBAA) and poly[(phenyl glycidyl ether)-co-formaldehyde] (PGEF), using MS and UV detection.

## Benefits

- ACQUITY UPC<sup>2</sup> can analyze thermally labile and higher molecular weight polymers than GC.
- ACQUITY UPC<sup>2</sup> is compatible with polar and non-polar polymers.
- Supercritical fluid mobile phase and sub-2- $\mu$ m particle stationary phases shorten retention times for large molecular weight compounds.
- ACQUITY UPC<sup>2</sup> reduces the use of toxic solvents in comparison with normal phase LC.
- MS provides complimentary information to UV data, which can be useful for characterization of individual oligomers, impurity determination, and formulation analysis.

## Introduction

The most common polymer analyses use gel permeation chromatography (GPC) to determine average molecular weight and polydispersity. However, when high resolution separations for individual oligomers are required to evaluate the material performance or understand polymer structure, other analytical techniques are used.<sup>1-4</sup> Low molecular weight polymers can be analyzed by liquid chromatography (LC), gas chromatography (GC), and supercritical fluid chromatography (SFC) among others. The choice of separation technique usually is defined by solubility, average molecular weight, and thermal stability of the polymer. Waters UltraPerformance Convergence Chromatography (UPC<sup>2</sup>), the next step in the evolution of SFC, offers several advantages for the separation of complex oligomeric materials. Due to the low viscosity of supercritical carbon dioxide in comparison with liquids, higher flow rates can be used, which results in shorter analysis times than LC. Convergence chromatography operates at lower temperatures than GC, which is beneficial for the analysis of thermally labile material. Furthermore, UPC<sup>2</sup> can separate higher mass, non-volatile oligomers than GC. Another advantage is the use of sub-2- $\mu$ m particle columns which provide more theoretical plates and better resolution than traditional SFC. If the polymer has a chromophore, UV detection can be used. If information about isomer molecular weight is needed, a mass spectrometer (MS) can be used as the detector. UPC<sup>2</sup> can be interfaced with both UV and MS detectors.

The simplest types of polymers are addition polymers. They are formed by sequential addition of monomer units without a loss of any molecules. Condensation polymers are formed in a condensation reaction between two or more different monomers, where individual molecules bind together and expel a by-product such as water. During polymerization reactions the individual molecules can attach to each other not only linearly but can also form branched isomers. Due to their ability to form various isomers, their separation and characterization can be challenging. In addition, degradation products and by-products can be formed under polymerization conditions which need to be characterized. The performance of polymeric materials can be affected by the isomer and oligomer distribution.

In this application note, we investigated various addition polymers, polystyrenes (PS) and polymethylmethacrylates (PMMA), to evaluate the separations range of UPC<sup>2</sup>. This knowledge was then applied to the analysis of condensation co-polymers, bisphenol A- formaldehyde condensation polymer (PBAA) and poly[(phenyl glycidyl ether)-co-formaldehyde] (PGEF), using MS and UV detection.

## Experimental

### Sample preparation

All polymer samples were dissolved in tetrahydrofuran (THF) at a concentration of 10 mg/mL.

#### UPC<sup>2</sup> conditions

System:	ACQUITY UPC <sup>2</sup> with PDA and ACQUITY SQD
Mobile phase A:	CO <sub>2</sub> (food grade)
Mobile phase B:	0.3% ammonium hydroxide in methanol
Column temp.:	60 °C
Injection volume:	1.0 µL
MS ionization:	ESI (+ or – depending on sample)

## UPC<sup>2</sup> conditions

MS scan range:	150 to 2000 <i>m/z</i>
Capillary:	1 kV
Cone:	25 V
Make-up solvent:	0.3% ammonium hydroxide in methanol
ABPR:	see specific sample
Flow rate:	see specific sample
Vials:	Clear Glass 12 x 32 mm Screw Neck Vial, 2-mL volume
Data management:	Empower 3 CDS

## Results and Discussion

Various molecular weight polystyrenes and PMMAs (Figure 1) were evaluated on sub-2- $\mu$ m particle size columns for UPC<sup>2</sup>. Figure 2 shows the separation of three different polystyrenes. Separation of all oligomers for PS-1000 and PS-1300 was achieved in less than 2.5 minutes. However, only partial separation for PS-2500 was attained. As molecular weight increases, the complexity of the polymer increases so that baseline resolution is no longer achieved.

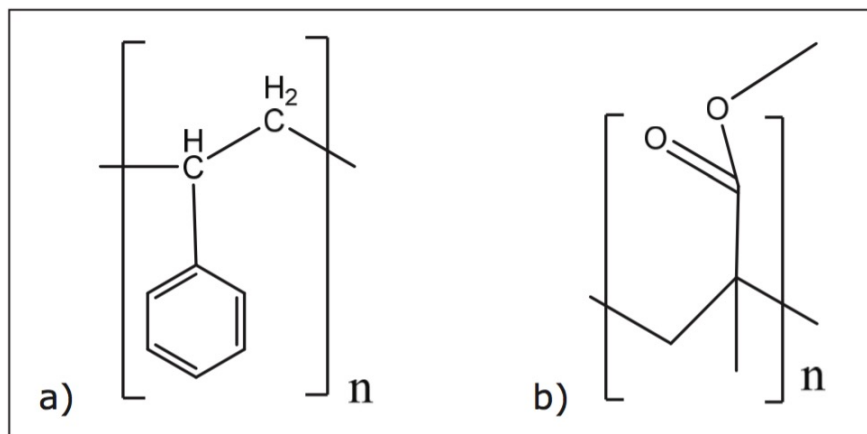


Figure 1. Structures for: a) polystyrene and b) PMMA.

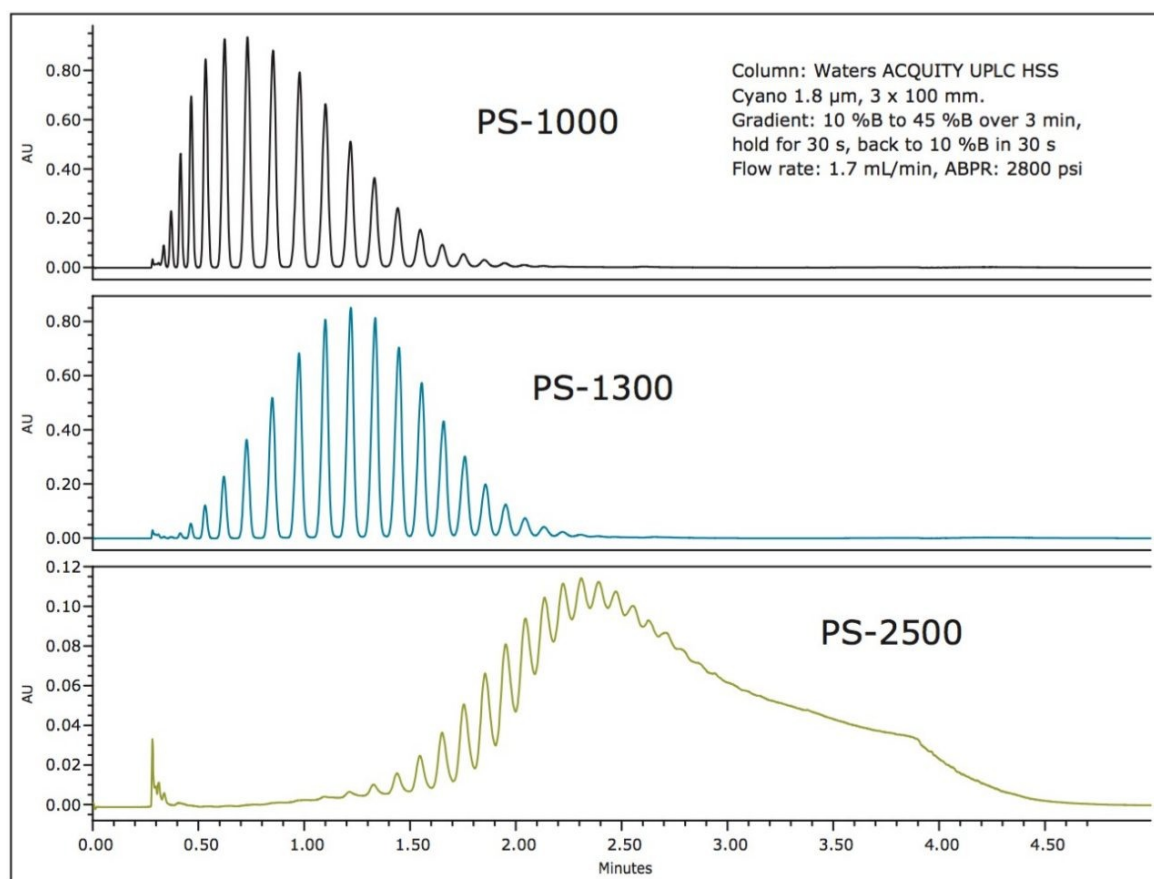


Figure 2. UV chromatograms showing separations for three different polystyrenes.

It was possible to resolve higher mass PMMA oligomers than polystyrene, as shown in Figure 3. As the average molecular weight of polymer increases, the retention time for complete elution goes up as well. The molecular weight range of polymers that can be analyzed by UPC<sup>2</sup> will depend on the solubility of the sample in CO<sub>2</sub>, the type of polymer, and the length of run time the analyst is willing to accept for the separation to be achieved.

Higher molecular weight polymers generally require a higher percentage of organic modifier to elute off the

column. However, increasing the organic modifier percentage increases the back pressure. Keeping the back pressure within an acceptable range requires the flow rate to be decreased, which ultimately can extend the run time.

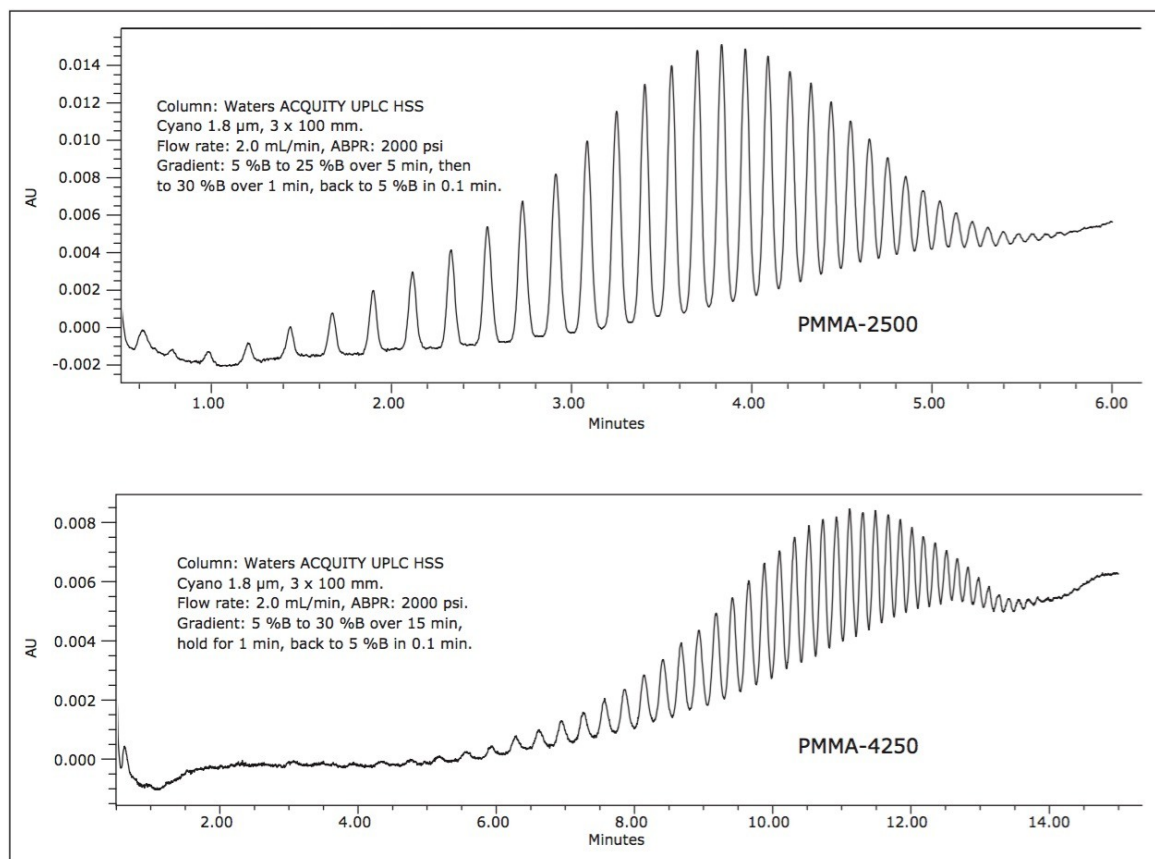


Figure 3. UV chromatograms for PMMA separation.

## Case study 1

Examples of the value of UPC<sup>2</sup> in polymer analysis are demonstrated in the following two case studies. The first one involves the analysis of bisphenol A- formaldehyde condensation co-polymer (PBAA), shown in Figure 4. The co-polymer is formed by addition of polybisphenol A to formaldehyde and expelling a water molecule in the process. Analyzing PBAA, the expected dimer, trimer, and subsequent oligomer peaks were observed. However, an additional large peak was observed at the retention time of 0.7 minutes with  $m/z$  227 (Figure 5). The starting compounds for the polymer were bisphenol A and formaldehyde.  $m/z$  227 (ESI-) corresponds to the bisphenol A molecular ion  $[\text{M-H}]^-$ .

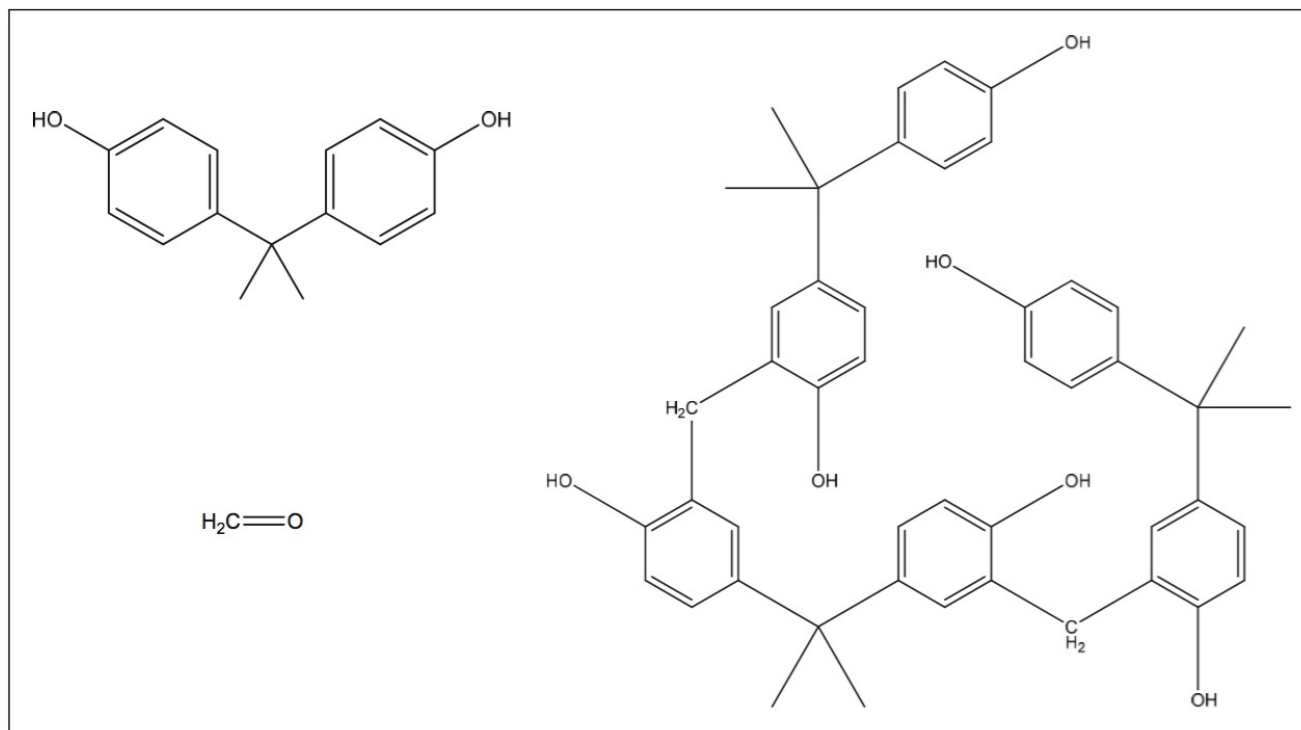


Figure 4. Structures for bisphenol A, formaldehyde, and a trimer of their co-polymer

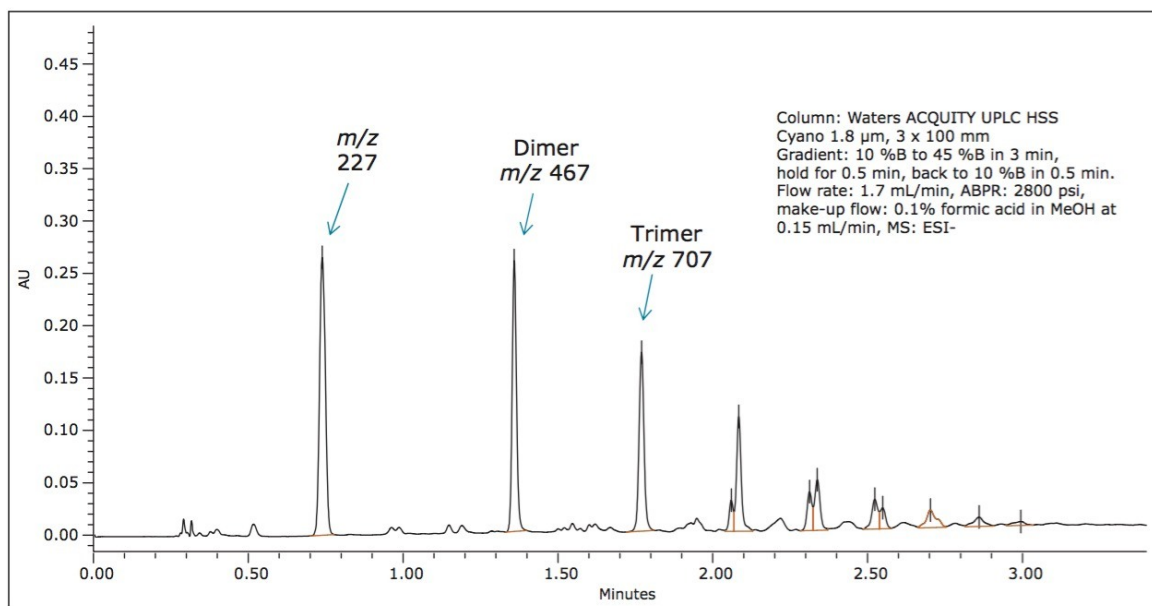


Figure 5. PBA separation.

Confirmation of the unreacted bisphenol A was performed by UV and MS detection with an authentic standard (Figure 6). The retention time of bisphenol A standard corresponded to the unknown peak in the polymer sample. Also, the MS spectrum of bisphenol A matched the spectrum for the peak of interest. Additional confirmation was provided by the formic acid adduct also seen in the mass spectrum.

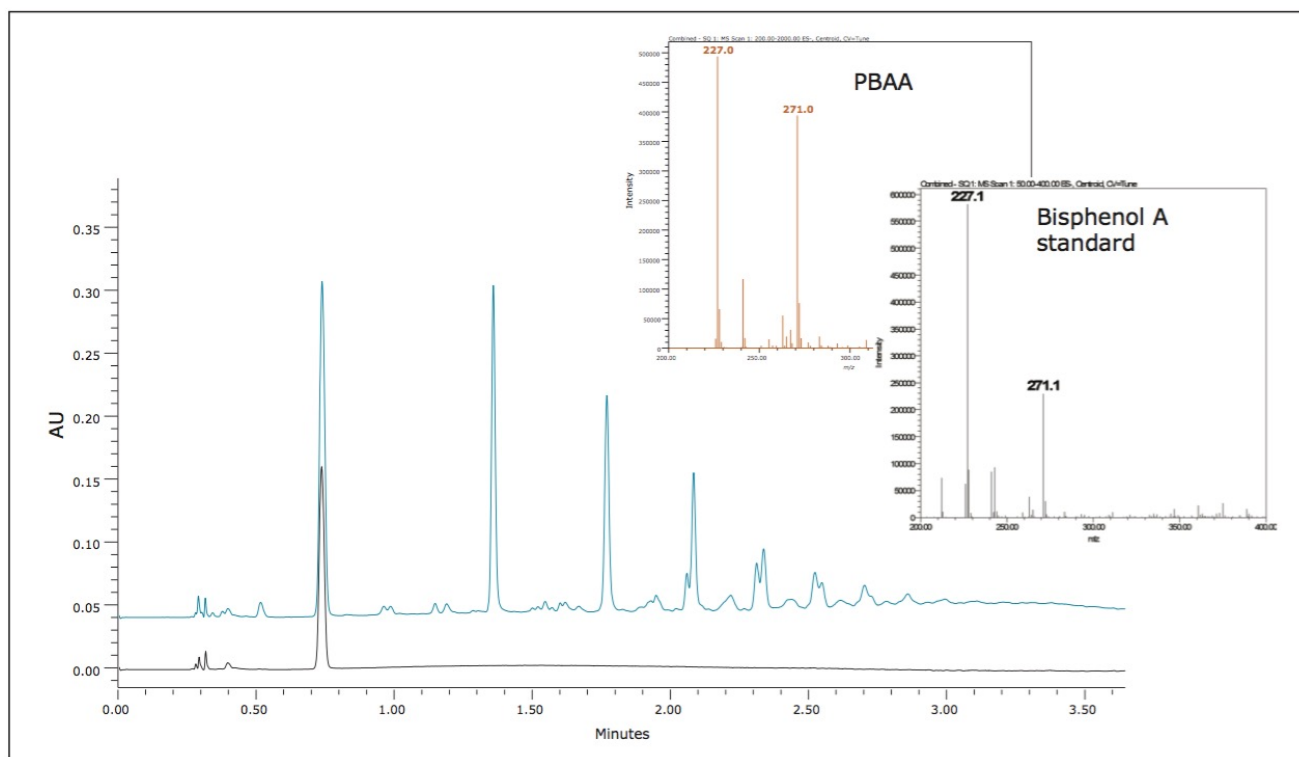


Figure 6. Confirmation for  $m/z$  227 at the retention time of 0.7 minute Black UV trace-bisphenol A standard 0.1 mg/mL, blue UV trace-polymer sample with unreacted bisphenol A.

In this case, MS provided valuable information about an unreacted starting material in the polymerization reaction. This analysis method could be used in reaction monitoring to ensure complete utilization of the starting materials.

## Case study 2

The second test case involved analysis of poly[(phenyl glycidyl ether)-co formaldehyde] (Figure 7). As shown in Figure 8, a separation for individual isomers was easily achieved for dimers. Based on the starting molecule's structure, there are three possible positions for the attachment of the next unit. For the dimer that means six different isomers can be present in the sample – only three were observed. For the trimers and subsequent oligomers, the possible structures increased exponentially. In the current separation, seven individual trimers were resolved.



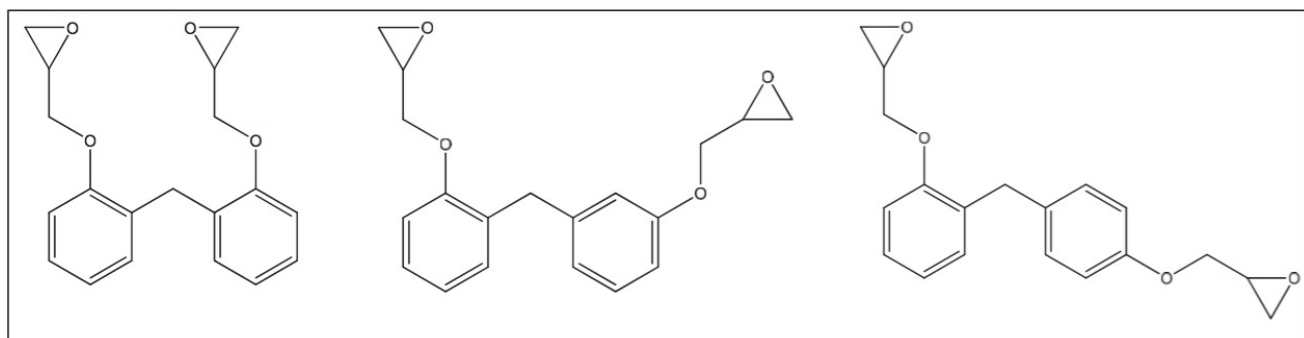


Figure 7. Representative structures for PGEF dimers

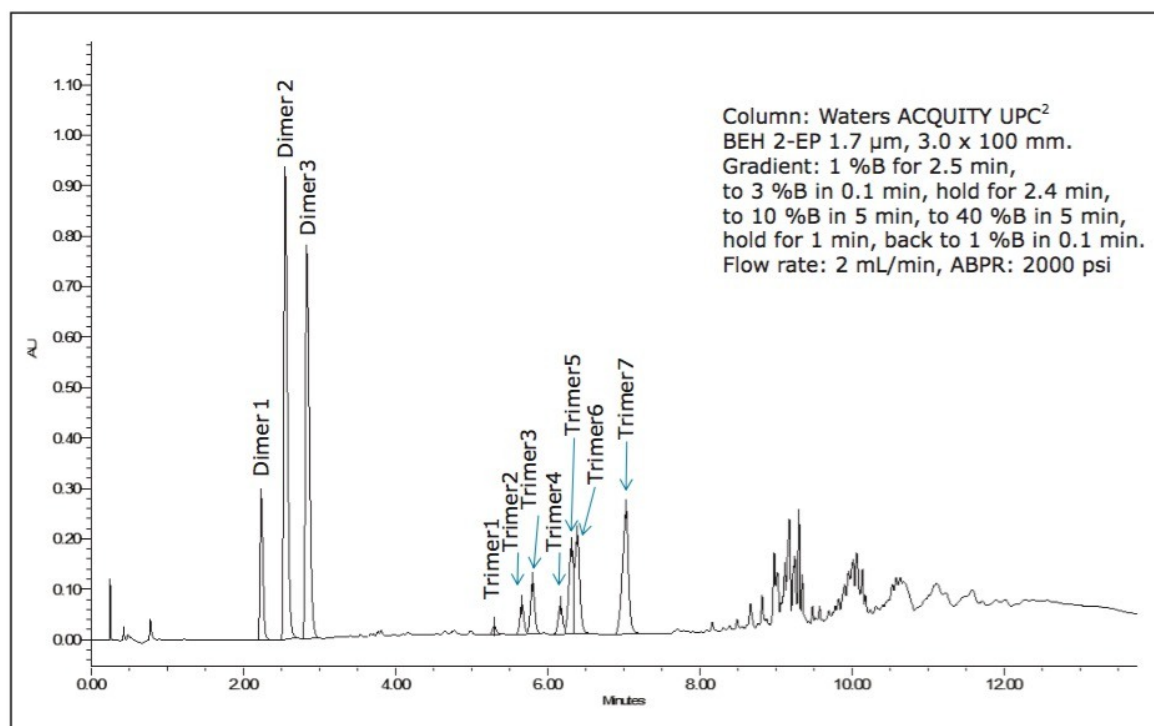


Figure 8. UV chromatograms for the separation of PGEF co-polymer.

When looking at the total ion chromatogram (TIC) from the same separation (Figure 9), additional peaks were seen between the dimers and trimers. The observed  $m/z$  ratios of the clusters were 404 and 402. These masses can result from changes in phenyl glycidyl ether with either loss of a glycidyl ether chain or with opening of an ether ring (Figure 10). Subsequent oligomers containing the described degradation units appeared between trimers and tetramers as well.

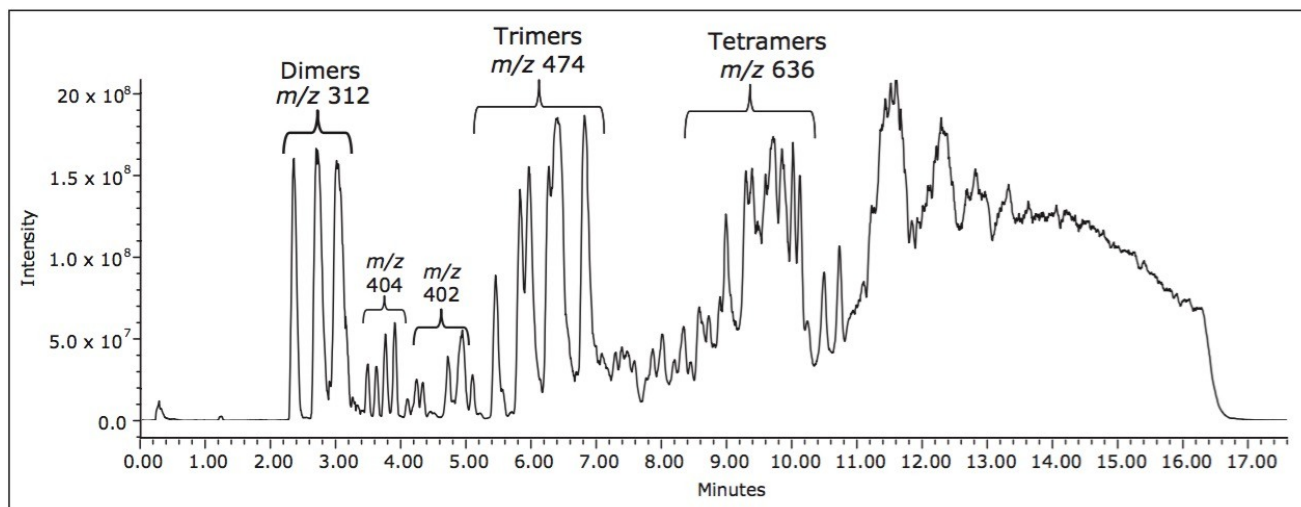


Figure 9. TIC for the separation of PGEF co-polymer.

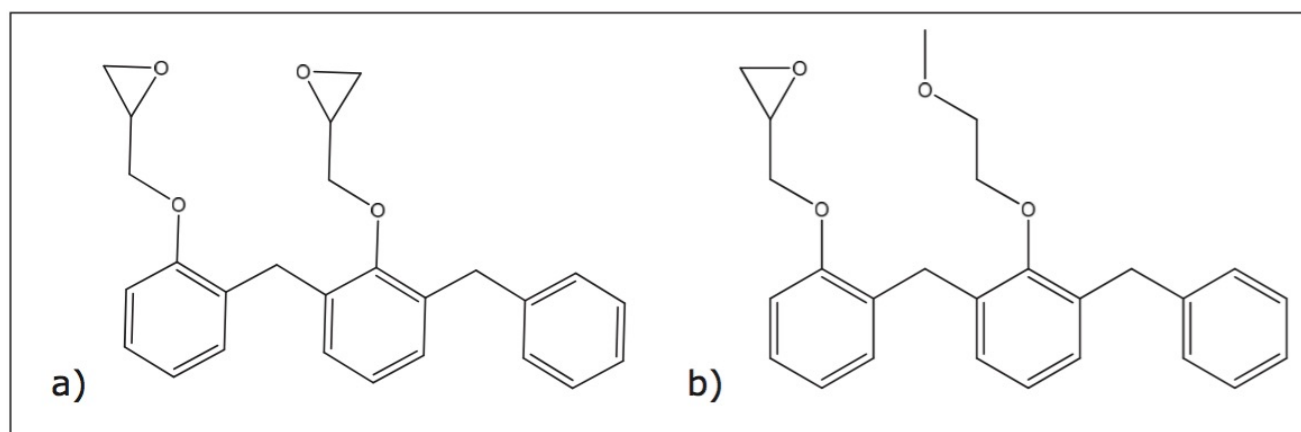


Figure 10. Proposed structures for a)  $m/z$  404 and b)  $m/z$  402.

## Conclusion

UPC<sup>2</sup>-MS is a very powerful tool for characterization of complex oligomeric materials. A wide selectivity space is beneficial for the separation of similar compounds like isomers of polymer oligomers. Additional advantages include compatibility with polar and non-polar polymers, lower analysis temperatures, and higher mass range than GC. The use of supercritical fluid mobile phase shortens retention times for large molecular weight compounds in comparison to LC. The addition of detection by MS provides complementary information to UV data, which can be useful for reaction monitoring, characterization of individual oligomers, impurity determination, and formulation analysis.

## References

1. Ibanez E, Senorans FJ. Tuning of mobile and stationary phase polarity for the separation of polar compounds by SFC. *J Biochem Biophys Methods*. 2000; 43: 25-43.
2. Hoffman BJ, Taylor LT, Rumbelow S, Goff L, Pinkston JD. Separation of derivatized alcohol ethoxylates and propoxylates by low temperature packed column supercritical fluid chromatography using ultraviolet absorbance detection. *J Chromatogr A*. 2004; 1034: 2007-212.
3. Hanton SD. Mass Spectrometry of Polymers and Polymer Surfaces. *Chem Rev*. 2001; 101: 527-569.
4. Takahashi K. Polymer analysis by supercritical fluid chromatography. *J Bioscience Bioeng*. In press, available online 5 March 2013.

## Featured Products

- [ACQUITY UPC2 System <https://www.waters.com/134658367>](https://www.waters.com/134658367)
- [ACQUITY UPLC PDA Detector <https://www.waters.com/514225>](https://www.waters.com/514225)
- [SQ Detector 2 <https://www.waters.com/134631584>](https://www.waters.com/134631584)
- [Empower 3 Chromatography Data Software <https://www.waters.com/10190669>](https://www.waters.com/10190669)

720004759, July 2013



©2019 Waters Corporation. All Rights Reserved.