

## Robustness of ionKey/MS System in the Analysis of Pharmaceutical Compounds in Biological Fluids

---

Paul D. Rainville, Catalin E. Doneanu

Waters Corporation



### Abstract

---

This work illustrates the robustness of the ionKey/MS System when analyzing complex biological samples produced by complex biological samples produced by common bioanalytical sample preparation techniques. We demonstrate the ionKey/MS System's robustness in the analysis of plasma samples prepared by crude protein precipitation (PPT), liquid-liquid extraction (LLE), and by immunoaffinity isolation followed by trypsin digestion (IA/TD).

## Benefits

This work illustrates the robustness of the ionKey/MS System when analyzing complex biological samples produced by common bioanalytical sample preparation techniques.

## Introduction

The robustness and reliability of pharmacokinetic (PK) data is an essential part of bioanalysis. LC-MS is the technique of choice in quantitative bioanalysis due to the high selectivity and sensitivity the technique offers. The use of reduced-bore chromatographic column dimensions such as 180–300  $\mu\text{m}$  I.D. for the analysis of biological fluids can be traced back to the late 1990's by Fraser et al. and has been shown as a means to increase assay sensitivity as well as reduce the amount of sample required to perform drug metabolism and pharmacokinetic (DMPK) analysis.<sup>1,2</sup> However, the implementation of these scales of chromatography require: specialized equipment, smaller tubing I.D.s, and connections-related dead volumes in order to achieve the best chromatographic performance. Plugging of the small scale chromatographic components by the relative dirtiness of biological samples can be a major concern when implementing this scale of chromatography. In this work, we present the robustness in the utilization of the novel ionKey/MS System for analysis of plasma samples prepared by crude protein precipitation (PPT), liquid-liquid extraction (LLE), and by immunoaffinity isolation followed by trypsin digestion (IA/TD).

## Experimental

### UPLC conditions

LC system:

ACQUITY UPLC M-Class

Separation device:	iKey Peptide BEH 130Å, 1.7 µm, 150 µm x 50 mm (p/n 186006764)
iKey temp.:	40 °C
Sample temp.:	4 °C
Injection vol.:	1 µL
Flow rate:	4 µL/min
Mobile phase A:	0.1 % Formic acid in water
Mobile phase B:	0.1% Formic acid in Acetonitrile
Gradient (PPT):	5% B to 95% B over 5 minutes
Gradient (LLE):	5% B to 95% B over 5 minutes
Gradient (IA/TD):	0% B to 30% B over 5 minutes

## MS conditions

MS system:	Xevo TQ-S
Ionization:	Positive ESI
Acquisition mode:	MRM
Capillary voltage:	3.2 kV

Cone voltage and collision energies were optimized for each compound.

## Data management

Chromatography software:

MassLynx 4.1

## Sample preparation

### Protein precipitation (PPT)

Human plasma was prepared by the addition of acetonitrile in a ratio of 2:1 (acetonitrile:plasma). The plasma sample was then vortex mixed for one minute and subsequently centrifuged at 5,000 relative centrifugal force (RCF) for five minutes. The supernatant was then removed, pipetted into an LC vial, and injected onto the LC-MS system. At regular intervals of fifty injections, a QC standard, consisting of dextromethorphan and propranolol, was monitored to assess chromatographic performance over the test period.

### Liquid-liquid extraction (LLE)

Human plasma was prepared by the addition of hexane in a ratio of 10:1. The plasma sample was then vortex mixed for one minute and subsequently centrifuged at 5,000 RCF for five minutes. The supernatant was then removed into a new vial. The sample was then dried down and reconstituted in one fifth the initial volume, and injected onto the LC-MS system. At regular intervals of twenty injections, a QC standard, consisting of dextromethorphan and propranolol, was monitored to assess chromatographic performance over the test period.

### Immunoaffinity isolation and tryptic digestion of a monoclonal antibody (IA/TD)

Samples were kindly obtained from Bristol Myers Squibb (BMS). Human plasma was spiked with a therapeutic monoclonal antibody (mAb) and immunoaffinity isolation, implemented in the magnetic bead format, was used for the isolation of the mAb from plasma. After denaturation, the mAb was digested with trypsin. Over 1,000 injections of the mAb digest were performed and two signature peptides were monitored in each LC-MS run to evaluate the chromatographic performance over the test period.

## Results and Discussion

Routine analysis within a bioanalytical laboratory usually consists of a batch of two to four 96-well sample plates that have been prepared by the bench chemist using protein precipitation, liquid-liquid extraction, or in the case of biopharmaceuticals such as monoclonal antibodies, immunoaffinity isolation followed by proteolytic digestion.

Of these techniques, protein precipitation is the most commonly utilized due to the speed and relative low cost of the technique for the analysis of small molecules.<sup>3</sup> However this technique is also the crudest in its ability to produce clean samples for analysis. Because of this fact, robustness testing of a novel 150  $\mu\text{m}$  iKey Separations Device was carried out under these conditions as they provided the most challenging of the samples preparation techniques of biological fluids. The ionKey/MS System, comprised of the Xevo TQ-S Mass Spectrometer, the ACQUITY UPLC M-Class, the ionKey Source, and the iKey Separation Device is shown in Figure 1. The iKey Separation Device consists of the ceramic-based separations device with an integrated emitter together in a single device that is placed directly into the source of the mass spectrometer.



Figure 1. ionKey/MS System: comprised of the Xevo TQ-S, the ACQUITY UPLC M-Class, the ionKey Source, and the iKey Separation Device.

Figure 2 illustrates the separation of dextromethorphan and propranolol, during a generic five minute LC gradient. Here we observe a peak width at 10% height for dextromethorphan and propranolol of one second. The calculated resolution at this peak height was 0.8. In Figure 2 we again show the injection of the QC standard after 1,000 injections (5 days) of continuous 1  $\mu\text{L}$  injections of the protein precipitated human plasma. It should be noted that a 1  $\mu\text{L}$  injection on the 150  $\mu\text{m}$  I.D. iKey Separation Device is equivalent to 200  $\mu\text{L}$  on a traditional 2.1 mm I.D. analytical column. The comparison of the data shown in Figure 2 indicates that excellent peak symmetry

and chromatographic resolution of 0.9 were maintained over the course of the testing period.

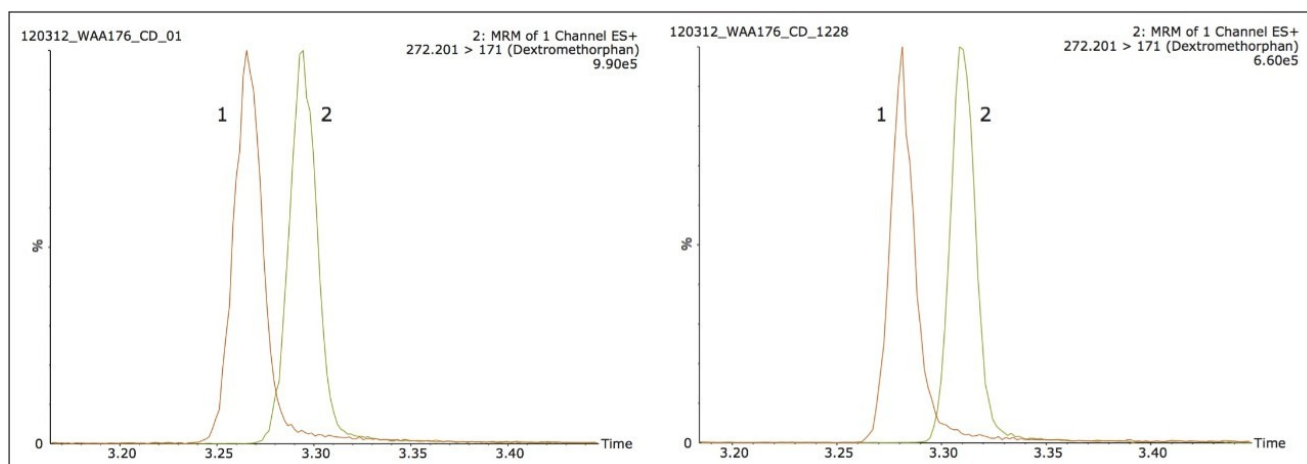


Figure 2. Resolution of propranolol (1) and dextromethorphan (2) at the start and finish of 1,000 continuous 1 µL injections of human plasma, prepared by protein precipitation.

The system pressure traces from the study are shown in Figure 3. Here we observe no discernible increases in the system pressure, indicating that none of the frits, tubing or connective fittings have been blocked over the course of the study. It should again be noted that this injection volume of 1 µL is analogous to injecting roughly 200 times this volume or roughly 200 µL onto a standard 2.1 mm I.D. column. Cleanliness of the MS source is often a key parameter in the continued acquisition of quality data over the course of a study. The inherent ability in the use of smaller volumes to achieve similar sensitivity results as with standard 2.1 mm I.D. scale LC-MS equates to less contamination of the MS source from the sample.

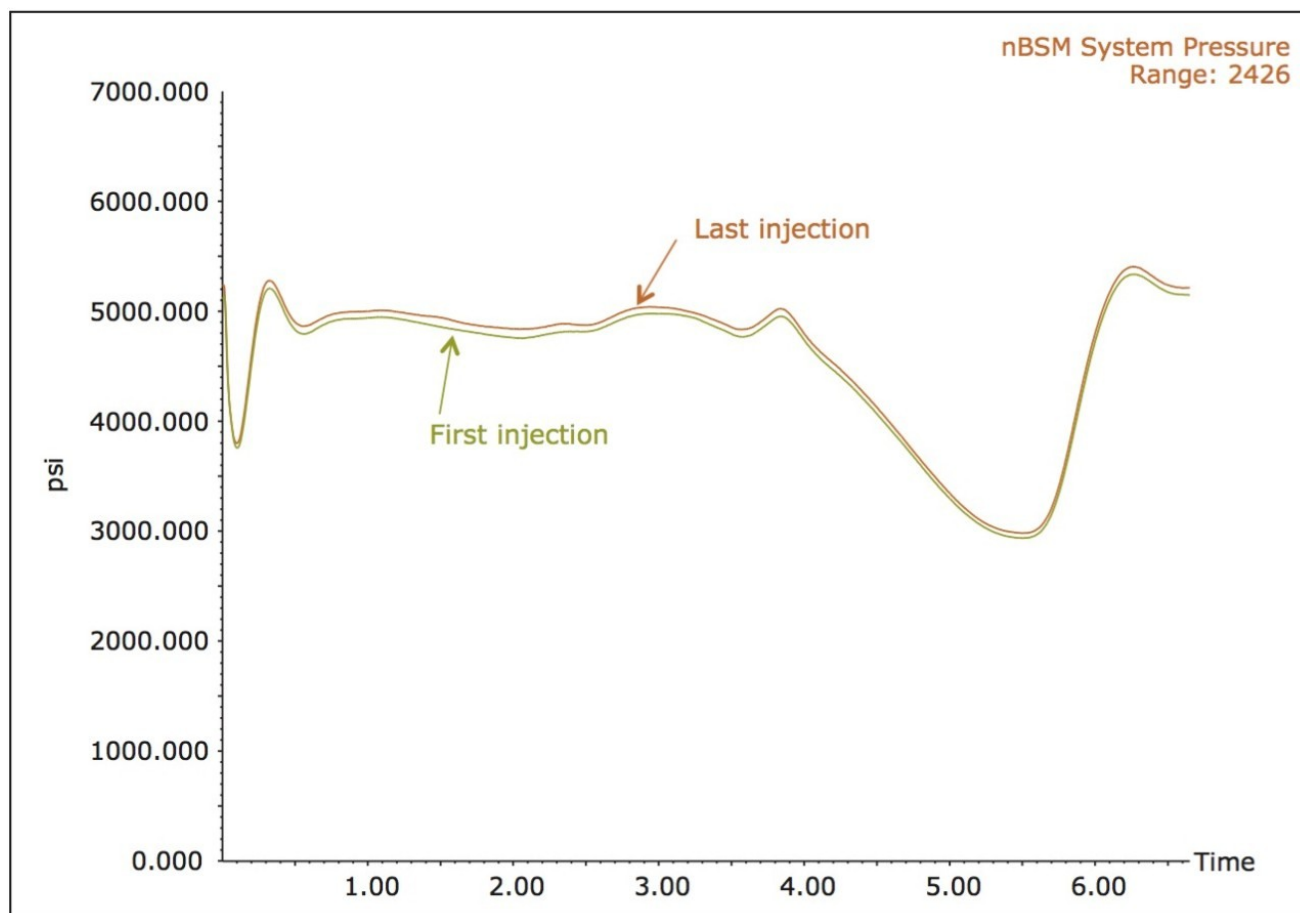


Figure 3. IonKey/MS System pressure traces at injection 1 and injection 1,000, after continuous injections of 1  $\mu$ L of protein precipitated human plasma.

Figure 4 illustrates the stability of the peak area from the QC sample over a course of 1000 injections of the protein precipitated human plasma sample. This data further illustrates the robustness, not only of the iKey Separations Device, integrated emitter, MS source but the entire system. The system was next challenged with human plasma samples prepared by LLE. This test utilized the same experimental conditions as with the previous example. Figure 5 shows the QC standard of propranolol and dextromethorphan at injection 1 and injection 1000 and illustrates that both the chromatographic peak shape and the resolution were maintained over the course of the study, much in the same manner as with the previous example. It should be further noted that again no discernible increase in system backpressure was observed.

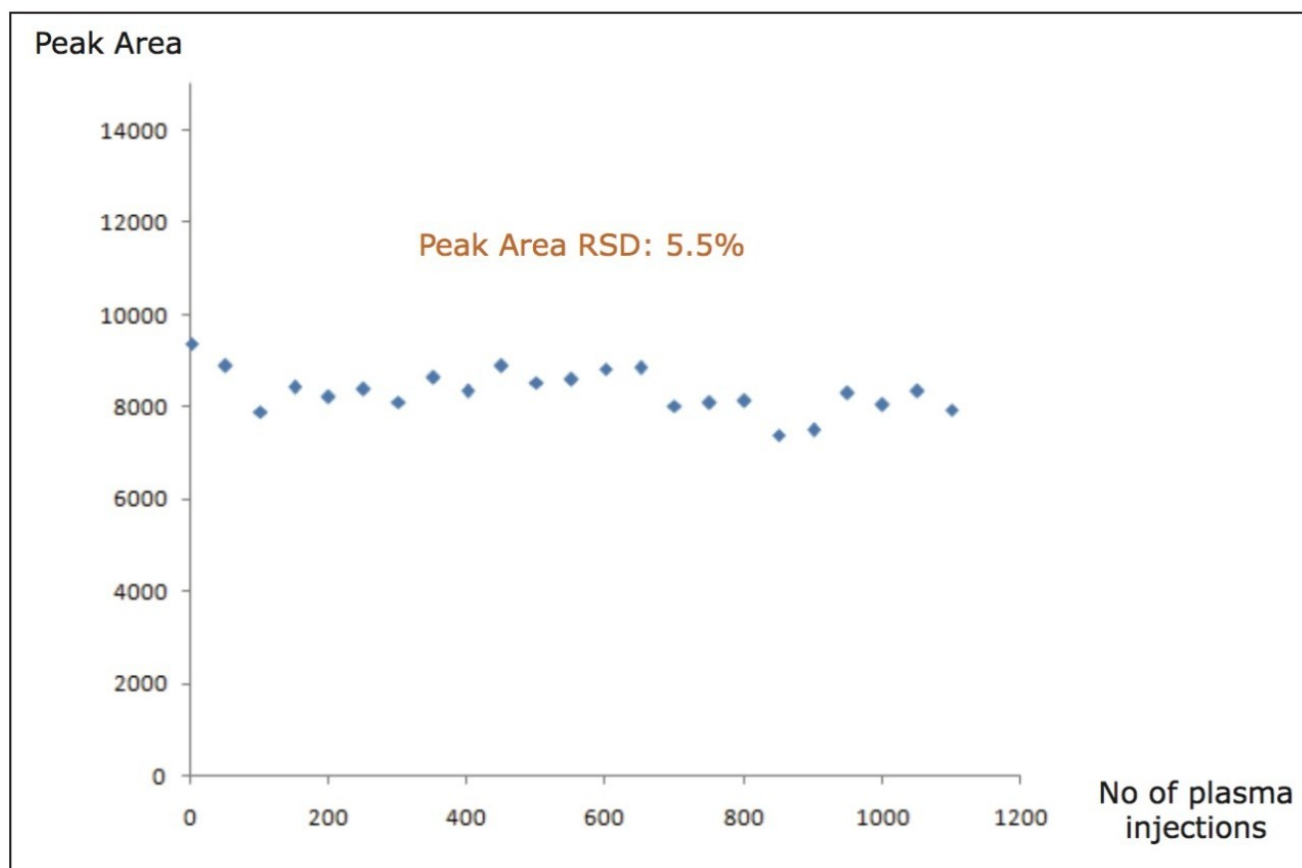


Figure 4. Dextromethorphan peak area counts from various test time points for the QC sample taken during the robustness testing with human plasma prepared by protein precipitation.



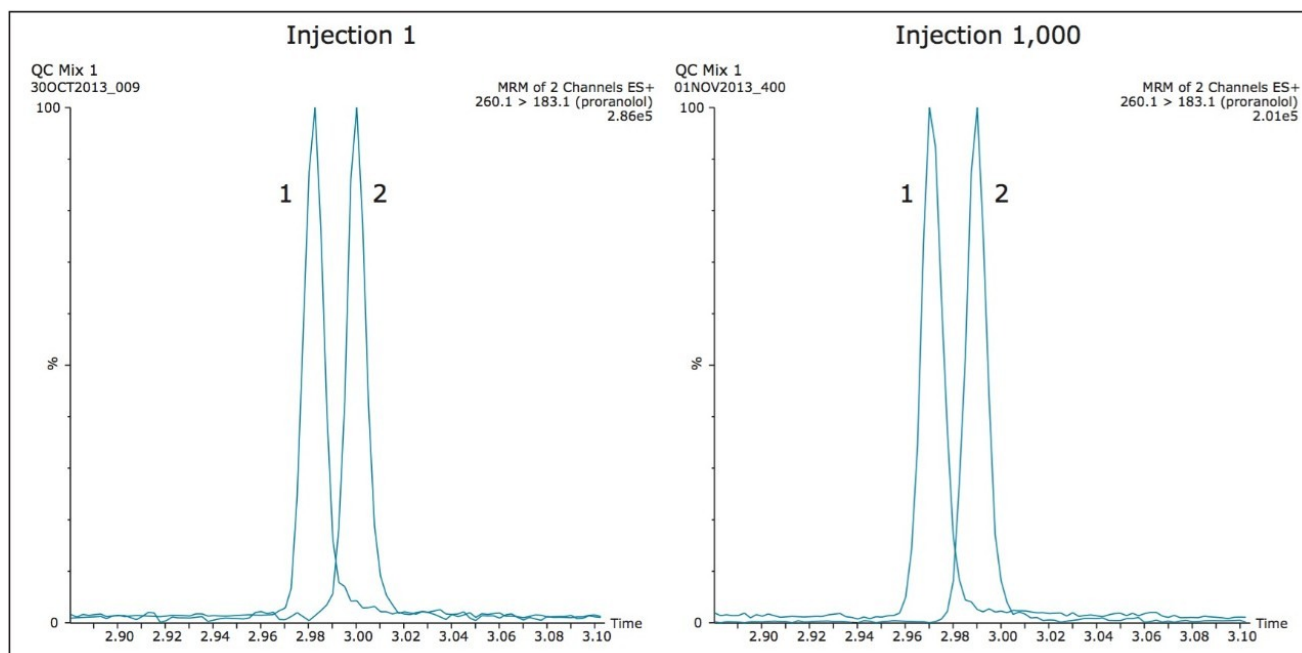


Figure 5. Resolution of propranolol (1) and dextromethorphan (2) at the start and finish of 1,000 continuous 1  $\mu$ L injections of human plasma, prepared by liquid-liquid extraction.

Due to the increase in the development of biopharmaceuticals, such as antibody therapeutics, the robustness of the ionKey/MS System was tested with a common sample preparation scheme of affinity isolation followed by digestion and subsequent analysis of peptides generated from the sample preparation. As with the previous studies chromatographic performance and system pressure were monitored. In this example, specific peptides produced by the trypsin digestion of the antibody were monitored over the course of the study. Figure 6 illustrates the peak shape at injection 1 and injection 1,000 for two of the signature peptides monitored during the study. The typical peptide peak widths (at 10% peak height) observed under the experimental conditions employed were 3–4 seconds.

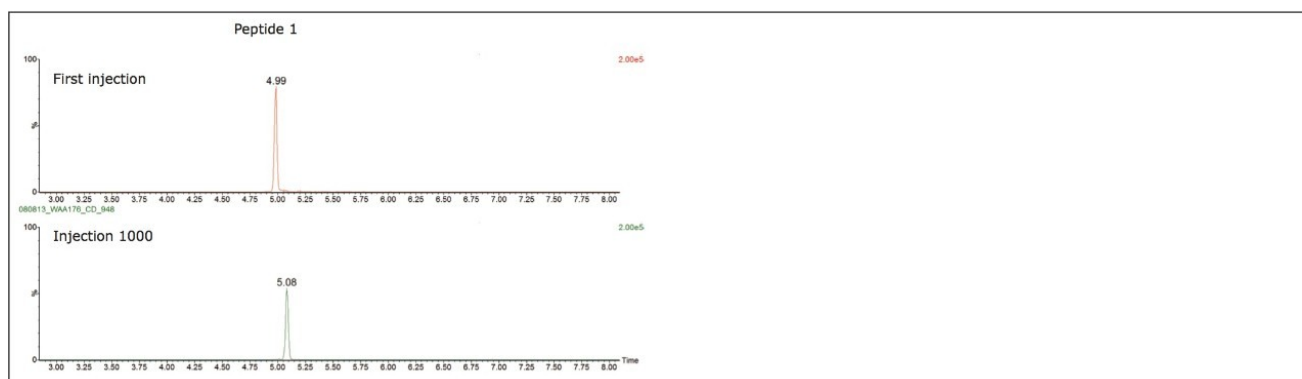


Figure 6. MRM chromatograms of two signature peptides from a therapeutic mAb digested with trypsin following immunoaffinity isolation from mouse serum.

As in the previous experiments, pressure traces were recorded throughout the entire study and the first and the last pressure traces are displayed in Figure 7.

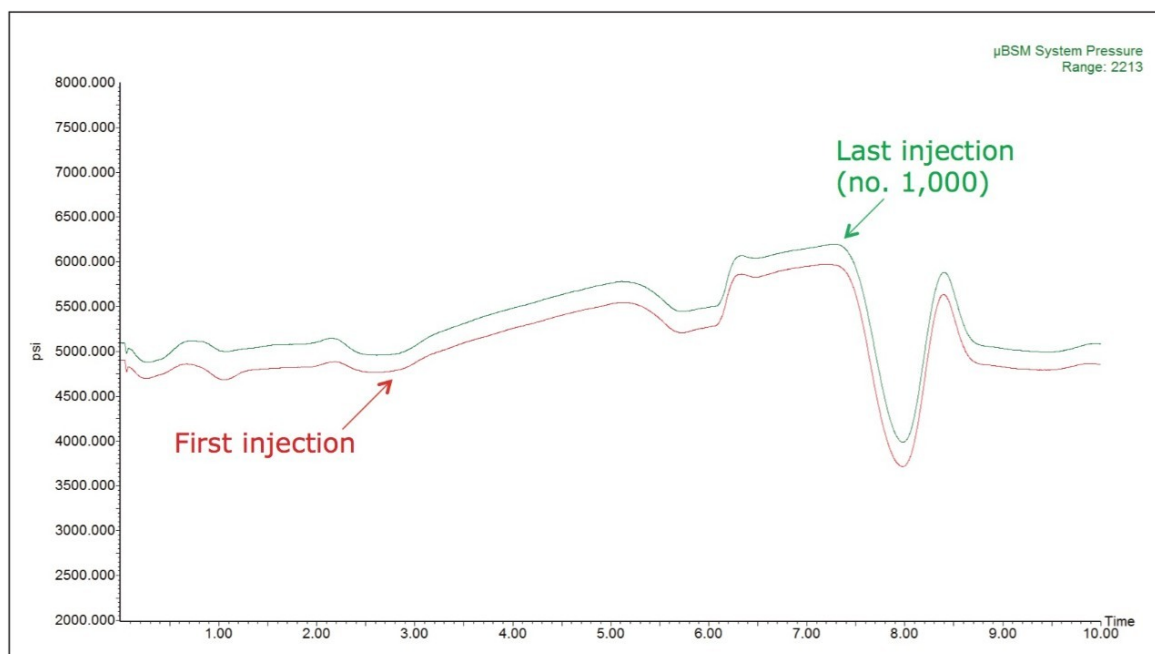


Figure 7. ionKey/MS System pressure traces recorded for the first and the last (1,000th) injection of a mAb isolated by immunoaffinity from mouse serum and digested with trypsin.

## Conclusion

The performance of the ionKey/MS System, using a novel 150  $\mu\text{m}$  I.D. iKey Separation Device packed with 1.7  $\mu\text{m}$  chromatographic particles, in the analysis of biological fluids was shown to be:

- Robust for over 1,000 injections of human plasma prepared by protein precipitation, liquid-liquid extraction, and by immunoaffinity isolation with trypsin digestion, using injection volumes of 1  $\mu\text{L}$ . This volume is roughly equivalent to a 200  $\mu\text{L}$  injection on a standard 2.1 mm I.D. column.
- Capable of maintaining excellent peak symmetry and resolution for a critical pair of small molecules over a continuous testing period of 5 days. The resolution and peak width measured at 10% peak height was 0.8 and 1 second, at injection 1 and injection 1,000 respectively.
- Capable of maintaining very good chromatographic performance in terms of peak shape and peak width over 1,000 injections of a complex mAb digest sample.
- Maintaining consistent system pressure over the course of the study, indicating no accumulation of plasma

components in the narrow I.D. tubing, or frits that could lead to poor chromatographic performance and system over-pressure.

## References

1. Fraser, I., et al., The use of capillary high performance liquid chromatography with electrospray mass spectrometry for the analysis of small volume blood samples from serially bled mice to determine the pharmacokinetics of early discovery compounds, *Rapid Communications in Mass Spectrometry*, 1999, 13, 2366–2375.
2. Dear, G., et al., Applications of capillary LC(-MS) to drug metabolism studies, *LCGC Europe*, 2001, 2–8.
3. Wells, D.A., High throughput bioanalytical sample preparation, 2003, Elsevier, St. Paul, Minnesota, USA.

## Acknowledgements

Waters would like to thank Eugene Ciccimaro, Bogdan Slecicka, Celia D'Arienzo, and Timothy Olah from BMS for the production and use of the digested antibody samples utilized in this study.

## Featured Products

- [ACQUITY UPLC M-Class System <https://www.waters.com/134776759>](https://www.waters.com/134776759)
- [ionKey/MS <https://www.waters.com/134782630>](https://www.waters.com/134782630)
- [Xevo TQ-S <https://www.waters.com/10160596>](https://www.waters.com/10160596)
- [MassLynx MS Software <https://www.waters.com/513662>](https://www.waters.com/513662)

720004953, February 2016



©2019 Waters Corporation. All Rights Reserved.