


Nota de aplicación

# Metabolite Identification of Complex Cyclic Peptides Using WebMetabase, Ion Mobility-Enabled DIA, and Product Ion Confirmation

---

Mark D. Wrona

Waters Corporation

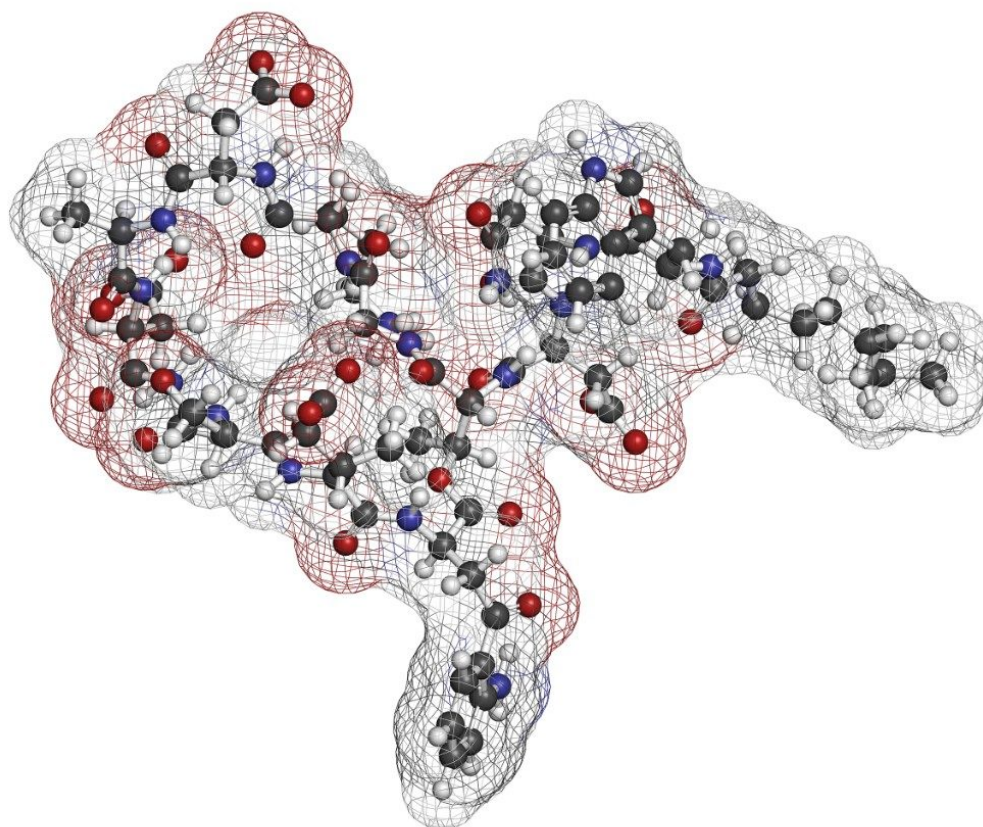

---

## Abstract

In this application note, we describe a solution that uses an information-rich data acquisition mode, which effectively deals with the complexity of modified peptides. This enables key metabolic data to be generated unambiguously while tackling complex peptide studies. The use of DIA combined with ion mobility (HDMS<sup>E</sup>) improves upon the selectivity provided by MS-only techniques. The addition of routine Collision Cross Section (CCS) measurements makes HDMS<sup>E</sup> a powerful identification acquisition mode, as well as an ideal separation technique.

## Benefits

- Ability to screen metabolism from complex cyclic peptides using comprehensive HDMS<sup>E</sup> combined with optional product ion confirmation scan (PICS, HS-MSMS/DDA scans)
- Use of QuanRecovery vials to minimize impact of non-specific losses
- Automatic CCS value generation to help track metabolites across studies
- Cloud based workflow to enable uploading of data to a central server for processing and web based review

---

## Introduction

The incorporation of unnatural amino acids, backbone modifications, conjugations, and cyclizations is a commonly used strategy to improve the efficacy and ADME profile of peptide therapeutics. Analyzing these complex peptides (often natural products which may be synthetically modified) requires overcoming a variety of analytical and technical challenges, including challenging chromatography, non-specific binding, limited or complex fragmentation, and the prediction and review of many catabolites and/or novel metabolites. To ensure safe and efficacious therapies, the metabolic fate of potential drug candidates needs to be thoroughly and rapidly investigated using LC-MS coupled with informatics capable of identifying and characterizing their profiles.

Current approaches for analyzing complex modified peptides suffer from challenges that include:

1. Rules developed for small molecule metabolism which are largely not applicable to or are not observed in

appreciable amounts for peptides.

2. Tracking and identifying all multiply charged species for all potential metabolites.
3. Data Dependant Acquisition (DDA) approaches generate high quality data, but MS-MS spectra may not be triggered on relatively low abundance peaks without multiple runs, longer run times, or detailed inclusion/exclusion rules. DDA may also have an insufficient survey scan frequency to characterize all closely eluting relevant peaks. Data Independent Acquisition (DIA) (all ions) approaches are comprehensive for identification, unlikely to miss precursor information, and enable retroactive analysis, but may not always be sufficient for generating fully resolved product ion spectra for all peaks.
4. Predictive approaches to generate inclusion lists can improve the quality of hits but may still miss peaks due to imperfect prediction of the precursor metabolite masses in earlier studies. Even modest peptide lengths may generate very long lists of potential targets.
5. Non-specific binding, leading to underreporting and loss of some metabolites.

Here, we describe a solution that uses an information-rich data acquisition mode, which effectively deals with the complexity of modified peptides. This enables key metabolic data to be generated unambiguously while tackling complex peptide studies. The use of DIA combined with ion mobility (HDMS<sup>E</sup>) improves upon the selectivity provided by MS-only techniques. The addition of routine Collision Cross Section (CCS) measurements makes HDMS<sup>E</sup> a powerful identification acquisition mode, as well as an ideal separation technique. For those projects requiring definitive product ion scans, we describe the use of simultaneous product ion confirmation scans (PICS, IMS MS/MS) for metabolites with complex product ion spectra to enable the highest confidence in elucidating structures. Loss of metabolites and signal response can be mitigated through the use of low binding protocols or products such as QuanRecovery vials.

All data for daptomycin, dalbavancin, and lanreotide were screened for metabolism using an ion mobility-enabled HDMS<sup>E</sup> acquisition workflow and processed using Mass-MetaSite (MMS) and WebMetabase (WMB) macromolecule software packages (Lead Molecular Design, S.L, Sant Cugat del Vallés, Spain). HDMS<sup>E</sup> and HDMS<sup>E</sup> + product ion data were reviewed using the UNIFI screening application.

---

## Experimental

---

Samples were incubated and processed according to methodology reported by Radchenko et al.<sup>1</sup> Briefly, chymotrypsin (1 mg/mL) (MilliporeSigma, St.Louis, MO) was dissolved in simulated intestinal fluid (SIF) containing pancreatin (Ricca Chemical, Arlington, TX). 50  $\mu$ L incubations were prepared in 300  $\mu$ L QuanRecovery vials (Waters Corporation, Milford, MA). One mM stock solutions of daptomycin, dalbavancin or lanreotide (MilliporeSigma, St.Louis, MO) were prepared. 0, 5, 15, 45, 120, and 300 min timepoints for daptomycin and 0 and 300 min for dalbavancin and lanreotide were prepared at a starting concentration of 10  $\mu$ M. Each timepoint was quenched with two volumes cold acetonitrile containing 1% ammonium hydroxide (by volume).

## Method conditions

### LC conditions

LC system:	ACQUITY UPLC I-Class PLUS with FTN
Vials:	QuanRecovery with MaxPeak
Column:	ACQUITY UPLC Peptide BEH C <sub>18</sub> 300 Å, 1.7 $\mu$ m
Column temp.:	60 °C
Sample temp.:	8 °C
Injection volume:	2 $\mu$ L
Flow rate:	0.5 mL/min
Mobile phase A:	Water, 0.1% formic acid
Mobile phase B:	Acetonitrile, 0.1% formic acid
Gradient:	2 to 60% B over 7 min; 60 to 90% B from 7 to 7.5 min; 90% B hold from 7.5 to 8.5 min; and re-equilibrate at 2% B from 8 to 10 min.

## MS conditions

MS system:	Vion IMS QToF Mass Spectrometer
Ionization mode:	ESI+
Acquisition range:	50–2000 <i>m/z</i>
Capillary voltage	1.5 kV
HDMS <sup>E</sup> collision energy:	4 (low energy) 15–70 eV ramp (high energy)
HS-MSMS collision energy:	12–40 eV ramp @ 50 <i>m/z</i> to 22–80 @ 2000 <i>m/z</i>
Cone voltage:	80 V

## Data management

UNIFI Scientific Information System, version 1.9 SR4 API enabled Mass-MetaSite 3.4.9, WebMetabase 4.03 (Lead Molecular Design, S.L, Sant Cugat del Vallés, Spain)

## Data processing

All data were acquired in UNIFI and processed using the Mass-MetaSite batch processor. For each peptide, time points with blanks and standards were processed automatically (in the US) and uploaded to a WebMetabase server (based in the UK), and reviewed via a web browser interface.

---


## Results and Discussion

### Selection of Test Set, Overall Catabolism of Peptides

The primary goal of this study was to verify that data for complex peptide structures could be rapidly acquired

using HDMS<sup>E</sup> acquisition on a Vion IMS QTof MS and subsequently processed for catabolism identification using Mass-MetaSite and WebMetabase directly using the UNIFI Application Programming Interface (API).<sup>2</sup>

A cross section of peptides were tested containing various non-natural substitutions, adducts, and increasingly complex backbones and cyclizations within the macromolecular structures (Figure 1). For all peptides, the top three metabolites were reported and for daptomycin, a larger discussion of metabolites and CCS values generated is provided. Table 1 shows the complexity observed in the peptides tested and summarizes the structural modifications that can be handled by Mass-MetaSite and WebMetabase. Simpler peptides shown in Table 1, such as GLP-1 and somatostatin, are not discussed here, but have been previously published by several labs.<sup>3,4</sup>

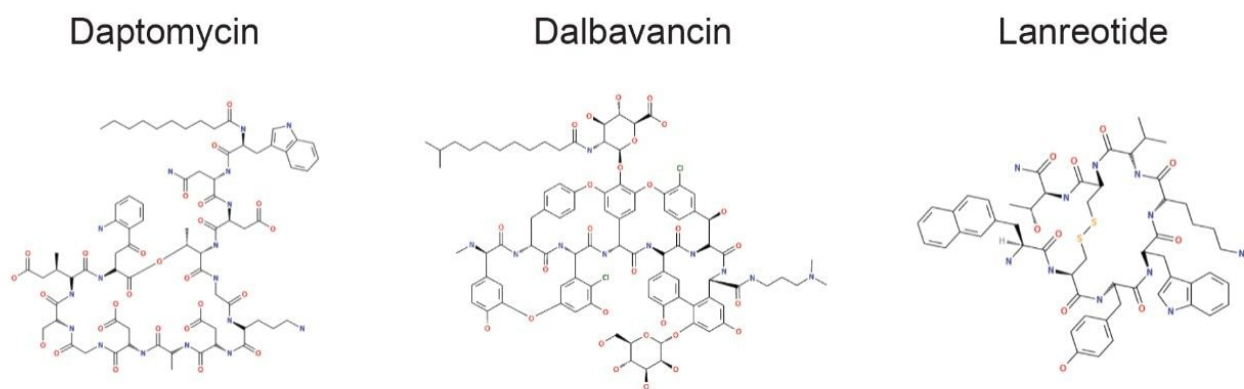

*Figure 1. Examples of semi-synthetic and/or natural product-derived therapeutic peptides.*

Modification	MMS/WMB compatibility	Selected examples
Simple linear (none)	Yes	GLP-1
non-natural/non-proteinogenic AA	Yes	Kynurenine, 3-methylglutamic acid (daptomycin)
Phenyl and chlorophenyl based units (dalbavancin)		
D-amino acids	Yes	Tryptophan (lanreotide)
Disulfide linked	Yes	Somatostatin, lanreotide
End-to-sidechain linked	Yes	Threonine-Kynurenine ester (daptomycin)
Non-natural linkages	Yes	Dalbavancin
		Decanoic acid (daptomycin)
		Napthalene (lanreotide)
Lipid modifiers/adducts	Yes	Methyl ester, glucopyranosyl, mannopyranosyl, dimethylaminopropyl, chlorophenyl based groups, glycolipid (dalbavancin)

Table 1. Peptide modification complexity.

## Turnover of Daptomycin, Dalbavancin, and Lanreotide

A small to moderate amount of catabolism was detected for all peptides. Daptomycin showed higher turnover after 300 min (>60%), while other peptides were found to be more stable and showed <10% turnover. The top three catabolites were reported for each peptide (Table 2). Daptomycin was chosen as a model compound to further demonstrate capabilities in DIA/DDA acquisition modes and software data processing.

Peptide	Turnover	Observed CCS (Å <sup>2</sup> )	Major catabolites		
Daptomycin	Med	406.1 +/- 1.6	+18 (hydrolysis)	-1130 (hydrolysis)	-986 (hydrolysis)
Dalbavancin	Low	456.7 +/- 1.8	-14 (demethylation)	-14 (demethylation)	-28 (2 x demethylation)
Lanreotide	Low	355.65 +/- 0.85	+18 (hydrolysis)	-197 (hydrolysis, loss of naphthalene amino group)	+18 (hydrolysis)

Table 2. Summary of turnover and major catabolites.

## Daptomycin catabolism

A summary of the major catabolites for daptomycin is reported in Figure 2, along with the summary trace at 300

min. Data were reviewed with filters applied that excluded singleton metabolites (those present only at one timepoint) and results that were  $> \pm 2$  mDa different from expected values. All other metabolites were reviewed for probable matches. Cluster plots (time courses) were reviewed to ensure profiles were consistent with turnover. Data were reviewed for plausible CCS values that corresponded to a given structure change. For example, the loss of an amino acid or a fatty acid tail was expected to result in a slightly smaller CCS at a given charge state, while ring opening was generally expected to decrease mobility and increase CCS. Consistency of mass accuracy, retention time, and CCS accuracy were also tracked by the software and used to help confirm metabolic profiles and further remove matrix false positives.



Figure 2. WebMetabase display of daptomycin incubation at 300 min. The top chromatogram is an overview of the substrate and major metabolites. The bottom table is a data summary for each observed peak.

The increase and/or disappearance of major species are displayed in the software as a cluster plot. An example of response versus time of major species for daptomycin is shown in Figure 3. WebMetabase automatically filters the data to show only metabolites that appear in more than one sample through the remove singletons option, further refining table results.



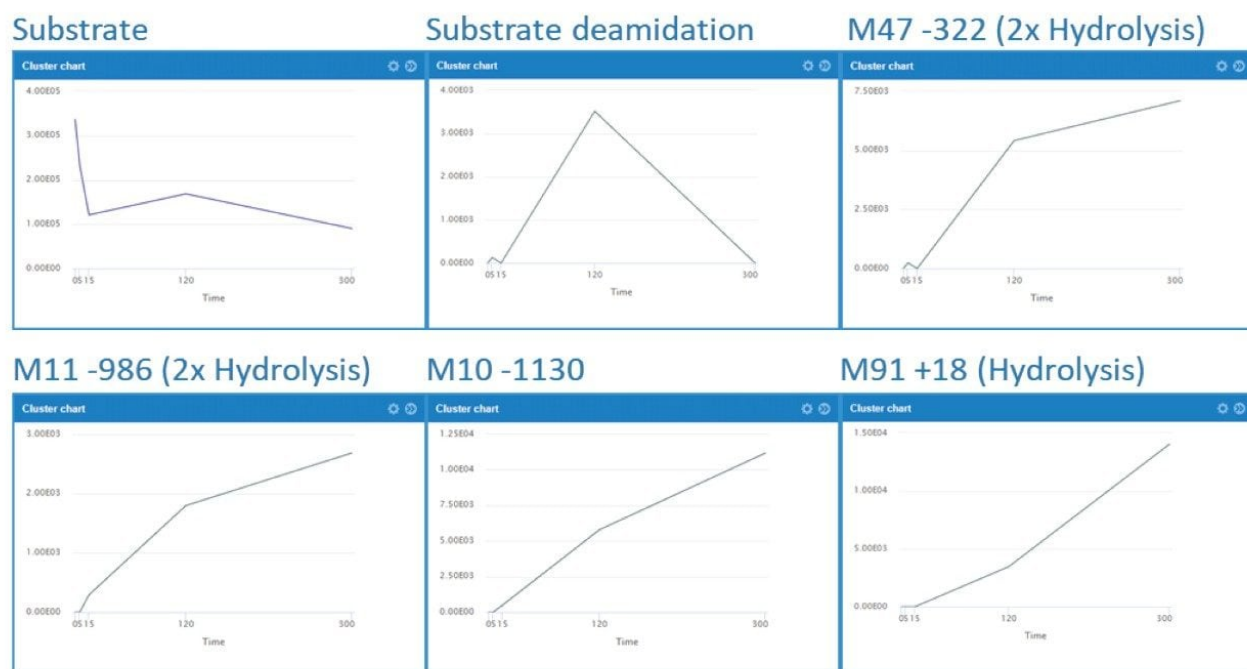


Figure 3. Daptomycin and major catabolic species. Cluster plots (time course) shown from 0–300 min.

Several peaks corresponding to amide hydrolysis were observed and putative assignments shown in Figure 4. For the major hydrolysis peak, M91, the site of biotransformation was localized to a cyclic cleavage product involving the hydrolysis of the ester, or a close by amide bond between the 3-anthraniloylalanine (kynurenine) and the 3-methylglutamic acid residues. Both sites have been previously reported as primary sites of metabolism.<sup>5</sup> Cleavage of the terminal fatty acid tail (M59-154) in isolation and in combination with the terminal tryptophan (M47-322) were also observed and reported by the software in these samples.

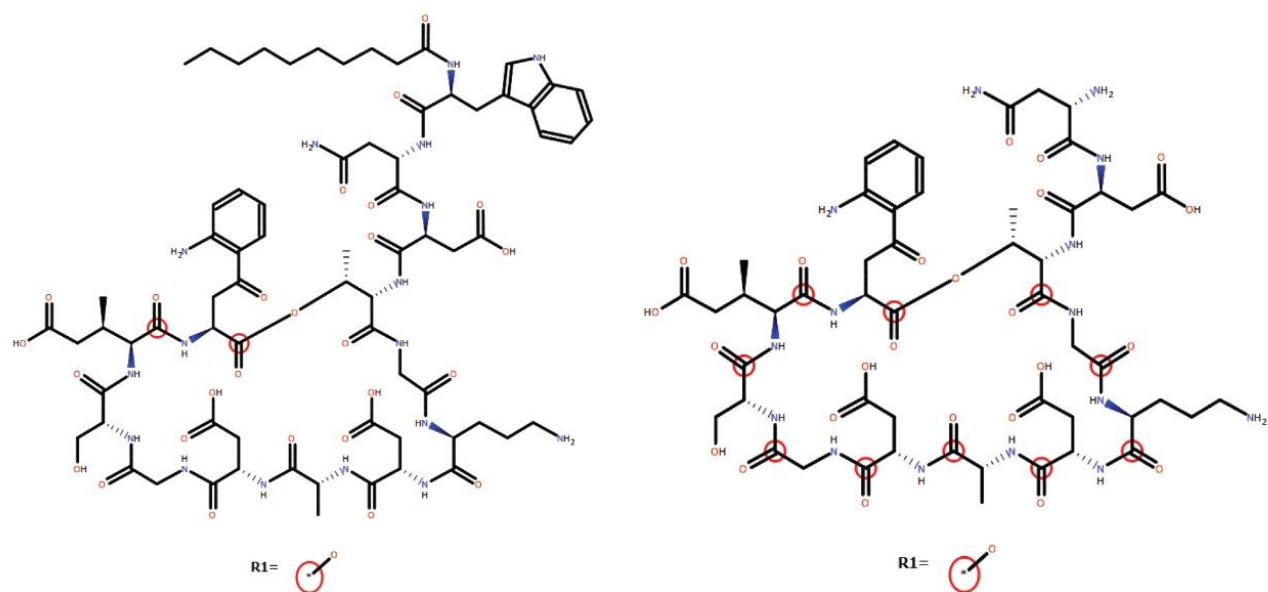


Figure 4. Examples of putative assigned structures in WebMetabase: Left/A) Major amide hydrolysis species, M91 and Right/B) loss of fatty acid/tryptophan tail section and ring hydrolysis, M47.

Experiments using QuanRecovery vials showed equivalent performance or better recovery for the API and metabolites. For this series of experiments, the incubation time was extended to 480 minutes. API recovery was found to be equivalent for polypropylene and QuanRecovery vials. For several metabolites such as M48 and M47, the recovery was consistently better using QuanRecovery vials with an average of 18% and 26% higher response, respectively.

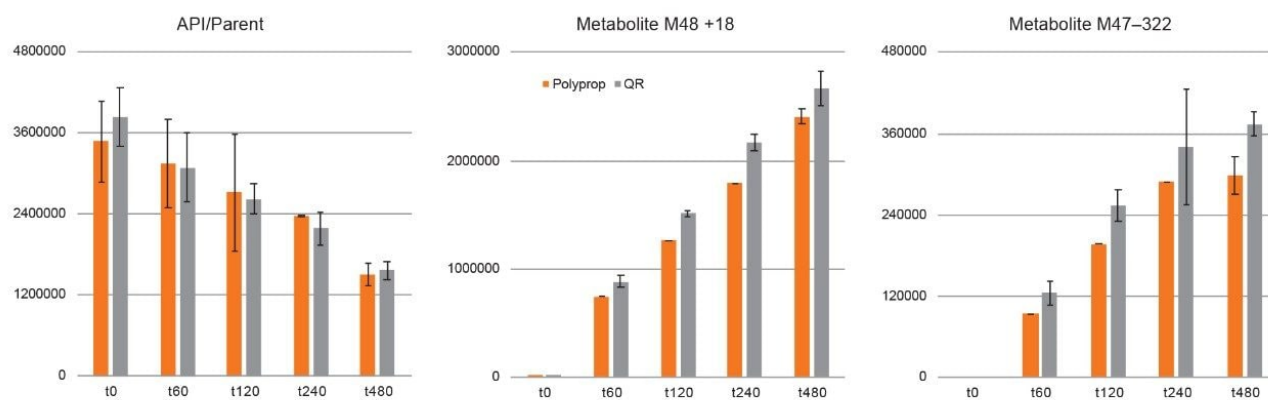


Figure 5. QuanRecovery performance for parent and hydrolysis products, M48 and M47.

## Using Product Ion Confirmation (HS-MSMS) With HDMSE

The addition of directed scanning functions to DIA methods enables full confirmation for ions of interest. Product ion confirmation scans (PICS) can be enabled for MS<sup>E</sup> or HDMS<sup>E</sup> analyses and spectra may be viewed in UNIFI. Figure 6 shows representative examples of data acquired without IMS processing, with IMS processing (HDMS<sup>E</sup>) and with HS-MSMS product ion spectra (ion mobility precursor selected MSMS) enabled. The data shows that IMS significantly enhances the spectrum by improving specificity and removing ions present in the unfiltered top trace, cleaning up noisy baselines and improving the quality of spectra. The use of drift time alignment can be used to combine multiple charge states into a single rich spectrum, containing all fragments produced by all major charge states. HS-MSMS data was optimized further (higher CE, mass dependent ramp) to generate a high-quality spectrum for the major (+2) charge state peaks. This results in a further refined spectrum with better peak quality, particularly in the mid region of the spectrum, which is more straightforward to interpret.

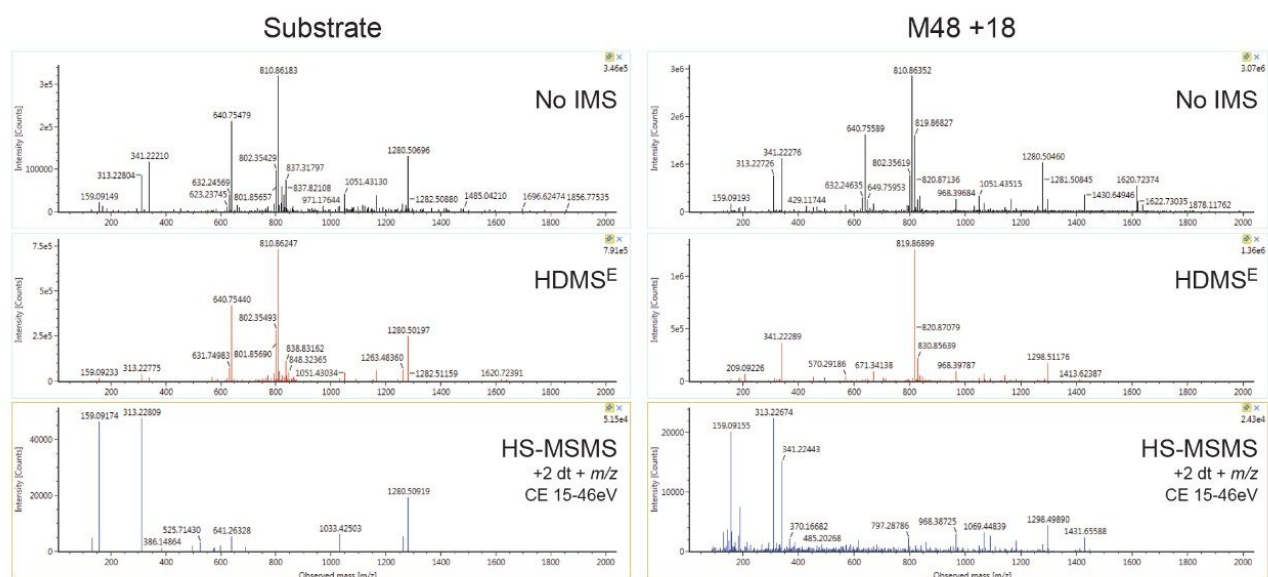


Figure 6. Substrate (left) and M48 (right) product ion scans using various acquisition modes: Top –  $MS^E$  (non IMS filtered DIA data), 15–70 eV ramp; middle –  $HDMS^E$  (IMS filtered DIA data), 15–70 eV ramp; bottom – targeted IMS PICS of  $m/z$  810, drift time 7.46 ms (Parent  $[M+2H]^{2+}$ ) and  $m/z$  819, drift time 7.67 ms (M48  $[M+2H]^{2+}$ ), 15–46 eV.

## Conclusion

Metabolites for 3 FDA-approved cyclic peptide drugs of varying structural complexity were identified and tracked using the WebMetabase macromolecule workflow. Ion mobility afforded additional resolution that was shown to help further discriminate, characterize, resolve, and track metabolites from interfering matrix ions having similar  $m/z$  but significantly different ion mobility. Ion mobility benefits both DIA and DDA approaches for generation of high-quality spectra. CCS enabled  $HDMS^E$  provided tools for tracking metabolites across multiple samples and instruments. Product ion confirmation scans (PICS) offer an easy way to build on the strength of the  $HDMS^E$  approach by providing a capability for adding true ion mobility and mass directed HS-MS/MS scans to DIA analysis.

The use of an ion mobility capable software package to process and review RT,  $m/z$ , MS-MS, and CCS information, along with cluster plots and integrated structural ID tools, help users quickly decide if peaks are real

and whether they warrant further investigation. Proper handling of peptide therapeutics is essential and the use of high-performance non-binding products such as QuanRecovery vials enables better and more consistent recovery of detected metabolites.

Combining ion-mobility enabled HDMS<sup>E</sup> data with a powerful software solution tailored for biotherapeutics enabled comprehensive metabolite identification of complex cyclic peptides needed for the development of efficacious biotherapeutics.

## Glossary - Ion mobility modes of acquisition and terminology

HDMS<sup>E</sup> – MS<sup>E</sup> is a full scan acquisition mode that uses alternating collision energies and aligns the precursor and fragment ion information for all species.<sup>6</sup> The low energy channel, is optimized for normal full scan transmission; and the high energy channel is optimized for precursor transmission. High Definition (HD) modes adds TWIMS (Travelling Wave Ion Mobility-based separation to both functions. Drift time information is captured by recording the drift time of all ions that pass through the ion mobility separation device and improves DIA data quality.

HS-MS/MS or HS-DDA MSMS – Once the drift time is known (determined from a routine HDMS scan mode), targeted MS/MS scans can select precursors exiting the ion mobility cell at a given time. HS (High Selectivity) scans further increase the specificity of precursor ions passed through for MS/MS using ion mobility. This type of scan is enabled when PICS is selected for an HDMSE analysis.<sup>7</sup>

CCS – Collision Cross Section – Represents the average cross section area for a molecule as it passes through a gas cell. Larger, extended structures and higher charge state species take longer to pass through a gas cell versus smaller more compact structures for the same given m/z creating an additional mechanism for ion separation.<sup>8</sup> For CCS on a Vion IMS QToF, drift times are calibrated against a known set of standards (using Major Mix, PN 186008113) during routine calibration. CCS values are calculated automatically for all analyses run using HD modes in UNIFI.

---

## References

1. Radchenko T, Brink A, Siegrist Y, Kochansky C, Bateman A, Fontaine F, Morettoni L, Zamora I. Software-Aided

Approach to Investigate Peptide Structure and Metabolic Susceptibility of Amide Bonds in Peptide Drugs Based on High Resolution Mass Spectrometry, 2017 PLoS ONE 12(11): e0186461.

2. Kirk J, Wrona M, Mortishire-Smith R, Zamora I, Fontaine F, Goshawk J, Lunt M, Doering K, Anderson N. Using Mass-MetaSite and WebMetabase to Process HDMSE Data Acquired on the Vion IMS QToF Mass Spectrometer. 2018. Waters Application Note. 720006362EN.
3. Sharma R, McDonald TS, Eng H, Limberakis C, Stevens BD, Patel S, Kalgutkar AS. In Vitro Metabolism of the Glucagon-Like Peptide-1 (GLP-1)-Derived Metabolites GLP-1(9–36) Amide and GLP-1(28-36) Amide in Mouse and Human Hepatocytes. Drug Metab Dispos, 2013 Dec; 41(12):2148–57.
4. Wrona MD, Kirk JM, Zamora I, Radchenko T, Escola A, Riera A, Mortishire-Smith RJ. Somatostatin Analogue Catabolite Screening and Identification Using Vion IMS Qtof with WebMetabase. 2019. Waters Application Note. 720006586EN.
5. d'Costa VM, Mukhtar TA, Patel T, Koteva K, Waglechner N, Hughes DW, Wright GD, De Pascale G. Inactivation of the Lipopeptide Antibiotic Daptomycin by Hydrolytic Mechanisms. Antimicrob Agents Chemother 2012 Feb; 56(2):757–64.
6. An Overview of the Principles of MSE. Waters White Paper 720004036EN.
7. Tomczyk N, Wallace A, Richardson K, Grzyb A, Wildgoose J. High Definition Data Directed Analysis. 2013. Waters Application Note. 720004737EN.
8. Pringle SD, Giles K, Wildgoose JL, Williams JP, Slade SE, Thalassinou K, Bateman RH, Bowers MT, and Scrivens JH. An Investigation of the Mobility Separation of Some Peptide and Protein Ions Using a New Hybrid Quadrupole/Travelling Wave IMS/oa-ToF instrument. Int J Mass Spectrom 2007 261(1): 1–12.

---

## Featured Products

[ACQUITY UPLC I-Class PLUS System <https://www.waters.com/134613317>](https://www.waters.com/134613317)

[Vion IMS QToF Ion Mobility Quadrupole Time-of-flight Mass Spectrometry <https://www.waters.com/134845751>](https://www.waters.com/134845751)

UNIFI Scientific Information System <<https://www.waters.com/134801648>>

720006850, April 2020



©2019 Waters Corporation. All Rights Reserved.

[Condiciones de uso](#) [Aviso de privacidad de Waters](#) [Marcas comerciales](#) [Empleo](#) [Avisos legales y de privacidad](#) [Cookies](#) [Preferencias de cookies](#)