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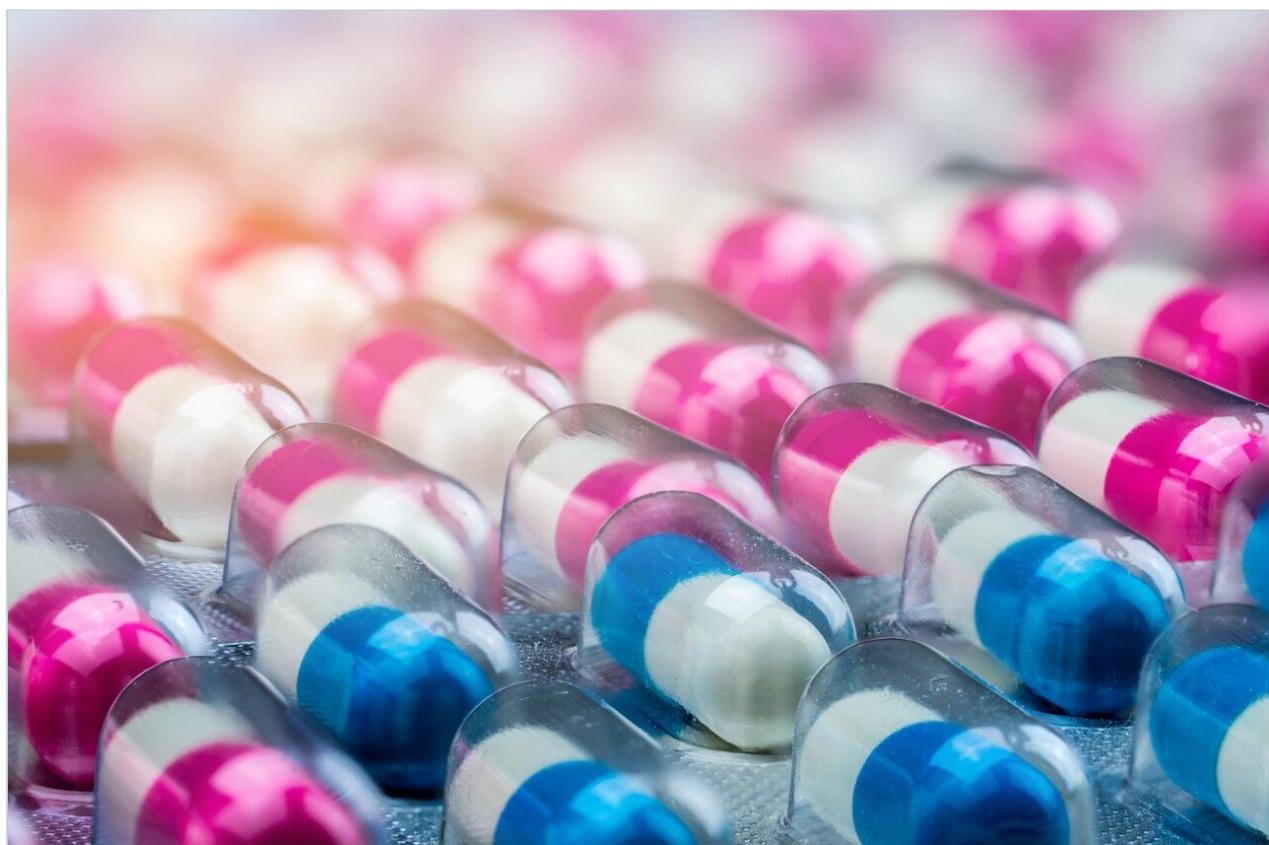
| Parameter | New HPLC method with MS compatible conditions | USP monograph method ⁶ |
|---|---|---|
| LC system | ACQUITY Arc | ACQUITY Arc |
| Detection | PDA (derived at 254 nm) and ACQUITY QDa | PDA (derived at 254 nm) |
| Column(s) | XSelect CSH C ₁₈ , 4.6 × 100 mm, 3.5 μm | Symmetry C ₁₈ , 4.6 × 250 mm, 5 μm |
| Column temp. | 40 °C | 30 °C |
| Injection volume | 10 μL | 20 μL |
| Flow rate | 1.5 mL/min | 1.0 mL/min |
| Mobile phase | Acetonitrile/10 mM ammonium formate (10/90) with 0.1% formic acid | Methanol/acetonitrile/water/phosphoric acid (100:100:800:2). Add 96 mg of sodium 1-pentanesulfonate in the resulting solution, filter |
| System suitability: standard and system suitability (SS) solution | <ul style="list-style-type: none"> Standard solution: 0.05 mg/mL of hydroxychloroquine (HCQ) in water SS solution: Chloroquine phosphate (CQ) and hydroxychloroquine sulfate (HCQ) at 0.05 mg/mL in water | <ul style="list-style-type: none"> Standard solution: 0.05 mg/mL of hydroxychloroquine (HCQ) in mobile phase SS solution: Chloroquine phosphate (CQ) and hydroxychloroquine sulfate (HCQ) at 0.05 mg/mL in mobile phase |

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| Parameter | USP requirement ⁶ | New HPLC method with MS compatible conditions | USP Method |
|--|------------------------------|---|---|
| Resolution between CQ and HCQ | Not less than (NLT) 1.8 | 3.7 | 2.5 |
| Relative standard deviation (RSD) for replicate injections of HCQ standard | Not more than (NMT) 1.5% | <ul style="list-style-type: none"> RSD of peak areas: 0.0% RSD of retention times: 0.1% | <ul style="list-style-type: none"> RSD of peak areas: 0.2% RSD of retention times: 0.2% |
| Peak tailing for CQ | N/A | 1.4 | 2.1 |
| Peak tailing for HCQ | N/A | 1.4 | 1.9 |

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1. Al-Bari, MA.Targeting endosomal acidification by chloroquine analogs as a promising strategy for the treatment of emerging viral diseases.*Pharmacol Res Perspect*.5(1):e00293 (2017).
2. Liu, J. et al.Hydroxychloroquine, a less toxic derivative of chloroquine, is effective in inhibiting SARS-CoV-2 infection in vitro.*Cell Discov*.6, 16 (2020).
3. Kupferschmidt, K. Big studies dim hopes for hydroxychloroquine.*Science*.368, 1166-1167 (2020).DOI: 10.1126/science.368.6496.1166
4. Who.int [Internet].World Health Organization; c2020 [cited 09 Jul 2020].Available from: <https://www.who.int/news-room/detail/04-07-2020-who-discontinues-hydroxychloroquine-and-lopinavir-ritonavir-treatment-arms-for-covid-19> <<https://www.who.int/news-room/detail/04-07-2020-who-discontinues-hydroxychloroquine-and-lopinavir-ritonavir-treatment-arms-for-covid-19>>
5. NIH.gov [Internet].Bethesda: U.S. Department of Health and Human Services; c2020.[cited 09 Jul 2020].Available from: <https://www.nih.gov/news-events/news-releases/nih-halts-clinical-trial-hydroxychloroquine> <<https://www.nih.gov/news-events/news-releases/nih-halts-clinical-trial-hydroxychloroquine>>
6. Tropmedres.ac [Internet].Thailand: Mahidol University; c2020 [cited 09 Jul 2020].Available from: <https://www.tropmedres.ac/covid-19/copcov> <<https://www.tropmedres.ac/covid-19/copcov>> .
7. USP Monograph, Hydroxychloroquine Sulfate Tablets, USP42-NF37, The United States Pharmacopeia Convention, official 23-March-2020.

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ACQUITY QDa 質量検出器 <<https://www.waters.com/134761404>>

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© 2021 Waters Corporation. All Rights Reserved.