# Waters<sup>™</sup>

#### 應用手冊

# Monitoring System Performance With System Suitability and Control Charts in Empower

Haining Pee

Waters Corporation

這是一篇應用簡報,不含詳細的實驗內容章節。

#### Abstract

System Suitability is a series of injections you can make to help you determine the suitability and effectiveness of your chromatographic system before use. System Suitability requirements must pass to guarantee the quality of the sample run results. System Suitability injections also enable you to monitor method performance over time.

### Introduction

Empower<sup>™</sup> Chromatographic Data System (CDS) can assess System Suitability standards interactively, checking the performance of standards against your System Suitability requirements in an automated, unattended manner. Upon failing the requirements, Empower CDS can be configured to continue the sample run, abort the sample run, or reinject the standards. Subsequently, System Suitability parameters can be plotted on control charts for data trending. Data trending allows users to monitor method performance and take corrective action before out-

of-specification products are produced.

The functionality of interactive System Suitability and control charts are highlighted through this case study below:

Six System Suitability standards are analyzed at the start of the sample run and, to ensure that the system is performing well, the analyte's Retention Time must fall within  $\pm 3\%$  of the target Retention Time.

Analyte	Target retention time (min)
Acesulfame K	0.650
Benzoic acid	1.004
Sorbic acid	1.450
Caffeine	1.700
Aspartame	2.000

*Table 1. Table of target Retention Time of the various analytes.* 

### **Results and Discussion**

#### In the Processing Method

First, System Suitability criteria are set up in your Empower Processing Method through the *Limits Tab*. Click on the target component and select the parameters that you wish to monitor in the Suitability Limits table. For this example, the target Retention Time and an error % of 3% for all components has been set.

Ac     Be     Soi     Cat     As     Ret	enzoic A orbic Aci- affeine spartame Field N etention	e K cid d e Vame Time	Flag Val. Na Target 0.650	me Error % 3.00	Lower E	rror Limit (Li			Suital	bility Co late Sui IV IV IV	emponent t Results	:S						Flag Out	tside Limits Iマ Iマ	\$	
Ac Be So Cat As Ref Weather Newton	enzoic A orbic Aci offeine spartame Field N etention	e K cid e Vame Time	Na Target 0.650	me Error % 3.00	Lower E	rror Limit (L			Suital	bility Co late Sui r r r	mponent t Results	S						Flag Out	tside Limits जि जि	<u>،</u>	
Ac Be So Cat As Ref Ref Verofilme	enzoic A orbic Aci- offeine spartame Field N etention	e K d a Vame Time	Na Target 0.650	Error % 3.00	Lower E	rror Limit (L			Calcul	late Sui	t Results							Flag Out	tside Limits 교 교 교		
Ac Be So Ca Asj Rel Rel	enzoic A orbic Aci affeine spartame Field N etention	e K d a Vame Time	Target 0.650	Error % 3.00	Lower E	rror Limit (L				র র র									হা হা		
Be So Ca As Ref Ref	enzoic A orbic Acia affeine spartame Field N etention	cid d e Vame Time	Target 0.650	Error % 3.00	Lower E	rror Limit (L				র র র									ম ম		
So Car As Rel Resulting	orbic Aci affeine spartame Field N etention	d e Vame Time	Target 0.650	Error % 3.00	Lower E	rror Limit (L				<u> </u>									v		
Cal As As Rel	affeine spartame Field N etention	e Name Time	Target 0.650	Error % 3.00	Lower E	rror Limit (L)				<b>K</b>									_		
As Rel	Field N	e Name Time	Target 0.650	Error % 3.00	Lower E	rror Limit (L				7					-				V		
Ret Ret	Field N	Name Time	Target 0.650	Error % 3.00	Lower E	rror Limit (L			~										<b>V</b>		
Re'	Field N	Name Time	Target 0.650	Error % 3.00	Lower E	rror Limit (L															
Rei megatan)	etention	Time	0.650	3.00		1-	CL) Uppe	r Error Limit	(UCL)	Warni	ng % L	ower W	Varni	ng Limit	Upp	er Warning L	imit	Ignore Blan	k Values	Check	Limits
Integration]						0.0	531		0.670											During Quar	ntitation
neepation) Neevallame Serial Acid Serial Acid Serial Acid																					
negation) Neevaltame Social Acid Social Acid Social Acid Social Acid																					
Serac Acie Serac Acie Cafeire Napatame	H Fag	Values Outside Law Rigense		S. Ce	tability Corroor tubility Corroor	ents As		Flag Cutside Linit	ti -	_	Anufarat	D RgV	Anni Dubid Iarine	a Lianka	-	Data Calcul	riky Compo da Duk Res	urda ulta	-	Flag Outside Linit	is
Soroc Acie Cafeire Nepatame	ekt				P			P		1	Ewizoic Acid									¥	
hapatane	•				8					4	2 Selscard 4 Caline			, i i i i i i i i i i i i i i i i i i i			-				
	(				P			P		-	Aspatiane									9	
		1			Sublicky Line	5		1				-	1			0.	tability Live				
Autorition 1	Time 1.00	04 3.00	0.074	1.3	4	Creative search Texa	chine arriving care	C	During Quart	Hation 1	Fadetian Tro	· 1.450	24	0	1407	1 434	waring is	Low Wanny Link	Caper Warring Link	Influence Billion Statuse	During Quantitation
Hegeter]	(troatingOth 19 Ray)	est   Dormonenty Value: Outville Link	treporty] Pare Ration :	MD too Famoro Dorteuro	and the first of the second	and brogs[Treed St	nigh) Outstatuty Linea	( horse and (off)			<u>).</u> megatan ji Sm	P Reg Va	Correan Ani Dunik	estallergunts ( Peer - Seite	(Rates) (M	ti on Rokos   Detaun Arte Sunat	ataituny (n Rey Compos	samed Groups) firmet Gro nems	ngn Suituolity Canta	Notive and Det()	
		Name		0	culate Suit Res	da		Flag Outside Line	is .	1	4	N	ane .			Cathyl	ee Sut Res	uts		Plag Dutside Link	h
Ackselfame Decesio Re-	w K		-		9			2		2	Acessifame K Bantaia Acid				-		-		-		
Sorbe Acid	6				8		-	8		3	Sotie Arid										
afere .					N.			ų		4	Caffeine						9				
Aspatane					9		_	P		5	Aquiane						P			9	
-					Subability Liev		1.8	1				1000			100	Se	tablity Line			10-1-12-12	
Field Na Retention 1		et Dror % S	we Dror Linit (LCL) 1.443	Upper Davar Limit (J.C 1.3	2) Warning Ni	Lower Warning Limit	Upper Warring Lini	I Sprore Dlank Yalves	Oving Quert	ints I	Field Nerve Retextion: Time	2000	Enu %	bree Enviro	4 (SCL) 1 540	Upper Dros Limit (UCL) 2 800	Warning %	Cover Warning Link	Upper Warning Linit	lignore Diank Volues	Check Links During Quantitation
	terre Targ	80 5.08			-																4

Figure 1. Click on each of the target component and select the parameters you wish to monitor in the Suitability Limits table.

#### In the Method Set

The Processing Method is added into the Method Set to allow Empower to process and evaluate your System Suitability criteria during the run.



Figure 2. Adding the Processing Method to the Method Set to allow for automated processing.

#### In the Run Samples Window

After your Sample Set is set up, the Run Mode and Suitability Mode are configured to enable interactive System Suitability. In the *Run Mode*, choose "Run Only", "Run and Process", or "Run and Report" for online data processing. "Run and Process" and "Run and Report" will result in the generation of results. Here, select "Run and Process".

	1	3	6	k	a 🖬 🍕 🐚	W III Ru	n Only 💌	Contin	ue on Fault	•
						Sample Set Method: Soi Bur	and Process	3		
E	Vial	Inj Vol (uL)	# of Injs	Label	SampleName	Function	Export Method	Label Reference	Processing	Run Time (Minutes)
1	1	5.0	1		Blank	Inject Samples	Soft Drink Analysis ACQ		Ignore Faults	3.00
2	2	5.0	6	SST	System Suitability Standard	Inject Samples	Soft Drink Analysis ACQ		Normal	3.00
3						Summarize Custom Fields			Normal	
4	3	5.0	1		Standard 1	Inject Standards	Soft Drink Analysis ACQ		Normal	3.00
5	4	5.0	1		Standard 2	Inject Standards	Soft Drink Analysis ACQ		Normal	3.00
6	5	5.0	1		Standard 3	Inject Standards	Soft Drink Analysis ACQ		Normal	3.00
7	6	5.0	1		Standard 4	Inject Standards	Soft Drink Analysis ACQ		Normal	3.00
8	7	5.0	1		Standard 5	Inject Standards	Soft Drink Analysis ACQ		Normal	3.00
9	8	5.0	1		Sample A	Inject Samples	Soft Drink Analysis ACQ		Ignore Faults	3.00
10	9	5.0	1		Sample B	Inject Samples	Soft Drink Analysis ACQ		Ignore Faults	3.00
11	10	5.0	1		Sample C	Inject Samples	Soft Drink Analysis ACQ		Ignore Faults	3.00

Figure 3. Select the Run Mode for online data processing.

The *Suitability Mode* allows you to configure Empower to take actions to continue the run, stop the run or reinject samples, in the event of System Suitability faults. Faults happen when the injection fails to meet the limits set. For this example, we will choose "Stop on Fault".

2	0	8	8	k	a 🖬 🚺 🐚	W III H Ru	n and Process 💌	Continue on Fault	]
						Sample Set Method: Soft Drin	k Analysis	Stop on Fault	-
Ē	Vial	Inj Vol (uL)	# of Injs	Label	SampleName	Function	Method Set / Report or Export Method	Reinject on Fault Re Next Sample on Fault	)
1	1	5.0	1		Blank	Inject Samples	Soft Drink Analysis ACQ	Ignore Faults 3.0	щ
2	2	5.0	6	SST	System Suitability Standard	Inject Samples	Soft Drink Analysis ACQ	Normal 3.0	00
3						Summarize Custom Fields		Normal	
4	3	5.0	1		Standard 1	Inject Standards	Soft Drink Analysis ACQ	Normal 3.0	00
5	4	5.0	1		Standard 2	Inject Standards	Soft Drink Analysis ACQ	Normal 3.0	00
6	5	5.0	1		Standard 3	Inject Standards	Soft Drink Analysis ACQ	Normal 3.0	00
7	6	5.0	1		Standard 4	Inject Standards	Soft Drink Analysis ACQ	Normal 3.0	00
8	7	5.0	1		Standard 5	Inject Standards	Soft Drink Analysis ACQ	Normal 3.0	00
9	8	5.0	1		Sample A	Inject Samples	Soft Drink Analysis ACQ	Ignore Faults 3.0	00
10	9	5.0	1		Sample B	Inject Samples	Soft Drink Analysis ACQ	Ignore Faults 3.0	00
11	10	5.0	1		Sample C	Inject Samples	Soft Drink Analysis ACQ	Ignore Faults 3.0	00

*Figure 4. Select the Suitability Mode to configure Empower to take specific actions in the event of a System Suitability failure.* 

Next, the System Suitability standards are analyzed. Empower continues your sample run if System Suitability limits pass or aborts the run if System Suitability limits fail.



Figure 5. Empower continues the sample run if System Suitability limits pass or aborts the run if

System Suitability limits fail. This allows for an unattended, hands-free operation.

Upon detecting a System Suitability fault in the second injection of the System Suitability standard, Empower aborts the run and indicates a suitability failure. You can review possible faults in the Review window.



Figure 6. Review window displaying System Suitability failure for Benzoic Acid.

System Suitability faults are automatically highlighted in red, bold, and italics for you. Here, it is observed that Benzoic Acid's Retention Time of 0.973 min falls outside the  $\pm 3\%$  error range of 0.974 to 1.034 min.

#### Data Trending in Publisher

Benzoic acid's Retention Time can be trended across multiple injections to determine if there is a drift towards an out-ofspecification result. To observe the data trending across multiple results, a Summary Report is required. This can be set up through Method Properties. Access Method Properties through the Edit Menu and select Method Properties. Here, we choose a Summary by All report.

Method Properties	×
Method Snap-To-Grid Legacy Report Type O Individual O Summary By All	Scaling Paper Size Vise Printer Defaults Size: 8.5x11
Margins Left: 0.50 Right: 0.50 Bottom: 0.50 Constant Constant C	Width: 8.50" Height: 11.00" Orientation Use Printer Defaults Portrait C Landscape
Scale to % of Paper Size: 100 Sign Off Allow this metho Note: This value will be be clear Suppress Errors Suppress data mismatch an	0.00 v d to be used in Sign Off red whenever a method is modified. d wrong chrom channel type errors
OK Cano	Apply Help

Figure 7. Method Properties in Report Publisher.



Figure 8. Report Publisher allows users to plot control charts via System Suitability groups.

On the right is the report method layout, while the list of report groups can be found on the left. System Suitability groups allow you to plot control charts. To observe the data trending for Retention Time, the SS Summary Plots are used. There are three display options for your Summary Plot, namely, Bar Plot, Line Plot, and Scatter Plot.

By default, Area will be trended for your Summary Plot. This can be changed in the Summary Plot Properties. Right click on the plot to access the Summary Plot Properties. There is an extensive selection of parameters – Empower built-in fields and custom fields inclusive – to choose from. We will select Retention Time to assess the trending of benzoic acid's Retention Time over six System Suitability standard injections.

Fonts		Colors			Name	
Sum Plot X-Axis Plot	Peaks	Scaling	Rotation	Legend	Order By	Bars/Points
Relative Response Relative RT Response Recention Time RF RF RT Ratio Slope Start Height Start Time TargetAmount Width PDA Peak Sys. Suit. Peak Sys. Suit. Peak Libraries Fraction	<ul> <li>Format</li> <li>Ba</li> <li>V</li> <li>I</li> <li>So</li> <li>Include</li> <li>I</li> <li< td=""><td>ar Plot he Plot catter Plot e in Plot bel Points w/L oper Error oper Warning ower Warning arget elative Limit 1 elative Limit 2 elative Target</td><td>Fit Parameter  Plot Fit  Fit Parameter  Plot Fit  Fit Parameter  Fit Parameter</td><td>rs r Fit ntial Notation ming ming</td><td></td><td></td></li<></ul>	ar Plot he Plot catter Plot e in Plot bel Points w/L oper Error oper Warning ower Warning arget elative Limit 1 elative Limit 2 elative Target	Fit Parameter  Plot Fit  Fit Parameter  Plot Fit  Fit Parameter  Fit Parameter	rs r Fit ntial Notation ming ming		
<		OK		ancel	Apply	Help

Figure 9. Summary Plot Properties allow users to set the parameter that they would like to trend.

The Summary Line Plot of benzoic acid's Retention Time over six injections revealed that the injections were mostly trending near the lower error limit. Investigations and preventive corrective actions would be necessary to ensure that the method continues to maintain its optimum performance.



Figure 10. Summary line plot of benzoic acid's Retention Time trending over six injections.

## Conclusion

Empower CDS allows for an assessment of your System Suitability standards in an automated manner. Laboratory productivity is increased as you do not have to manually assess System Suitability criteria. Upon failing the requirements, sample runs can be configured to continue, abort, or reinject the standards, allowing for a true hands-off system operation. Control charts provide you graphical representations of the result performance trends over time, allowing you to easily review and identify out-oftrend results.

Empower CDS's interactive System Suitability and control charts work in tandem to help assure that your laboratory continuously generates valid results while maintaining laboratory productivity.

# Featured Products

Empower Chromatography Data System (CDS) <https://www.waters.com/nextgen/global/products/informaticsand-software/chromatography-software/empower-software-solutions/empower-cds.html>

720007010, October 2020

 $\wedge$ 

© 2024 Waters Corporation. All Rights Reserved. 使用條款 隱私權政策 商標 就業機會 法律和隱私權聲明 Cookie Cookie偏好設定