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1. <http://www.pharmtech.com/enhancing-drug-development-applying-lc-ms-ms-cleaning-validation-manufacturing-equipment>. <<http://www.pharmtech.com/enhancing-drug-development-applying-lc-ms-ms-cleaning-validation-manufacturing-equipment>>
2. Sharma A, Vaghela J.S, Sharma P, Sharma B, Agrawal M, Jagiwala S. Cleaning Validation in Pharmaceutical Industry.A Comprehensive Approach; pharमतutor-art-1755.
3. Robert Fredrick.The Basic Facts of Cleaning Validation.Canada: Pharmainfo.net; 2004.Available from <http://www.pharmainfo.net/reviews/basic-facts-cleaning-validation>.
4. LeBlanc DA.Establishing Scientifically Justified Acceptance Criteria for Cleaning Validation of Finished Drug Products.Pharm Tech.1998; 22(10): 136–148.
5. Fourman GL, Mullen MV.Determining Cleaning Validation Acceptance Limits for Pharmaceutical Manufacturing Operations.Pharm Tech. 1993; 17: 54–60.
6. M.B. Boca, Z. Apostolides, E. Pretorius.A Validated HPLC Method for Determining Residues of a Dual Active Ingredient Anti-Malarial Drug on Manufacturing Equipment Surfaces.J. Pharm.Biomed.Anal. 37 (3), 461–468 (2005).
7. M.J. Nozal, et al. Development and Validation of an LC Assay for Sumatriptan Succinate Residues on Surfaces in the Manufacture of Pharmaceuticals.J. Pharm.Biomed.Anal. 30 (2), 285-291 (2002).
8. J. Lambropoulos, G.A. Spanos, N.V. Lazaridis.Development and Validation of an HPLC Assay for Fentanyl, Alfentanil, and Sufentanil in Swab Samples.J. Pharm.Biomed.Anal. 23 (2-3), 421–428 (2000).
9. T. Mirza, et al. Cleaning Level Acceptance Criteria and a HighPressure Liquid Chromatography Procedure for the Assay of Meclizine Hydrochloride Residue in Swabs Collected from Pharmaceutical Manufacturing Equipment Surfaces. J. Pharm.Biomed Anal. 19 (5), 747-756 (1999).
10. D.A. LeBlanc.Establishing Scientifically Justified Acceptance Criteria for Cleaning Validation of APIs.Pharm Technol. 24 (10), 160-168 (2000).
11. D.A. LeBlanc.Establishing Scientifically Justified Acceptance Criteria for Cleaning Validation of Finished Drug Products.Pharm Technol. 22 (10), 136-148 (1998).

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