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---

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1. Tran K, Young M, Organtini K, Twohig M, Hudalla C. Determination of Pesticide and Mycotoxin Residues in Dried Cannabis Flower: LC-MS/MS and GC-MS/MS Methodology to Meet the Recommended AOAC Regulatory Requirements for US States and Canada. Waters Application Note [720006780EN](https://www.waters.com/webassets/cms/library/docs/720006780en.pdf) <<https://www.waters.com/webassets/cms/library/docs/720006780en.pdf>> 2020 Feb.
2. Canellas E, Vera P, Domeño C, Alfaro P, Nerín C. Atmospheric Pressure Gas Chromatography Coupled to Quadrupole-Time of Flight Mass Spectrometry as a Powerful Tool for Identification of Non-Intentionally Added Substances in Acrylic Adhesives used in Food Packaging Materials. Journal of Chromatography A .2012 Apr 27;1235:141–8.
3. Bureau of Cannabis Control. California Code of Regulations Title 16 Division 42. Adopted June 2018. [https://bcc.ca.gov/law\\_regs/readopt\\_text\\_final.pdf](https://bcc.ca.gov/law_regs/readopt_text_final.pdf) <[https://bcc.ca.gov/law\\_regs/readopt\\_text\\_final.pdf](https://bcc.ca.gov/law_regs/readopt_text_final.pdf)> .
4. United States Pharmacopeia and National Formulary. USP <467> Residual Solvents. Official December 1, 2020. [https://www.uspnf.com/sites/default/files/usp\\_pdf/EN/USPNF/revisions/gc-467-residual-solvents-ira-20190927.pdf](https://www.uspnf.com/sites/default/files/usp_pdf/EN/USPNF/revisions/gc-467-residual-solvents-ira-20190927.pdf) <[https://www.uspnf.com/sites/default/files/usp\\_pdf/EN/USPNF/revisions/gc-467-residual-solvents-ira-20190927.pdf](https://www.uspnf.com/sites/default/files/usp_pdf/EN/USPNF/revisions/gc-467-residual-solvents-ira-20190927.pdf)> .
5. European Medicines Agency. ICH Guideline Q3C (R6) on Impurities: Guideline for Residual Solvents. August 9, 2019. [https://www.ema.europa.eu/en/documents/scientific-guideline/international-conference-harmonisation-technical-requirements-registration-pharmaceuticals-human-use\\_en-33.pdf](https://www.ema.europa.eu/en/documents/scientific-guideline/international-conference-harmonisation-technical-requirements-registration-pharmaceuticals-human-use_en-33.pdf) <[https://www.ema.europa.eu/en/documents/scientific-guideline/international-conference-harmonisation-technical-requirements-registration-pharmaceuticals-human-use\\_en-33.pdf](https://www.ema.europa.eu/en/documents/scientific-guideline/international-conference-harmonisation-technical-requirements-registration-pharmaceuticals-human-use_en-33.pdf)> .

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