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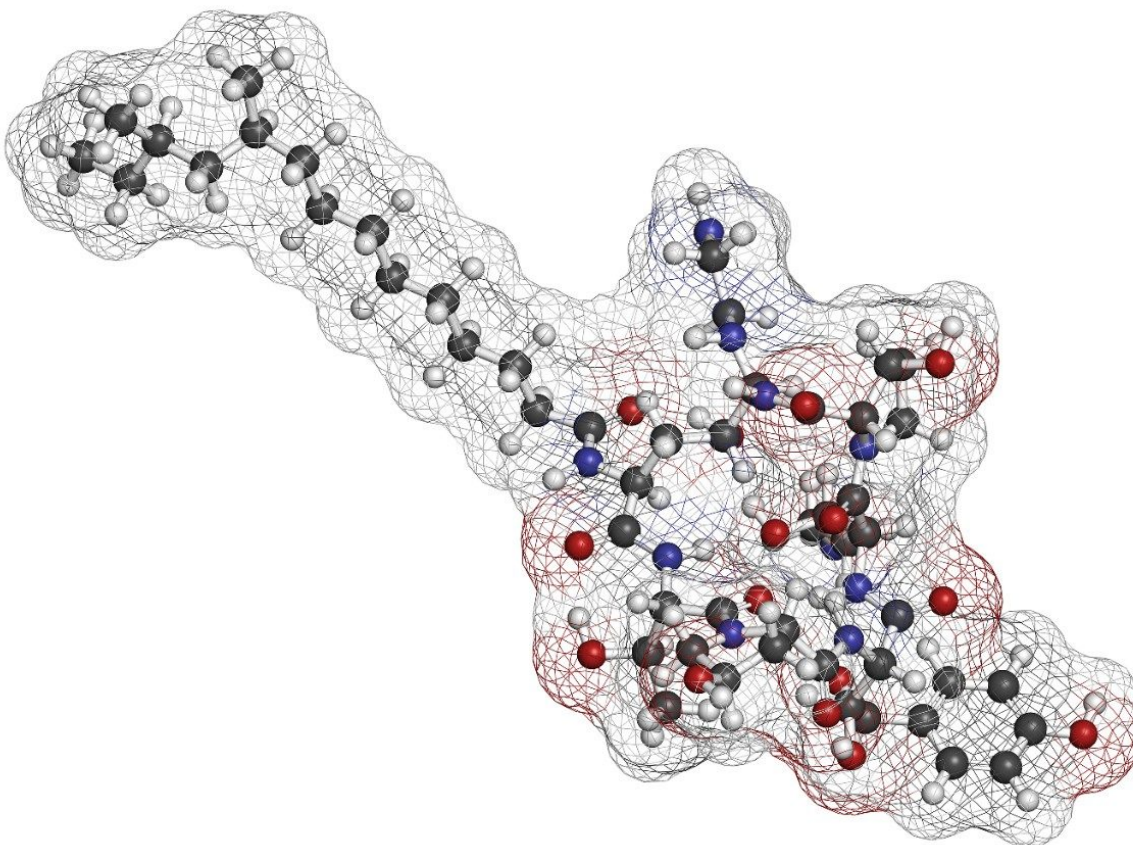
1. Higel F, *et al.* N-glycans of Complex Glycosylated Biopharmaceuticals and their Impact on Protein Clearance. *European Journal of Pharmaceutics and Biopharmaceutics*.2019; 139, 123–131.
2. Zhang P, *et al.* Challenges of Glycosylation Analysis and Control: An Integrated Approach to Producing Optimal and Consistent Therapeutic Drugs.*Drug Discovery Today*. 2016; 21(5), 740–765.
3. Yu, Y. Applying a Novel Glycan Tagging Reagent, *RapiFluor*-MS, and An Integrated UPLC-FLR/QToF MS System for Low Abundant N-Glycan Analysis.Waters Application Note, 2015.[720005383EN < https://www.waters.com/nextgen/us/en/library/application-notes/2015/applying-glycan-tagging-reagent-rapi-fluor-ms-and-integrated-uplc-flr-qtof-ms-system-for-low-abundant-n-glycan-analysis.html> .](https://www.waters.com/nextgen/us/en/library/application-notes/2015/applying-glycan-tagging-reagent-rapi-fluor-ms-and-integrated-uplc-flr-qtof-ms-system-for-low-abundant-n-glycan-analysis.html)
4. Zhang X, Reed C, Shion H, Alley W, Birdsall R, and Yu Y. Increasing Productivity and Confidence for N-linked Glycan Analysis of Biosimilars Using the BioAccord System.Waters Application Note, 2019.[720006545EN < https://www.waters.com/nextgen/us/en/library/application-notes/2019/increasing-productivity-and-confidence-for-n-linked-glycan-analysis-of-biosimilars-using-the-bioaccord-system.html> .](https://www.waters.com/nextgen/us/en/library/application-notes/2019/increasing-productivity-and-confidence-for-n-linked-glycan-analysis-of-biosimilars-using-the-bioaccord-system.html)
5. DeLoff M. A Comparison of Premier Columns with MaxPeak HPS Technology Versus PEEK-lined Column Hardware.Waters Application Note, 2021.[720007210EN < https://www.waters.com/nextgen/us/en/library/application-notes/2021/a-comparison-of-premier-columns-with-maxpeak-hps-technology-versus-peek-lined-column-hardware.html> .](https://www.waters.com/nextgen/us/en/library/application-notes/2021/a-comparison-of-premier-columns-with-maxpeak-hps-technology-versus-peek-lined-column-hardware.html)
6. Isaac G. and Plumb R. ACQUITY Premier LC Technology Significantly Improves Sensitivity, Peak Shape, and Recovery for Phosphorylated and Carboxylate Lipids.Waters Application Note, 2021.[720007092EN < https://www.waters.com/nextgen/us/en/library/application-notes/2021/acquity-premier-lc-technology-significantly-improves-sensitivity,-peak-shape-and-recovery-for-phosphorylated-and-carboxylate-lipids.html> .](https://www.waters.com/nextgen/us/en/library/application-notes/2021/acquity-premier-lc-technology-significantly-improves-sensitivity,-peak-shape-and-recovery-for-phosphorylated-and-carboxylate-lipids.html)
7. Tanna N. and Plumb R. Advantages of using ACQUITY Premier System and Columns with MaxPeak HPS

Technology for Bioanalysis of Gefitinib – an EGFR Inhibitor. Waters Application Note, 2021. [720007122EN <https://www.waters.com/nextgen/us/en/library/application-notes/2021/advantages-of-using-acquity-premier-system-and-columns-with-maxpeak-hps-technology-for-bioanalysis-of-gefitinib-an-egfr-inhibitor.html>](https://www.waters.com/nextgen/us/en/library/application-notes/2021/advantages-of-using-acquity-premier-system-and-columns-with-maxpeak-hps-technology-for-bioanalysis-of-gefitinib-an-egfr-inhibitor.html) .

8. Liu X. and Lauber M. High Performance Surface Technology Improves HILIC Profiling of Released N-Glycans. Waters Application Note, 2021.
9. Kang J.F., *et al.* Lysosomal Targeting Enhancement by Conjugation of Glycopeptides Containing Mannose-6-phosphate Glycans Derived from Glyco-engineered Yeast. *Scientific Reports*. 2018; 8, 8730.
10. Yu Y. N-linked Glycan Characterization and Profiling: Combining the Power of Accurate Mass, Reference Glucose Units, and UNIFI Software for Confident Glycan Assignments. Waters Application Note, 2016. [720004845EN <https://www.waters.com/nextgen/us/en/library/application-notes/2016/n-linked-glycan-accurate-mass-reference-glucose-units-glycan-assignments0.html>](https://www.waters.com/nextgen/us/en/library/application-notes/2016/n-linked-glycan-accurate-mass-reference-glucose-units-glycan-assignments0.html) .

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© 2025 Waters Corporation. All Rights Reserved.

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