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Discovery Elemental Composition ChemSpider Fragment Match

Results (10 found) ^

Component Name	Elemental Composition	i-FIT Confidence (%)	Common Name	Fragment Matches	Citations	Predic
1 Candidate Mass 284.2709...	C16H33N3O	100.00	MFCD02671276	12	7	

Assign

Information ^

MFCD02671276

Synonyms

- Decanamide, N-[2-(1-piperaziny)ethyl]
- N-[2-(1-Piperaziny)ethyl]decanamide
- N-[2-(1-Piperaziny)ethyl]decanamid

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1. USP-NF/PF, Assessment of Drug Product Leachables Associated with Pharmaceutical Packaging/Delivery Systems.1664. https://doi.usp.org/USPNF/USPNF_M7127_03_01.html <
https://doi.usp.org/USPNF/USPNF_M7127_03_01.html>
2. USP-NF/PF, Assessment of Extractables Associated with Pharmaceutical Packaging/Delivery Systems.1663. https://doi.usp.org/USPNF/PNF_M7126_03_01.html <https://doi.usp.org/USPNF/PNF_M7126_03_01.html>
3. Norwood D., Paskiet D., Ruberto M., Feinberg T., Schroeder A., Poochikian G., Wang Q., Deng T., Degrazio F., Munos M., Nagao L. Best Practices for Extractables and Leachables in Orally Inhaled and Nasal Drug Products: An Overview of the PQRI Recommendations.*Pharmaceutical research*.25.727–39, 2008.
4. NIST/EPA/NIH Mass Spectral Library Compound Scoring: Match Factor, Reverse Match Factor, and Probability.Jordi Labs.
5. Stevens D. M., Cabovska B., Bailey A. E. Detection and Identification of Extractable Compounds from Polymers.Waters Application Note [720004211](#).January 2012.
6. Sanig R., Kirk J., Cojocariu C. Screening for Extractables and Leachables in Nasal Spray Devices Using Liquid Chromatography and High-Resolution Mass Spectrometry With a Data Independent Informatics Strategy.Waters Application Note [720008000](#), August 2023.

7. Atmospheric Pressure GC (APGC).Waters White Paper, 720004771, August 2013.
8. Reid L., Pickles D. Improved Transmission of Labile Species on the Xevo G3 QTof Mass Spectrometer with the StepWave™ XS.Waters Application Note [720007794](#), November 2022.
9. McCullagh M., Mortishire-Smith R. J., Goshawk J., Cooper J., Obkircher M., Sprecher H., Koehling R., Nold m., Jacobsen J., Sanig R., Extractables, Leachables, and Contact Materials: The Invaluable Benefit of Ion Mobility-Enhanced Mass Spectrometry Libraries.Waters Application Note [720007655](#), June 2022.
10. ISO 10993–18:2020 Biological Evaluation of Medical Devices—Part 18: Chemical Characterization of Medical Device Materials Within a Risk management process, <https://www.iso.org/standard/64750.html> <
<https://www.iso.org/standard/64750.html>>
11. Cabovska B. Screening Workflow for Extractable Testing Using the UNIFI Scientific Information System, Technical Note [720005688](#), April 2016.

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© 2025 Waters Corporation. All Rights Reserved.

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