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1. National Pesticide Information Center, DDT General Fact Sheet, accessed 14 February 2024, [LINK <http://npic.orst.edu/factsheets/ddtgen.pdf>](http://npic.orst.edu/factsheets/ddtgen.pdf) .
2. United Nations Environment Programme, The 12 initial POPs under the Stockholm Convention, accessed 14 February 2024, [LINK <http://chm.pops.int/TheConvention/ThePOPs/The12InitialPOPs/tabid/296/Default.aspx>](http://chm.pops.int/TheConvention/ThePOPs/The12InitialPOPs/tabid/296/Default.aspx) .
3. Van Bavel, Bert, *et al.* Atmospheric-Pressure Chemical Ionization Tandem Mass Spectrometry (APGC/MS/MS) an Alternative to High-Resolution Mass Spectrometry (HRGC/HRMS) for the Determination of Dioxins. *Analytical Chemistry* 87.17: 9047–9053, 2015.
4. De Wit, Cynthia A., Katrin Vorkamp, and Derek Muir. Influence of Climate Change on Persistent Organic Pollutants and Chemicals of Emerging Concern in the Arctic: State of Knowledge and Recommendations for Future Research. *Environmental Science: Processes & Impacts* 24.10: 1530–1543, 2022.
5. Karasek, F. W., and F. I. Onuska. Trace Analysis of the Dioxins. *Analytical Chemistry* 54.2: 309–324, 1982.
6. Yost, Richard A. The Triple Quadrupole: Innovation, Serendipity and Persistence. *Journal of Mass Spectrometry and Advances in the Clinical lab* 24: 90, 2022.
7. Commission Regulation (EU) No.589/2014 Laying Down Methods of Sampling and Analysis for the Control of Levels of Dioxins, Dioxin-Like PCBs and Non-Dioxin-Like PCBs in Certain Foodstuffs and Repealing Regulation (EU) No.252/2012, [LINK <https://www.fao.org/faolex/results/details/en/c/LEX-FAOC134439/>](https://www.fao.org/faolex/results/details/en/c/LEX-FAOC134439/) .

8. Converting Semivolatile GC-MS/MS Methods from Helium to Nitrogen Carrier Gas with APGC, an Atmospheric Pressure Ionization Source, March 2023, [LINK <https://www.researchgate.net/publication/373921828_Converting_Semivolatile_GC-MSMS_Methods_from_Helium_to_Nitrogen_Carrier_Gas_with_APGC_an_Atmospheric_Pressure_Ionization_Source>](https://www.researchgate.net/publication/373921828_Converting_Semivolatile_GC-MSMS_Methods_from_Helium_to_Nitrogen_Carrier_Gas_with_APGC_an_Atmospheric_Pressure_Ionization_Source) .
9. Smalling, Kelly L., and Kathryn M. Kuivila. Multi-Residue Method for the Analysis of 85 Current-Use and Legacy Pesticides in Bed and Suspended Sediments. *Journal of Chromatography A* 1210.1: 8–18, 2008.

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© 2024 Waters Corporation. All Rights Reserved.

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