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1. Hansen K J, Clemen L A, Ellefson M E., Johnson H O. Compound-Specific Quantitative Characterisation of Organic Fluorochemicals in Biological Matrices.*Environ.Sci.Technol.* 35, 766–770, 2001.
2. Giesy J P and Kannan K Global Distribution of Perfluorooctane Sulfonate in Wildlife.*Environ.Sci.Technol.* 35: 1339–1342, 2001.
3. Kelly B C, Ikonomou M G, Blair J D, Surridge B, Hoover D, Grace R, Gobas F A P C. Perfluoroalkyl Contaminants in an Arctic Marine Food Web: Trophic Magnification and Wildlife Exposure.*Environ.Sci.Technol.* 43 (11), 4037–4043, 2009.
4. EFSA report.Results of the Monitoring of Perfluoroalkylated Substances in Food in the Period 2000–2009.February 2011.
5. United States Environmental Protection Agency.Analysis of Per and Polyfluoroalkyl Substances (PFAS) in Aqueous, Solid, Biosolids, and Tissue Samples by LC-MS/MS.4th Draft method 1633, July 2023.
6. Organtini K, Rosnack K, Hancock P. Analysis of Per and Polyfluoroalkyl Substances (PFAS) in Accordance with EPA 1633 Part 1: Establishing and Assessing the Method.Waters Application note [720008117](#), 2023.
7. Khoury-Hollins H., Reid L., Adams S. Direct Injection Screening Method for Per and Polyfluoroalkyl Substances (PFAS) in Drinking Water Using the High Resolution Time of Flight Mass Spectrometer, the Xevo™ G3 QTof.Waters Application note.[720008198](#), 2024.
8. Commission Recommendation (EU) 2022/1431, on the Monitoring of Perfluoroalkyl Substances in Food.L 221/105, 26.8.2022.
9. Travis D Evidence That There are Two Forms of Fluoride in Human Serum.*Nature* 214: 1050–1051, 1968.
10. Lau C, Perfluorinated Compounds: An Overview.In: DeWitt, J.C. (Ed.), Toxicological Effects of Perfluoroalkyl and Polyfluoroalkyl Substances.Springer, pp.1–21, 2015.
11. Dickman R A, Aga D S. A Review of Recent Studies on Toxicity, Sequestration, and Degradation of Per- and Polyfluoroalkyl Substances (PFAS) *Journal of Hazardous Material.* 436: 129120, 2022.
12. Trudel D, Horowitz L, Wormuth M, Scheringer M, Cousins IT, Hungerbühler K. Estimating Consumer Exposure

to PFOS and PFOA. *Risk Anal.* 28:251–69, 2008.

13. De Silva A O, Armitage J M, Bruton T A, Dassuncao C, Heiger-Bernays W, Hu X C, Kärrman A, Kelly B, Ng C, Robuck A, Sun M, Webster T F, Sunderland E M. PFAS Exposure Pathways for Humans and Wildlife: A Synthesis of Current Knowledge and Key Gaps in Understanding Environ *Toxicol Chem.* 40: 631–657, 2021.
14. Schrenk D, Bignami M, et al, EFSA Panel on Contaminants in the Food Chain (EFSA CONTAM Panel), Risk to Human Health Related to the Presence of Perfluoroalkyl Substances in Food. *EFS2*;18(9) 2020.
15. EURL for Halogenated POPs in Feed and Food (2022): Guidance Document on Analytical Parameters for the Determination of Per- and Polyfluoroalkyl Substances (PFAS) in Food and Feed, version 1.2 of 11 May 2022. Available online under (https://eurl-pops.eu/core-working-groups#_pfas <https://eurl-pops.eu/core-working-groups#_pfas>)
16. Adams S, Dreolin N, Organtini K L, Hancock P. Analysis of Per- and Polyfluoroalkyl Substances (PFAS) in Animal Products with an Enhanced Sensitivity LC-MS/MS Method using Fish Reference Materials as a Case Study. *Waters Application note*. [720008108](#), 2023.
17. Twohig M, Fujimoto G, Mohan A, Organtini K L, Rosnack K J, Hird S. Approaches to Non-targeted Analyses of Per- and Polyfluoroalkyl Substances (PFAS) in Environmental Samples. *Waters Application note*. [720007184](#), 2021.
18. U.S. Environmental Protection Agency. CompTox Chemicals Dashboard. <https://comptox.epa.gov/dashboard/chemical-lists/pfasmaster> <<https://comptox.epa.gov/dashboard/chemical-lists/pfasmaster>> accessed November 2023.
19. Suspect List of Possible Per- and Polyfluoroalkyl Substances (PFAS). <https://data.nist.gov/od/id/mds2-2387> <<https://data.nist.gov/od/id/mds2-2387>> Accessed March 2023.
20. Bugsel B, Zweigle J, Zwiener C. Nontarget Screening Strategies for PFAS Prioritization and Identification by High Resolution Mass Spectrometry: A Review. *Trends in Environmental Analytical Chemistry.* 40: e00216, 2023.
21. Charbonnet J A, McDonough C A, Xiao F, Schwichtenberg T, Cao D, Kaserzon S, Thomas K V, Dewapriya P, Place B J, Schymanski E L, Field J A, Helbling D E, Higgins C P. Communicating Confidence of Per- and Polyfluoroalkyl Substance Identification via High-Resolution Mass Spectrometry. *Environ Sci Technol Lett.* 9: 473–481, 2022.

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© 2025 Waters Corporation. All Rights Reserved.

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