

Fast and Simple Bioanalytical Plasma Sample Extraction Using Oasis® PRiME HLB SPE for High Analyte Recovery

Mary Trudeau, Margot Lee

Waters Corporation

Ce document est une note d'application et ne contient pas de section détaillée concernant l'expérimentation.

Abstract

This application describes a fast and simple method for bioanalytical sample preparation from plasma. Utilizing Oasis PRiME HLB solid-phase extraction (SPE), the protocol requires no method development and achieves high recovery and low matrix effects for several small molecule pharmaceuticals.

Experimental

Oasis PRiME HLB SPE Protocol
96-well plate, 3 mg sorbent per well,
P/N 186008052



Sample pretreatment:

200 µL 4% H₃PO₄ to 200 µL plasma sample

Load:

400 µL pre-treated sample

Wash:

200 µL 95:5 Water: Methanol

Elute:

2x 25 µL 90:10 Acetonitrile/Methanol

Dilute:

50 µL H₂O

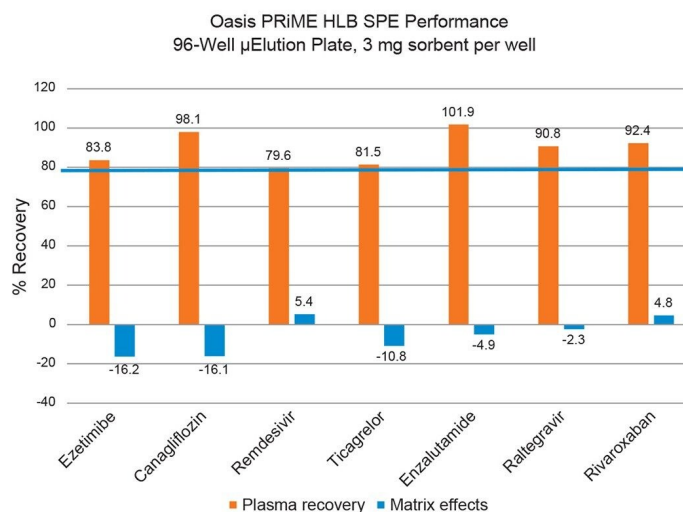


Figure 1. Demonstration of Oasis PRiME HLB SPE bioanalytical performance, requiring no optimization of the starting protocol, with high recovery (> 79%) and low matrix effects (-16,2 to 5.4%) for several pharmaceuticals extracted from plasma.

Results and Discussion

High Analyte Recovery with low matrix from plasma is achieved Using Oasis PRiME HLB SPE for top-selling small molecule pharmaceuticals in 2023.¹

References

1. Williams, RE and Leatherwood HM [Top 200 Small Molecule Drugs by Retail Sales in 2023](https://bpb-us-e2.wpmucdn.com/sites.arizona.edu/dist/9/130/files/2024/05/2023Top200SmallMoleculePosterV5.pdf) <<https://bpb-us-e2.wpmucdn.com/sites.arizona.edu/dist/9/130/files/2024/05/2023Top200SmallMoleculePosterV5.pdf>> .
(accessed 21 June 2024).

Featured Products

720008423, July 2024



© 2024 Waters Corporation. All Rights Reserved.

[Conditions d'utilisation](#) [Politique de confidentialité](#) [Marques](#) [Carrières](#) [Mentions légales et déclaration de confidentialité](#) [Cookies](#) [Préférences de cookies](#)