

Simple and Selective Bioanalytical Plasma Sample Extraction Using Oasis® MCX SPE for High Analyte Recovery

Mary Trudeau, Margot Lee

Waters Corporation

这是一份应用简报，不包含详细的实验部分。

Abstract

This application demonstrates the simple and selective bioanalytical plasma sample extraction using mixed-mode ion-exchange solid phase extraction (SPE), requiring no protocol method development, for high extraction MCX SPE recovery of several tyrosine kinase inhibitors (TKI) used in the treatment of leukemia.

Experimental

Oasis MCX SPE Protocol
96-well plate, μ Elution Plate, 2 mg
sorbent per well, 30 μ m
P/N 186001830BA



Sample Pretreatment:
200 μ L 4% H_3PO_4 to 200 μ L plasma sample
Load:
400 μ L pre-treated sample
Wash:
200 μ L 2% Formic acid in H_2O
Elute 1 (Neutral):
2x 25 μ L 100% MeOH
Elute 2 (M-M Fraction):
2x 25 μ L 5% NH_4OH in 100% MeOH
Dilute:
50 μ L 100% H_2O

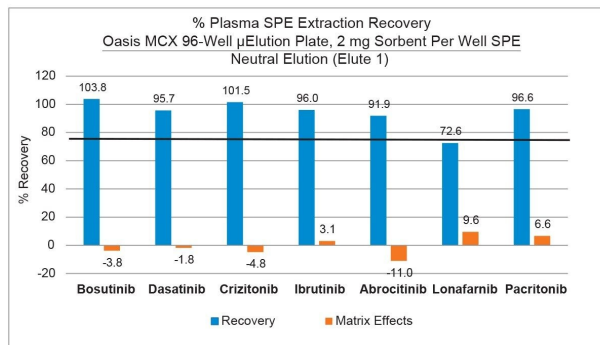


Figure 1. Demonstration of Oasis MCX SPE bioanalytical extraction performance of several TKI pharmaceuticals, requiring no optimization of the starting protocol, resulting in > 72.6% recovery and matrix effects between -3.8 and -11.0% from plasma.

Results and Discussion

High analyte recovery and selectivity from plasma using Oasis MCX SPE for top-selling small molecule tyrosine kinase inhibitor pharmaceuticals in 2023.¹

References

1. Williams, RE and Leatherwood HM Top 200 Small Molecule Drugs by Retail Sales in 2023. <https://bpb-us-e2.wpmucdn.com/sites.arizona.edu/dist/9/130/files/2024/05/2023Top200SmallMoleculePosterV5.pdf>

> (accessed 21 June 2024).

Featured Products

720008424, July 2024



© 2024 Waters Corporation. All Rights Reserved.

[使用条款](#) [隐私策略](#) [商标](#) [招聘](#) [法律和隐私声明](#) [危险化学品生产经营许可证](#) [Cookie](#) [Cookie 设置](#)

[沪ICP备06003546号-2](#) [京公网安备 31011502007476号](#)