



ACQUITY UPLC BEH AMIDE COLUMNS

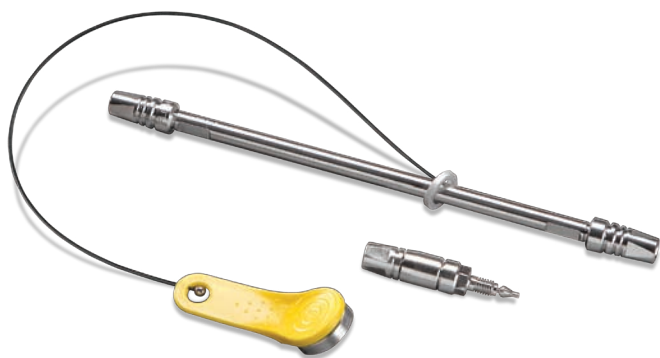

Acquity
 UltraPerformance LC®

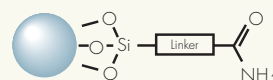

Particle Type:	Ethylene Bridged Hybrid [BEH]
Ligand Type:	Trifunctional Amide
Particle Size:	1.7 μm
Ligand Density:	7.5 $\mu\text{mol}/\text{m}^2$
Carbon Load:	12%
Endcap Style:	None
pH Range:	2-11
Pressure Tolerance:	1000 bar

[HYDROPHILIC INTERACTION CHROMATOGRAPHY]


Since 2003, Waters has developed innovative stationary phases for Hydrophilic Interaction Chromatography [HILIC] to overcome the challenge of retaining and separating extremely polar compounds. Based on Waters novel ethylene bridged hybrid [BEH] particle technology, New ACQUITY UPLC® BEH Amide columns utilize a chemically stable, trifunctionally-bonded amide phase, enabling a new dimension in stability and versatility for HILIC separations.

Designed to retain polar analytes and metabolites that are too polar to retain by reversed-phase [RP] chromatography, ACQUITY UPLC BEH Amide columns facilitate the use of a wide range of mobile phase pH [2 –11] to facilitate the exceptional retention of polar analytes spanning a wide range in polarity, structural moiety and pK_a .

In addition to enhanced retention of polar compounds, ACQUITY UPLC BEH Amide columns provide increased mass spectrometry response, direct compatibility with sample preparation eluates [PPT, LLE and SPE] as well as orthogonal selectivity compared to reversed-phase materials, making HILIC an attractive alternative separation technique.

Separation of Water Soluble Vitamins

Column:	ACQUITY UPLC BEH Amide , 1.7 μm 2.1 x 50 mm	
Part Number:	186004800	
Mobile Phase A:	50/50 ACN/H ₂ O with 10 mM CH ₃ COONH ₄ and 0.04% NH ₄ OH, pH 9.0	
Mobile Phase B:	90/10 ACN/H ₂ O with 10 mM CH ₃ COONH ₄ and 0.04% NH ₄ OH, pH 9.0	
Flow Rate:	0.5 mL/min	
Gradient:	Time (min)	Profile (%A %B)
	0.00	0.1 99.9
	3.50	70.0 30.0
	3.51	0.1 99.9
	7.50	0.1 99.9
Injection Volume:	5.0 μL (PICO)	
Sample Diluent:	75/25 ACN/MeOH with 0.2% HCOOH	
Column Temp.:	30 °C	
Weak Ndl. Wash:	ACN/H ₂ O 95/5	
Detection:	UV @ 265 nm	
Sampling Rate:	20 points/sec	
Filter Time Constant:	Normal	
Instrument:	Waters ACQUITY UPLC with ACQUITY UPLC PDA Detector	

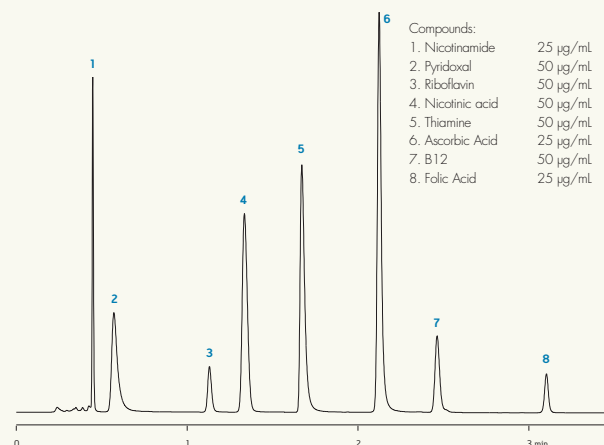

[CARBOHYDRATE ANALYSIS]

One of the most abundant and diverse classes of compounds analyzed by HILIC are carbohydrates [e.g., monosaccharides, disaccharides, oligosaccharides and polysaccharides].

The ACQUITY UPLC BEH Amide column is an exceptional tool for the analysis of carbohydrates, providing the following benefits:

- The 1.7 μm particle size enables high resolution, high speed analysis of carbohydrates in complex sample matrices while maintaining or improving chromatographic resolution.
- Increased chemical stability enables the use of high pH and high temperature to collapse reducing sugar anomers, shorten analysis times and improve MS detection.
- BEH particle technology, in combination with a trifunctionally-bonded amide phase, provides exceptional column lifetime, thus improving assay robustness.
- Unlike amine-based columns used for carbohydrate analysis, the ACQUITY UPLC BEH Amide column is not susceptible to Schiff base formation, thus improving quantitation accuracy.
- High pH mobile phase compatibility facilitates MS detection without the need for derivatization, post column addition or complexing with a metal cation, thus dramatically simplifying sample pre-treatment before LC/MS analysis and improving assay sensitivity.

ACQUITY UPLC BEH Amide Columns

Chemistry	Particle Size	Dimension	Part Number 1 pack	Part Number 3 pack	Part Number Method Validation Kit
BEH Amide	1.7 μm	1.0 x 50 mm	186004848	176001914	—
BEH Amide	1.7 μm	1.0 x 100 mm	186004849	176001915	—
BEH Amide	1.7 μm	1.0 x 150 mm	186004850	176001916	—
BEH Amide	1.7 μm	2.1 x 30 mm	186004839	176001906	—
BEH Amide	1.7 μm	2.1 x 50 mm	186004800	176001907	186004807
BEH Amide	1.7 μm	2.1 x 100 mm	186004801	176001908	186004808
BEH Amide	1.7 μm	2.1 x 150 mm	186004802	176001909	—
BEH Amide	1.7 μm	3.0 x 30 mm	186004803	176001910	—
BEH Amide	1.7 μm	3.0 x 50 mm	186004804	176001911	186004809
BEH Amide	1.7 μm	3.0 x 100 mm	186004805	176001912	186004810
BEH Amide	1.7 μm	3.0 x 150 mm	186004806	176001913	—
BEH Amide	1.7 μm	2.1 x 5 mm*	—	186004799	—


* VanGuard Pre-column 3-pack

UPLC/MS Analysis of Saccharides

Column: **ACQUITY UPLC BEH Amide**, 1.7 μm 2.1 x 50 mm
 Part Number: 186004800
 Mobile Phase A: 30/70 MeCN/H₂O with 0.10% NH₄OH
 Mobile Phase B: 80/20 MeCN/H₂O with 0.10% NH₄OH
 Flow Rate: 0.17 mL/min
 Gradient:

Time (min)	%A	%B
Initial	0	100
5.00	60	40
5.01	0	100

Injection Volume: 0.7 μL (PICO)
 Sample Diluent: 50/50 MeCN/H₂O
 Column Temp.: 35 °C
 Needle Wash: ACN/H₂O 75/25
 Instrument: Waters ACQUITY UPLC with ACQUITY UPLC TGD
 Ionization Mode: ES-
 Capillary: 2.8 kV
 Cone Voltage: 25 V
 Source Temp.: 120 °C
 Desolv. Temp.: 350 °C
 Desolv. Gas Flow: 500 L/Hr
 S/R m/z: 179.2 [fructose, glucose]; 341.3 [sucrose, maltose]; 503.4, 665.5, 827.6, 989.7, 1115 [maltooligosaccharides (n=1 to 5)]
 Dwell Time: 0.08 s

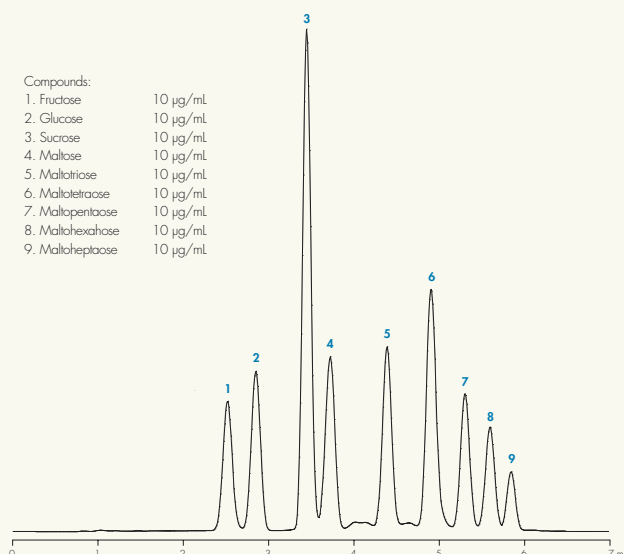

Waters

THE SCIENCE OF WHAT'S POSSIBLE.™



© 2009 Waters Corporation. Waters, The Science of What's Possible, ACQUITY UltraPerformance LC and ACQUITY UPLC are trademarks of Waters Corporation. All other trademarks are the property of their respective owners.

October 2009 720003122EN IH-PDF

Waters Corporation
 34 Maple Street
 Milford, MA 01757 U.S.A.
 T: 1 508 478 2000
 F: 1 508 872 1990
www.waters.com